

Oct 20, 2017

Version 1

0.5 YTSS Agar Plates (1.5%) V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.ehnbb5e>



Ben Temperton¹

¹University of Exeter

Temperton Lab



Ben Temperton

University of Exeter

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ehnbb5e>

Protocol Citation: Ben Temperton 2017. 0.5 YTSS Agar Plates (1.5%). **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.ehnbb5e>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: February 04, 2016

Last Modified: March 23, 2018

Protocol Integer ID: 2318

Keywords: ytss agar plate



- 1 16 g of Sigma Sea Salts
3.2 g of yeast extract
2 g of tryptone
12 g (1.5%) of bacto agar
800 mL MilliQ
- 2 Autoclave for 30 mins @ 121 °C
Allow to cool to ~ 60 °C
- 3 Pour into plates in a flow hood and leave for 30 mins to solidify
- 4 Stack plates and wrap them - store at RT or 4 °C