

Apr 06, 2019 Version 1

1. RNA isolation for tissue V.1



Forked from [1. RNA isolation for tissue](#)

DOI

dx.doi.org/10.17504/protocols.io.zu7f6zn

Sze-Xian Lim¹, Chin Yee Tan¹

¹Duke University



Sze-Xian Lim

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.zu7f6zn

Protocol Citation: Sze-Xian Lim, Chin Yee Tan 2019. 1. RNA isolation for tissue. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.zu7f6zn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: April 06, 2019

Last Modified: April 06, 2019

Protocol Integer ID: 22143



- 1 Homogenization in TriZol
Add 1mL of Trizol reagent per 30mg of tissue and homogenize using handheld homogenizer
- 2 Incubate at RT for 5 mins to allow nucleoprotein complexes to dissociate
- 3 Add 1/5 the volume of trizol of chloroform carefully, and vortex to mix well
- 4 Spin down at max speed, chilled centrifuge for 15 minutes
- 5 Carefully remove the top aqueous phase and transfer to a new Eppendorf tube
The interphase and bottom organic phase can be saved for DNA and protein respectively
- 6 To the aqueous phase, add 500 microliters of 100% isopropanol, mix by inversion and incubate at -20C for a minimum for 2hrs
- 7 Spin down at max speed for 30 minutes to precipitate RNA
- 8 Remove supernatant, and add 1mL of 75% Ethanol to wash the pellet
- 9 Spin down at max speed for 15 minutes and remove supernatant
- 10 Resuspend pellet in appropriate volume of nuclease free water