



Nov 08, 2019 Version 3

🌐 10X Genomics Single-Nucleus RNA-Sequencing for Transcriptomic Profiling of Adult Human Tissues V.3

DOI

dx.doi.org/10.17504/protocols.io.86khzcw

Sarah Urata¹, Blue Lake¹, Dinh Diep¹, Masato Hoshi², Sanjay Jain³, Kun Zhang¹

¹University of California, San Diego; ²Washington University, Saint Louis;

³Washington University School of Medicine, Saint Louis

KPMP

Human BioMolecular Atl...



Blue Lake

UCSD

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.86khzcw

External link: <http://genome-tech.ucsd.edu/ZhangLab/>

Protocol Citation: Sarah Urata, Blue Lake, Dinh Diep, Masato Hoshi, Sanjay Jain, Kun Zhang 2019. 10X Genomics Single-Nucleus RNA-Sequencing for Transcriptomic Profiling of Adult Human Tissues. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.86khzcw>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 08, 2019

Last Modified: November 08, 2019

Protocol Integer ID: 29612

Keywords: Kidney, KPMP, Single Nucleus, RNA Sequencing, 10X Genomics



Abstract

10X Genomics Single Cell 3' (v3) RNA sequencing is a microdroplet-based method that permits the effective capture and sequencing of the mRNA and pre-mRNA molecules from single nuclei [1]. RNA molecules are transcribed and processed within the nucleus before exporting to ER for translation into proteins. As such, nuclear RNA is a mixture of nascent transcripts, partially or fully processed mRNA, and various non-coding RNA molecules. The total RNA content within the nucleus is roughly 10% of the RNA content in a whole cell, but has been found to accurately represent whole cell expression values in adult human tissues [2,3] including the kidney [4]. Nuclei can be readily isolated from frozen tissues with a combination of chemical and physical treatments that can effectively circumvent the non-uniform or incomplete dissociation of solid tissues into single cells, as well as RNA degradation or artefacts (such as stress response) during dissociation. Here we present a modified version of the published 10X protocol [1] that we have adapted for the processing of adult human kidney nuclei.

References

1. Chromium Single Cell 3' Reagent Kits v3 User Guide (Rev A) CG000183, support.10xgenomics.com.
2. Lake et al. (2016). Science, doi:10.1126/science.aaf1204.
3. Lake et al. (2018). Nature Biotechnology, doi:10.1038/nbt.4038.
4. Lake et al. (2019). Nature Communications, doi:10.1038/s41467-019-10861-2.

Guidelines

Full protocol is from 10X Genomics. All modifications are to the original protocol (Chromium Single Cell 3' Reagent Kits v3 User Guide (Rev A) CG000183, support.10xgenomics.com)



Materials

MATERIALS

⊗ Chromium™ Single Cell 3 GEM Library & Gel Bead Kit v3 **10x Genomics Catalog #1000075**

⊗ Chromium Chip B Single Cell Kit **10x Genomics Catalog #1000074**

⊗ Chromium i7 Multiplex Kit **10x Genomics Catalog #120262**

Nuclease-free water

Ethanol (200 proof)

Dynabeads MyOne Silane

SPRI select reagent set

50% glycerol

10% Tween 20

Low TE Buffer (10mM Tris-HCl, pH 8.0, 0.1 EDTA)

Qiagen buffer EB

Qubit dsDNA HS Assay Kit




PCR strip tubes with flat cap

LoBind 1.5 ml tubes

Qubit Assay tubes



Isolate Nuclei

- 1 Prepare nuclei according to the protocol "Isolation of single nuclei from solid tissues" steps 1-14. [dx.doi.org/10.17504/protocols.io.ufketkw](https://doi.org/10.17504/protocols.io.ufketkw)
- 2 Resuspend nuclei in  100 μL to  1 mL of PBS + 0.1% RNase Inhibitor (volume depends on target concentration)
- 3 Count nuclei (e.g. BioRad T20 Cell Counter)
- 4 Check nuclei integrity under fluorescent microscope using DAPI channel. Nuclei should appear distinct, have rounded borders and the majority occurring as singlets. High clumping rates would indicate damaged nuclei and would require re-filtering using 30- μm CellTrics filter or exclusion.
- 5 Dilute nuclei stock to be in the 700-1200 nuclei per μL range in  50 μL total volume using the dilution guide

Note

If nuclei concentrations fall below target range, then centrifugation can be performed to increase nuclei concentrations to be within range


- Start with nuclei stock in PBS with 0.1% BSA
- Spin 900xg for 10 minutes
- Resuspend in appropriate volume PBS
- Use 10 μL to count; 50 μL for loading
-

QC cutoff: minimum of 10,000 nuclei (for 200 nuclei per μL in 50 μL total volume)

Safety information


Caution: BSA is necessary to avoid clumping and prevent nuclei loss from sticking to the tube during the spin down. DO NOT USE more than .01% BSA in nuclei stock

GEM Generation and Barcoding











- 6 Prepare Reagents for use
 1. Equilibrate gel beads to room temperature for  00:30:00 .

2. Thaw, vortex, and centrifuge RT Reagent, Template Switch Oligo, and Reducing Agent B. Verify no precipitate in Reducing Agent B.
3. Centrifuge RT Enzyme C before adding to Master Mix.

7 Prepare Master Mix

1. Prepare on ice. Pipette mix 15x and centrifuge briefly.
2. Add  33.4 μL Master Mix to each tube of a PCR 8-tube strip on ice.

8 Load Chromium Single Cell B Chip

1. Assemble Chromium Chip B in a 10X Chip Holder.
2. Must load chip in order according to row label for microfluidic channels to work properly: Row 1 \rightarrow Row 2 \rightarrow Row 3
3. Make sure NO bubbles are introduced while loading chip.
4. Dispense 50% glycerol solution into unused Chip Wells (if <8 samples used per chip)
 - a.  75 μL into unused wells in Row 1
 - b.  40 μL into unused wells in Row 2
 - c.  280 μL into unused wells in Row 3
5. Use Cell Suspension Volume Calculator Table to add the appropriate volume of nuclease-free water and corresponding volume of nuclei stock to make working nuclei suspension
6. Nuclei Loading
 - a. Minimum: load 800 nuclei \rightarrow target recovery 500 nuclei
 - b. Maximum: load 1600 nuclei \rightarrow target recovery 10000 nuclei
7. Add  46.6 μL working Nuclei Suspension to the Master Mix in PCR tubes on ice for  80 μL total volume
8. Load  75 μL Master Mix and Nuclei Suspension into the bottom center of each well in Row 1 without introducing bubbles.
9. Snap the Gel Bead strip into a 10X Vortex Adapter. Vortex  00:00:30 . Remove the Gel Bead strip and flick sharply downward to ensure maximum recovery. Confirm there are no bubbles at the bottom of the tubes and liquid levels look even.
10. Load  40 μL gel beads in Row 2. Only puncture the foil seal for gel bead tubes being used. Dispense slowly and without introducing bubbles.
11. Load  280 μL partitioning oil into each Row 3 by pipetting two aliquots of  140 μL .
12. Attach 10X gasket. Align the top-notch. Ensure gasket holes are aligned with the wells. Avoid touching the smooth gasket surface. Do not press down on the gasket.

9 Run the Chromium Controller

1. Press the eject button on the controller to eject.
2. Place the assembled chip with the gasket in the tray. Press the button to retract the tray.
3. Confirm the program on screen. Press the play button.
4. At completion of the run (~ 00:07:00), the controller will chime. Proceed immediately to the next step.

10 Transfer GEMs

1. Chill strip tubes on ice.
2. Press the eject button to remove the chip.
3. Discard the gasket. Open the chip holder. Fold the lid back until it clicks to expose the wells at 45 degrees.
4. Check the volume in rows 1-3. Abnormally high volume in any well indicates a clog.
5. Slowly aspirate 100 uL GEMs from the lowest points of the Recovery Wells without creating a seal between the pipette tips and the wells.

Note

GEMs should appear opaque and uniform across all channels. Excess partitioning oil (clear) in the pipette tips indicates a potential clog.

6. Slowly dispense (~ 00:00:20) GEMs into the tube strip on ice with the pipette tips against the sidewalls of the wells.

11 GEM-RT Incubation





1. Incubate in a thermocycler to complete reverse transcription.
2. Store at 4 °C for up to 72:00:00 or at -20 °C for up to a week. Or proceed to the next step.

12 Post GEM-RT Cleanup & cDNA Amplification - Prepare Reagents for use



1. Thaw, vortex, and centrifuge a tube of Reducing Agent B and cDNA primers.
2. Maintain Amp Mix on ice after vortex and centrifuge
3. Thaw Cleanup Buffer at 65 °C for 00:10:00 with shaking at max rpm then cool to room temperature.

13 Post GEM RT-Cleanup --Dynabead

1. Add 125 µL Recovery Agent to each sample at room temperature. DO NOT MIX.
Wait 00:01:00 .

2. Carefully aspirate  125 μL of pink oil phase from the bottom of the tube. DO NOT aspirate any aqueous sample.
3. Prepare Dynabeads Cleanup Mix
4. Vortex and centrifuge mixture and add  200 μL to each sample. Pipette 5x to mix.
5. Incubate at room temperature for  00:10:00 ; pipette mix again  00:05:00 into incubation.
6. Prepare Elution Solution I. Vortex and centrifuge briefly.
7. Place sample on magnetic separator and remove the supernatant.
8. Wash twice with freshly prepared 80% ethanol.
9. Elute with 35.5 μL Elution Solution I.
10. Transfer 35 μL to a new strip tube.






14 cDNA Amplification

1. Prepare cDNA Amplification Mix on ice. Vortex and centrifuge briefly.
2. Add  65 μL cDNA Amplification Reaction Mix to  35 μL sample.
3. Pipette mix 15x to mix and centrifuge briefly.
4. Incubate in a thermocycler to amplify cDNA.

Note

- Use cycle number optimization table for total number of cycles
- When using nuclei, increase the optimized number of cycles by one cycle
- Maximum: load 1600 nuclei - perform 12 cycles

15 cDNA Cleanup --SPRIselect

1. Perform a 0.6X beads purification with SPRIselect reagent
2. Wash twice with freshly prepared 80% ethanol
3. Elute with  40.5 μL Qiagen Buffer EB.
4. Transfer  40 μL to a new strip tube.
5. Store at  4 $^{\circ}\text{C}$ for up to  72:00:00 or at  -20 $^{\circ}\text{C}$ for up to 4 weeks. Or proceed to the next step.



16 cDNA QC & Quantification

1. Run TapeStation to obtain concentration and size
2. Expected size range for amplified cDNA: 400 to 2500 bp
3. cDNA total yield range: 80 to 1000 ng

Note


QC Cutoff: Minimum cDNA total yield of 80 ng

3' Gene Expression Library Construction

17 Prepare Reagents for use



1. Thaw, vortex, and centrifuge Fragmentation Buffer, Adaptor Oligos, Ligation Buffer, SI Primer.
2. Maintain on ice Fragmentation Enzyme, DNA Ligase, and Amp Mix.

18 Fragmentation, End Repair & A-tailing



1. Prepare Fragmentation Mix on ice. Pipette mix and centrifuge briefly.
2. Transfer  10 μ L purified cDNA to a new strip tube





Note

Note that 10 μ L (25%) cDNA sample is sufficient for generating 3' Gene Expression Library. The remaining 30 μ L (75%) cDNA sample can be stored at 4 C for up to 72 hours or at -20 C for up to 4 weeks for generating additional 3' Gene Expression Libraries.


3. Add  25 μ L Buffer EB to each sample
4. Add  15 μ L Fragmentation Mix to each sample
5. Pipette mix, centrifuge briefly, and transfer to pre-cooled block. "SKIP" hold step to initiate the protocol

19 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection --SPRIselect



1. Vortex to resuspend the SPRIselect reagent. Add  30 μ L SPRIselect reagent (0.6X) to each sample and pipette mix 15x.
2. Incubate  00:05:00 at room temperature.
3. Centrifuge briefly
4. Place sample on magnetic separator. DO NOT discard supernatant.

5. Transfer  75 μL supernatant to a new tube strip tube.
6. Vortex to resuspend the SPRIselect reagent. Add  10 μL SPRIselect reagent (0.8X) to each sample and pipette mix 15x.
7. Place sample on magnetic separator and remove supernatant. DO NOT discard any beads.
8. Wash twice with freshly prepared 80% ethanol
9. Elute with  50.5 μL Qiagen Buffer EB.
10. Transfer  50 μL to a new strip tube.



20 Adaptor Ligation

1. Prepare Adaptor Ligation Mix on ice. Pipette mix and centrifuge briefly.
2. Add  50 μL Adaptor Ligation Mix to sample. Pipette mix 15x. Centrifuge briefly.
3. Incubate in a thermocycler to ligate adaptor.

21 Post Ligation Cleanup --SPRIselect

1. Perform a 0.8X beads purification with SPRIselect reagent
2. Wash twice with freshly prepared 80% ethanol
3. Elute with  30.5 μL Qiagen Buffer EB.
4. Transfer  30 μL to a new strip tube.

22 Sample Index PCR

1. Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run.
2. Prepare Sample Index PCR Mix on ice. Pipette mix and centrifuge briefly.
3. Add  60 μL Sample Index PCR Mix to the sample
4. Add  10 μL of individual Chromium i7 Sample Index (from SI-GA plate) to each well. Record well assignment. Mix and centrifuge briefly.
5. Incubate in a thermocycler to incorporate sample indices.

Note

Use cycle number recommendation table for total number of cycles; recommended number of cycles is based on cDNA input



6. Store at 4 °C for up to 72:00:00 . Or proceed to the next step.

23 Post Sample Index PCR Double Sided Size Selection --SPRIselect

1. Vortex to resuspend the SPRIselect reagent. Add 60 µL SPRIselect reagent (0.6X) to each sample and pipette mix 15x.
2. Incubate 00:05:00 at room temperature.
3. Centrifuge briefly
4. Place sample on magnetic separator. DO NOT discard supernatant.
5. Transfer 150 µL supernatant to a new tube strip tube.
6. Vortex to resuspend the SPRIselect reagent. Add 20 µL SPRIselect reagent (0.8X) to each sample and pipette mix 15x.
7. Incubate 00:05:00 at room temperature.
8. Centrifuge briefly
9. Place sample on magnetic separator and remove supernatant. DO NOT discard any beads.
10. Wash twice with freshly prepared 80% ethanol xi. Elute with 35.5 µL Qiagen Buffer EB.
11. Transfer 35 µL to a new strip tube.
12. Store at 4 °C for up to 72:00:00 or at -20 °C for long-term storage.

Note

QC Post Library Construction:

- Quantify Library (e.g. using Qubit dsDNA HS Assay)
- Estimate Library size range (e.g. using TapeStation or BioAnalyzer) - expected size range of 300 to 800 bp, average: 475 bp

Sequencing

24 MiSeq Sequencing - **QC for estimation of library quality and number of nuclei captured**

1. Paired End, Single Indexing
 - a. Read 1: 28 cycles
 - b. i7 Index: 8 cycles
 - c. i5 Index: 0 cycles
 - d. Read 2: 91 cycles
2. Library Loading
 - a. 10X recommended Loading concentration: 11 pM
 - b. Optional: 1% PhiX
3. Output
 - a. 22-25 million reads

25 NovaSeq Sequencing (target - 25,000-50,000 reads per nucleus)


1. Paired End, Single Indexing
 - a. Read 1: 28 cycles
 - b. i7 Index: 8 cycles
 - c. i5 Index: 0 cycles
 - d. Read 2: 91 cycles
2. Library Loading
 - a. 10X recommended Loading concentration: 300 pM
 - b. Optional: 1% PhiX
3. Output
 - a. SP: 650–800 million reads
 - b. S1: 1.3–1.6 billion reads
 - c. S2: 3.3 –4.1 billion reads
 - d. S4: 8-10 billion reads

Cell Ranger Mapping and Analysis Pipeline

26 Generate pre-mRNA reference

Note

IMPORTANT: for nuclei experiments, we need to use the reference files with pre-mrna in order to count intronic reads in the UMI counts

 Capture.PNG

- b. Generate Sample Sheet




Note

Use sample sheet generator provided by 10X Genomics to generate a "SampleSheet.csv": **Sample Sheet Generator**


c. Generate fastq files

Note

Use **mkfastq** command

 Capture.PNG

d. Run count for each sample

 Capture.PNG