



Mar 26, 2025

Version 2

10x Protocols: Chromium Next GEM Single Cell 5' -- University of Minnesota TMCs (CG000331 Rev F) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvokn89l4o/v2>

Laura J Niedernhofer¹, David A Bernlohr¹

¹University of Minnesota, Minneapolis, MN

Cellular Senescence Net...



Allie Abdellatif

MIBAM, University of Minnesota, Minneapolis, MN

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvokn89l4o/v2>

Protocol Citation: Laura J Niedernhofer, David A Bernlohr 2025. 10x Protocols: Chromium Next GEM Single Cell 5' -- University of Minnesota TMCs (CG000331 Rev F). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5qpvokn89l4o/v2> Version created by [Allie Abdellatif](#)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 25, 2025

Last Modified: March 26, 2025

Protocol Integer ID: 125027

Keywords: minnesota genomics center, 10x genomic, chromium next gem single cell, barcoded cdna, cdna, university of minnesota tmc, nuclei isolation, cell, minnesota tmc, following protocol, protocol, nuclei, tmc, microfluidic, cell dissociation, university of minnesota genomics center, 10x protocol

Funders Acknowledgements:

NIH

Grant ID: 5U54AG076041-03

NIH

Grant ID: 5U54AG079754-02

Abstract

Following single-cell dissociation or nuclei isolation, the Next GEM 5' assay uses microfluidics to partition and assign cell- or nuclei-specific barcodes to transcript cDNA at the 5' end. Barcoded cDNA is prepared with adaptors for sequencing by synthesis (SBS). The following protocol has been used at the University of Minnesota TMCs in collaboration with the University of Minnesota Genomics Center. The following protocol has been adapted from protocols developed by 10x Genomics and Illumina to be used at the University of Minnesota TMCs in collaboration with the University of Minnesota Genomics Center. These protocols are owned by their respective companies and are subject to periodic revision.



Tissue Preparation

- 1 Complete single cell or nuclei isolation prior to starting this protocol

Library Preparation

- 2  CG000331_Chromium_Next_GEM_S... 4.7MB

Note

Sequence with the read format 29,10,10,89

Note

Sequencers used at UMN Genomics Center:

- Illumina NextSeq 2000
- Illumina NovaSeq 6000
- Illumina NovaSeq X Plus

FASTQ Generation

- 3 BCL data from Illumina sequencer is demultiplexed and converted into FASTQ format using bcl2fastq version 2.20.0