

May 17, 2024

2D differentiation of NPCs to neurons & astrocytes

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzjy9rlx1/v1>

Jessie Buth¹

¹UCLA



Jessie Buth

UCLA

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzjy9rlx1/v1>

Protocol Citation: Jessie Buth 2024. 2D differentiation of NPCs to neurons & astrocytes . protocols.io

<https://dx.doi.org/10.17504/protocols.io.rm7vzjy9rlx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: May 09, 2024



Last Modified: May 17, 2024

Protocol Integer ID: 99474

Keywords: astrocytes in 2d culture, 2d differentiation of npc, neural progenitor cell, astrocyte, 2d differentiation, neuron, cell, 2d culture, differentiation, npc

Abstract

Protocol to differentiate neural progenitor cells to mature neurons and astrocytes in 2D culture.

Protocol materials

 Corning® Matrigel® Basement Membrane Matrix, LDEV-free, 10 mL **Corning Catalog #354234**

 Corning® Matrigel® Basement Membrane Matrix, Phenol Red-free, LDEV-free, 10 mL **Corning Catalog #356237**



Overview

1 Typical workflow:

d(-14) thaw the NPCs (in ROCKi)

- maintain in +EGF+FGF

d(-7) passage (in ROCKi)

- maintain in +EGF+FGF

d0 passage (in ROCKi)

d1 full media change to start differentiation to neurons/astrocytes

d7 passage (without ROCKi)

d14 passage (without ROCKi)

d42 Mitomycin C treatment (**for neurons only**)

~d60-100 replate neurons/astrocytes for experiments

Note

If more cells are needed, the NPCs can be maintained additional passages in +EGF+FGF before differentiation, but can't be refrozen and thawed again. If frozen/thawed again, they don't proliferate or differentiate well.

Media Recipe

2 **NPC Base Media (500mL):**

240mL 50% DMEM/F12 (Hyclone VWR Cat # SH30023.02)

240mL 50% Neurobasal (FisherSci Cat # 21103-049)

2.5mL 0.5X N2 (100X; Invitrogen Cat # 17502048)

5mL 0.5X B27 without VA (50X; Invitrogen, 12587010)

5mL 1X Glutamax (100x Invitrogen Cat # 35050061)

5mL 1X NEAA (Invitrogen Cat # 11140050)

*250 uL 5 ug/mL Human Insulin Solution (250mg per 500mL) (Sigma Cat # I9278-5ML)

1mL 1X Primocin (stock 500X, InvivoGen ant-pm-2)

50uL 1µg/mL Heparin (10 mg/ml stock Sigma H3419-100KU)

**Note**

*Look up the lot number of human insulin & make aliquots of 2.5mg.
The concentration is typically 9-11 mg/mL -> 2.5 mg will be ~250uL.

For NPC maintenance:

+20 ng/mL bFGF (stock 100 ug/mL): 10uL/50mL NPC base

+20 ng/mL EGF (stock 200 ug/mL): 5uL/50mL NPC base

For Neuron differentiation:

d0-d80+

+20 ng/mL BDNF (stock 200 ug/mL): 5uL/50mL NPC base

+20 ng/mL GDNF (stock 100 ug/mL): 10uL/50mL NPC base

+1 uM cAMP (stock 10mM): 5uL/50mL NPC base

~d42 (can be d42-d49)

1hr pulse with 5 ug/mL Mitomycin C ready made solution (stock 10 mg/mL Sigma Cat# M5353)

^^ this is enrich for neurons, dividing cells die off or stop proliferating over the next 4-7 days

For Astrocyte differentiation:

d0-d14

+10 ng/mL hLIF (stock 100 ug/mL): 5uL/50mL

+10 ng/mL EGF (stock 200 ug/mL): 2.5uL/50mL

d14-d80+

+10 ng/mL hLIF (stock 100 ug/mL): 5uL/50mL

+10 ng/mL CNTF (stock 100 ug/mL): 5uL/50mL

Thawing Neural Progenitor Cells (NPCs)

- 3 See protocol for thawing hematopoietic progenitor cells (HPCs).
<https://www.protocols.io/view/thawing-frozen-hematopoietic-stem-cells-hpcs-dddt226n>

It is the same procedure for NPCs, except:

- the plates are coated with **0.5 mg/mL standard matrigel**
- the cells are maintained in **NPC base +EGF+FGF+10uM ROCKi** (ROCKi is removed 24hr after thawing)



- the cells are plated at **250-300K/well** in a **6w plate**

Passaging NPCs (every 6-7 days - volumes listed at for 1w of a 6w plate)

- 4 Before starting, coat 6-well plates with 0.5 mg/mL standard matrigel (Set aside for later)

30m

Note

Use standard matrigel (regular or phenol-free):

 Corning® Matrigel® Basement Membrane Matrix, LDEV-free, 10 mL **Corning Catalog #354234**

 Corning® Matrigel® Basement Membrane Matrix, Phenol Red-free, LDEV-free, 10 mL **Corning Catalog #356237**

To make aliquots of matrigel:

- The concentration of a lot of matrigel can be looked up here in "quality certificate lookup" section. (<https://www.corning.com/worldwide/en/products/life-sciences/resource-library.html>).
- To keep coating consistent across lots, I check the concentration and make aliquots. However, if you don't have the lot number information, 0.5mg/mL is close to a 1:20 dilution.

To coat plates:

- Thaw an aliquot of matrigel at  4 °C or  On ice
- Dilute with the appropriate amount of ice cold DMEM/F12
 -  6.5 mL for 6.5-wells = 3.25 mg matrigel
- Aliquot  1 mL per well in a cold 6-well plate
- Shake/tilt the plate until the matrigel evenly coats the bottom
- Incubate  37 °C for  00:30:00
- After incubating, aspirate the matrigel
- Add  2 mL per well DMEM/F12
- Place the plate in the incubator at  37 °C

Note

Plates can be stored in the incubator with DMEM/F12 for up to 1 week before use

- 4.1 Pre-treat the cells with **[1M] 10 micromolar (μM)** ROCK inhibitor (ROCKi)

1h 30m



- Pick up the media in the well to a 15mL conical tube.
- Add [M] 10 micromolar (μM) ROCKi ( 2 μL per well) and mix well.
- Gently return the media to the dish.
- Incubate the cells with ROCKi for  00:30:00 -  01:00:00 at  37 °C

4.2 Lift the cells with Accutase.

- Aspirate the media.
- Wash with 2mL/well PBS (without Ca/Mg).
- Aspirate the PBS.
- Add 1mL/well Accutase.
- Incubate 5 minutes at 37C.
- Pick up the cells in accutase with a glass pipette and blow the cells off the dish.
- Transfer the cells to a 15mL conical tube.
- Add 6mL NPC base media to the dish and blow the remaining cells off the dish.
- Transfer the remaining cells to the 15mL conical tube.
- Pipette up/down 3-5x in the conical tube to break apart clumps. To break apart clumps, pipette the cells up, and then press the pipette against the side of the tube and blow out quickly.
- Pick up the entire cell mixture and filter through a 100um cell strainer over a 50mL conical tube to remove any remaining large clumps.
- Transfer the cell mixture back to the 15mL conical tube.
- Centrifuge at 180g for 5 minutes room temperature.
- Aspirate the media.
- Close the conical tube lid and tap the pellet to loosen the cell pellet.
- Resuspend in 2-3mL of NPC base media.
- Pipette up/down 3-5x or until the cell solution is cloudy.

4.3 Count the cells and aliquot to new plates.

- Count 10uL cells + 10uL trypan blue with a hemocytometer.
- Determine the #uL for 1 million cells.
- Add 1 million cells/well in  2 mL per well NPC+EGF+FGF+ROCKi in new 0.5 mg/mL matrigel coated plates.

NPC+EGF+FGF+ROCKI (10mL):

10mL NPC base media

1uL 20ng/mL EGF (stock 200ug/mL)

2uL 20ng/mL FGF (stock 100ug/mL)

10uL 10uM ROCKi (stock 10mM)

- Gently place the plate in the incubator and shake evenly (4x forward/back, 4x side-to-side, 4x forward/back).



Note

If the cells were confluent, 1 million cells/well is typically around a 1:2-1:4 passage dilution.

Expansion of NPCs

- 5 d1 after passaging NPCs (see STEP 3)
- Perform a full media change and replace with NPC maintenance media without ROCKi (NPC base +EGF+FGF)

NPC+EGF+FGF (10mL):

10mL NPC base media

1uL 20ng/mL EGF (stock 200ug/mL)

2uL 20ng/mL FGF (stock 100ug/mL)

d3

- Perform a half media change. Remove 900uL/well, and add a fresh 1mL/well NPC base +EGF+FGF.

d5

- Perform a half media change. Remove 900uL/well, and add a fresh 1mL/well NPC base +EGF+FGF.

d6-7

- Passage to new dishes coated with 0.5mg/mL standard matrigel in NPC base +EGF+FGF.
 - for 6w plates → 1 million cells/well in 2mL NPC base +EGF+FGF
 - for 5cm cellBIND dishes → 2.4 million cells/dish in 6mL NPC base +EGF+FGF

Note

The high plating density is important, the cells need to be very dense to maintain progenitor identity.

I've stained the cells after maintaining in +EGF+FGF at high density up to 5x passages after thawing and > 90% of the cells remain PAX6+SOX2+NESTIN+. If passaged at lower density, the cells will stop dividing and/or differentiate.

Differentiation of NPCs to Neurons

- 6 d1 after passaging NPCs (see STEP 3)
- Perform a full media change and replace with neuron differentiation media (NPC base +BDNF+GDNF+cAMP)
NPC+BDNF+GDNF+cAMP (10mL):
10mL NPC base media
1uL 20 ng/mL BDNF (stock 200 ug/mL)
2uL 20 ng/mL GDNF (stock 100 ug/mL)
1uL 1 uM cAMP (stock 10mM)
- d3
- Perform a half media change. Remove 900uL/well, add a fresh 1mL/well NPC base +BDNF+GDNF+cAMP.
- d5
- Perform a half media change. Remove 900uL/well, add a fresh 1mL/well NPC base +BDNF+GDNF+cAMP.
- d7 [*differentiation passage 1 - no ROCKi during this step*]
- Passage to new dishes coated with 0.5mg/mL standard matrigel in NPC base +BDNF+GDNF+cAMP.
for 6w plates → 1 million cells/well in 2mL media
for 5cm cellBIND → 2.4 million cells/dish in 6mL media
for 10cm cellBIND dishes → 5.8 million cells/dish in 12mL media
- d9
- Perform a full media change and replace with neuron differentiation media (NPC base +BDNF+GDNF+cAMP) to remove any dead cells.
- d11
- Perform a half media change. Remove 900uL/well, add a fresh 1mL/well NPC base +BDNF+GDNF+cAMP.
- d13 [*differentiation passage 2 - no ROCKi during this step*]
- Passage to new dishes coated with 0.5mg/mL standard matrigel in NPC base +BDNF+GDNF+cAMP.
for 6w plates → 400K cells/well in 2mL media
for 5cm cellBIND → 1 million cells/dish in 6mL media
for 10cm cellBIND dishes → 2.5 million cells/dish in 12mL media
- d16+
- Perform a half media change every 3-4 days until 90-95% confluent. When the cells are 90-95% confluent, passage at the same density as d13 and continue with half media changes every 3-4 days until 90-95% confluent again.



for 5cm cellBIND → remove ~2-2.5mL (depending how much media evaporated, leave enough media to cover the bottom), and add a fresh 3mL/well NPC base +BDNF+GDNF+cAMP.
for 10cm cellBIND → remove ~4.5-5mL (depending how much media evaporated, leave enough media to cover the bottom), and add a fresh 6mL/well NPC base +BDNF+GDNF+cAMP.

Note

Seeding at the lower density and feeding the cells less often (every 4 days) results in more progenitor differentiation, plan to have around the number of dishes you want to end with after the 2nd differentiation passage.

~d42

- To enrich for neurons, around d42 (or close to this date when the dishes are 95% confluent again), treat the cells with 5ug/mL Mitomycin C for 1hour at 37C.
- After 1hr, perform 3x washes with 6mL PBS, and then replace with 6mL (5cm dishes) or 12mL (10cm dish) of fresh NPC base +BDNF+GDNF+cAMP.

Note

If the cells were ~95% confluent before Mitomycin C treatment, 1x 10cm typically yields around 3-6 million cells.

Continue performing half media changes until used for experiments. After Mitomycin C treatment, the cells should not expand further, so they likely won't need to be passaged again until plating for experiments.

I've been using these for experiments ~d60-100.

Differentiation of NPCs to Astrocytes

- 7 d1 after passaging NPCs (see STEP 3)
- Perform a full media change and replace with astrocyte differentiation media #1 (NPC base +hLIF+EGF)
NPC+hLIF+EGF (10mL):
10mL NPC base media
1uL 10 ng/mL hLIF (stock 100 ug/mL)
0.5uL 10 ng/mL EGF (stock 200 ug/mL)

d3

- Perform a half media change. Remove 900uL/well, and add a fresh 1mL/well NPC base +hLIF+EGF.

d5

- Perform a half media change. Remove 900uL/well, and add a fresh 1mL/well NPC base +hLIF+EGF.

d6-7 [*differentiation passage 1 - no ROCKi during this step*]

- Passage to new dishes coated with 0.5mg/mL standard matrigel in NPC base +hLIF+EGF.

for 6w plates → 1 million cells/well in 2mL media

for 5cm cellBIND → 2.4 million cells/dish in 6mL media

for 10cm cellBIND dishes → 5.8 million cells/dish in 12mL media

~d9

- Perform a full media change and replace with neuron differentiation media (NPC base +hLIF+EGF) to remove any dead cells.

~d11

- Perform a half media change. Remove 900uL/well, and add a fresh 1mL/well NPC base +hLIF+EGF.

~d13-14 [*differentiation passage 2 - no ROCKi during this step*]

- Passage to new dishes coated with 0.5mg/mL standard matrigel in **NPC base +hLIF+CNTF**.

for 6w plates → 400K cells/well in 2mL media

for 5cm cellBIND → 1 million cells/dish in 6mL media

for 10cm cellBIND dishes → 2.5 million cells/dish in 12mL media

NPC+hLIF+CNTF (10mL):

10mL NPC base media

1uL 10 ng/mL hLIF (stock 100 ug/mL)

1uL 10 ng/mL CNTF (stock 100 ug/mL)

~d16+

- Perform a half media change every 3-4 days with **NPC base +hLIF+CNTF** until 95-100% confluent.
- When the cells are 95-100% confluent, passage at the same density as ~d13 and continue with half media changes with **NPC base +hLIF+CNTF** every 3-4 days until 95-100% confluent again.

for 5cm cellBIND → remove ~2-2.5mL (depending how much media evaporated, leave enough media

to cover the bottom), and add a fresh 3mL/well **NPC base +hLIF+CNTF**.



for 10cm cellBIND → remove ~4.5-5mL (depending how much media evaporated, leave enough media to cover the bottom), and add a fresh 6mL/well **NPC base +hLIF+CNTF**.

Note

If the cells were ~95% confluent before passaging, 1x 10cm typically yields around 9-14 million cells.

The cells will expand slower with each passage, avoid passaging until 95-100% confluent. Continue performing half media changes until used for experiments, around d50-60 astrocytes with more complex branching should appear and increase in number.

I've been using these for experiments ~d60-100.

After replating for experiments or if maintained past d100, I change the media to NPC +BDNF+GDNF+cAMP to prevent the astrocytes from overgrowing and detaching from the plate.