

May 14, 2024

Version 3

## 4% PFA + 2.5% Glutaraldehyde Fixative V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.dm6gpr15dvzp/v3>

Allen Institute for Brain Science<sup>1</sup>

<sup>1</sup>Allen Institute

Allen Institute for Brain Science  
Tech. support  
[Click here to message tech. support](#)



Allen Institute

Allen Institute

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.dm6gpr15dvzp/v3>

**Protocol Citation:** Allen Institute for Brain Science 2024. 4% PFA + 2.5% Glutaraldehyde Fixative. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.dm6gpr15dvzp/v3> Version created by **Allen Institute**

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** May 14, 2024

**Last Modified:** May 14, 2024

**Protocol Integer ID:** 99811

**Keywords:** RP0229, tissue, electrophysiology, reagent preparation, tissue after electrophysiology recording, pfa, electrophysiology recording,

## Abstract

This protocol is used to prepare 4% Paraformaldehyde (PFA) + 2.5% Glutaraldehyde Fixative. This is utilized to fix the tissue after electrophysiology recordings.

**Note:** Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

## Attachments



RP0229\_4\_PFA\_2\_5\_Glu.



63KB

