

May 28, 2015

Adaptor Ligation Mixture (E7370)

DOI

dx.doi.org/10.17504/protocols.io.cg6tzd



New England Biolabs (NEB)

Tech. support phone: +1(800)632-7799 email: info@neb.com



New England Biolabs

New England Biolabs

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.cg6tzd>

External link: <https://www.neb.com/protocols/2012/11/15/adaptor-ligation-e7370>

Protocol Citation: New England Biolabs: Adaptor Ligation Mixture (E7370). [protocols.io](https://www.protocols.io)
<https://dx.doi.org/10.17504/protocols.io.cg6tzd>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

Created: September 29, 2014

Last Modified: January 17, 2018

Protocol Integer ID: 254

Keywords: adaptor ligation, adaptor, e7370, mixture

Troubleshooting



1 Blunt/TA Ligase Master Mix 15 μ l

2 NEBNext Adaptor for Illumina 2.5 μ l

Note

The NEBNext adaptor is provided in NEBNext Singleplex (**NEB #E7350**) or Multiplex (**NEB #E7335, #E7500**) Oligos for Illumina.

Note

If DNA input is < 100 ng, dilute the NEBNext Adaptor for Illumina 1:10 in sterile water. Use immediately to a final concentration of 1.5 μ M.

3 Ligation Enhancer 1 μ l