

Mar 31, 2016 Version 3

# Determining Genome Targeting Efficiency using T7 Endonuclease I (M0302) V.3



DOI

[dx.doi.org/10.17504/protocols.io.etbbein](https://dx.doi.org/10.17504/protocols.io.etbbein)

New England Biolabs (NEB)

Tech. support phone: +1(800)632-7799 email: [info@neb.com](mailto:info@neb.com)



**New England Biolabs**

New England Biolabs

OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.etbbein](https://dx.doi.org/10.17504/protocols.io.etbbein)

**External link:** <https://www.neb.com/protocols/2014/08/11/determining-genome-targeting-efficiency-using-t7-endonuclease-i>

**Protocol Citation:** New England Biolabs: Determining Genome Targeting Efficiency using T7 Endonuclease I (M0302). [protocols.io https://dx.doi.org/10.17504/protocols.io.etbbein](https://dx.doi.org/10.17504/protocols.io.etbbein)

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**Created:** March 31, 2016

**Last Modified:** March 27, 2021

**Protocol Integer ID:** 2627

**Keywords:** genomic targeting efficiency, genomic target, T7 Endonuclease for Gene targeting, determining genome targeting efficiency by digesting annealed PCR with T7 endonuclease, digest with T7 endonuclease I, editing efficiency, t7e1 assay, t7 endo assay

## Abstract

**T7 Endonuclease I** recognizes and cleaves non-perfectly matched DNA. This protocol describes how to determine genome targeting efficiency by digesting annealed PCR products with T7 Endonuclease I. In the first step PCR products are produced from the genomic DNA of cells whose genomes were targeted using Cas9, TALEN, ZFN etc. In the second step, the PCR products are annealed and digested with T7 Endonuclease I. Fragments are analyzed to determine the efficiency of genome targeting.

## Guidelines

### Overview:

**T7 Endonuclease I** recognizes and cleaves non-perfectly matched DNA. This protocol describes how to determine genome targeting efficiency by digesting annealed PCR products with T7 Endonuclease I. In the first step PCR products are produced from the genomic DNA of cells whose genomes were targeted using Cas9, TALEN, ZFN etc. In the second step, the PCR products are annealed and digested with T7 Endonuclease I. Fragments are analyzed to determine the efficiency of genome targeting.

### Required Materials:

- **Q5® Hot Start High-Fidelity 2X Master Mix (M0494S)**
- T7 Endonuclease I (M0302S)
- 0.25 mM EDTA
- Purified genomic DNA from targeted cells
- PCR primers to amplify a ~1kb region containing the target site
- The target site should be offset from the center of the amplicon so that digestion produces easily resolvable DNA fragments
- PCR primer design is critical. Please visit NEB's [Tools and Resources](#) page to optimize your primer design using the [NEB Tm Calculator](#)
- A PCR thermocycler with programmable temperature ramp rate
- DNA purification system - we recommend Ampure XP beads
- Apparatus to quantitate DNA - spectrophotometer or fluorometer
- Apparatus to analyze DNA fragments – e.g. Agilent Bioanalyzer, Qiagen Qiaxcel, or standard agarose gel electrophoresis

## PCR

### Cycling Conditions

Step	Temperature	Time
Initial Denaturation	98°C	30 seconds
35 cycles	98°C	5 seconds
*50-72°C	10 seconds	
72°C	20 seconds	
Final Extension	72°C	2 minutes
Hold	4-10°C	

**\*Use of the [NEB Tm Calculator](#) is highly recommended.**

**Note: Q5 Hot Start High-Fidelity 2X Master Mix does not require a separate activation step. Standard Q5 cycling conditions are recommended.**

### T7 Endonuclease I digestion:

**Anneal the PCR products in a thermocycler using the following conditions:**

**Hybridization Conditions**

Step	Temperature	Ramp Rate	Time
Initial Denaturation	95°C		5 minutes
Annealing	95-85°C	-2°C/second	
85-25°C	-0.1°C/second		
Hold	4°C		Hold

**REFERENCES:**

Guschin, D.Y., et. al.(2010) [A rapid and general assay for monitoring endogenous gene modification](#). Methods Mol Biol, 649, 247–256.

**Materials**

MATERIALS

☒ Q5 Hot Start High-Fidelity 2X Master Mix - 100 rxns **New England Biolabs Catalog #M0494S**

☒ T7 Endonuclease I - 250 units **New England Biolabs Catalog #M0302S**

## PCR

- 1 Set up a **50 µl** PCR reaction using ~100 ng of genomic DNA as a template. For each amplicon set up 3 PCR reactions (positive, negative, no-template control)

Component	50 µl reaction	Final Concentration
Q5® Hot Start High-Fidelity 2X Master Mix (M0494S)	25 µl	1X
10 µM Forward Primer	2.5 µl	0.5 µM
10 µM Reverse Primer	2.5 µl	0.5 µM
Template DNA	variable	100 ng total
Nuclease-free water	To 50 µl	

### Note

For each amplicon set up 3 PCR reactions using the following templates:

- gDNA from targeted cells (e.g. Cas9, or TALEN transfected cells)
- gDNA from negative control cells (e.g. non-specific DNA transfected cells)
- water (i.e. no template control)

### Protocol



NAME

**Q5 PCR Mixture for M0302**

CREATED BY

New England Biolabs

**PREVIEW**

- 1.1 Q5® Hot Start High-Fidelity 2X Master Mix (M0494S)


25 µL

Q5 Hot Start High-Fidelity 2X Master Mix - 100 rxns **New England Biolabs** Catalog #M0494S

- 1.2 10 µM Forward Primer, **2.5 µl**

3 µL

1.3 10  $\mu$ M Reverse Primer, 2.5  $\mu$ l

 3  $\mu$ L

1.4 Template DNA

1.5 Nuclease-free water to **50  $\mu$ l**

2 Gently mix the reaction.

3 Collect all liquid to the bottom of the tube by a quick spin if necessary.

4 Transfer PCR tubes to a PCR machine and begin thermocycling.

Note

Hot Start High-Fidelity 2X Master Mix does not require a separate activation step. Standard Q5 cycling conditions are recommended.

5 Analyze a small amount of the of the PCR product to verify size and appropriate amplification.

6 Purify the PCR reaction using 90  $\mu$ l of Ampure XP beads following the manufacturer's recommendations.

Note

Other PCR purification systems (e.g. Qiagen MinElute PCR Purification Kit PCR purification columns, or Zymo DNA Clean and Concentrator™) are acceptable.

7 Elute PCR products in **30  $\mu$ l** of water, recovering **25  $\mu$ l**.

8 Measure the concentration of the purified PCR products.

## T7 Endonuclease I digestion



9 Assemble reactions as follows:

### Protocol



NAME

**Digestion Mixture for M0302 (v2)**

CREATED BY

New England Biolabs

**PREVIEW**

9.1 DNA, **200 ng**

9.2 10X NEBuffer 2, **2 µl**

2 µL

9.3 Nuclease-free Water to **19 µl**

10 Anneal the PCR products in a thermocycler (see [guidelines](#) for conditions)

11 Add **1 µl** of the T7 Endonuclease I to the annealed PCR products

1 µL

12 Incubate at 37°C for 15 minutes

00:15:00

13 Stop the reaction by adding **1.5 µl** of 0.25 M EDTA.

2 µL

14 Purify the reaction using 36 µl of Ampure XP beads according to the manufacturer's suggestion.

### Note

This step is optional since 1 µl of the reaction will not interfere with analysis on an Agilent Bioanalyzer using DNA1000 reagents.



15 Elute the DNA fragments in **20  $\mu$ l** of water, recovering **15  $\mu$ l**.

## Analysis

16 Analyze the fragmented PCR products and determine the percent of nuclease-specific cleavage products (fraction cleaved)

17 Calculate the estimated gene modification using the following formula:  
% gene modification =  $100 \times (1 - (1 - \text{fraction cleaved})^{1/2})$