



Jan 21, 2015

Ethanol precipitation of nucleic acids (Eppendorf tubes)

DOI

<https://dx.doi.org/10.17504/protocols.io.cdrs55>



Openwetware

OpenWetWare

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.cdrs55>

External link: http://openwetware.org/wiki/Ethanol_precipitation_of_nucleic_acids

Protocol Citation: Openwetware: Ethanol precipitation of nucleic acids (Eppendorf tubes). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.cdrs55>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: August 12, 2014

Last Modified: November 09, 2017

Protocol Integer ID: 145

Keywords: ethanol precipitation of nucleic acid, nucleic acid precipitation, ethanol precipitation, nucleic acid, fact that nucleic acid, soluble in alcohol, solubility, water dipole, eppendorf tube, alcohol

Abstract

Nucleic acid precipitation is used to concentrate and/or purify nucleic acids. The below protocol is based on the fact that nucleic acids are less soluble in alcohol than in more polar water. Addition of salt further decreases solubility by competing for water dipoles; as does low temperature. Please see the [OpenWetWare](#) website for more details.

Citation: OpenWetWare contributors, 'Ethanol precipitation of nucleic acids', *OpenWetWare*, , 1 July 2012, 10:48 UTC, <http://openwetware.org/index.php?title=Ethanol_precipitation_of_nucleic_acids&oldid=611420> [accessed 21 January 2015]

Materials

STEP MATERIALS

 ammonium acetate P212121

 ammonium acetate P212121

Protocol materials

 ammonium acetate P212121



- 1 Add 1/10 volume of 3M sodium acetate, pH 5.2 or 1/2 volume of 5M ammonium acetate.

ammonium acetate P212121

- 2 Add 2-3 volumes of 100% Ethanol.

- 3 Mix and freeze overnight in -20.

Note

In general, the time you need to incubate in the freezer depends on how much nucleic acid you have, how big it is and the volume it is in. My general protocol is to freeze for 20 min to 1 hr at -80 °C. This seems to work well for most things, but you may want to freeze longer if you have only a small concentration of nucleic acid or if it is small in size(<15 nucleotides). (Kathleen)

Note

If you are in a hurry, you can also dip you epi shortly into liquid nitrogen. If you added enough ethanol, the mix won't freeze. Careful with isopropanol - it freezes more quickly. This works well for me and saves me a lengthy incubation in the fridge. (Jasu)

- 4 Spin at full speed in a standard microcentrifuge at 4 degrees for 30 minutes.

00:30:00

- 5 Decant (or carefully pipet off) the supernatant.

- 6 Dry the pellet.

Note

For this you can air dry (tubes open, ~15 min) or dry in a speedvac. DNA and RNA (if you don't have RNases in your sample) are typically hearty enough for you to air dry at 37 °C, if desired.

Note

Overdrying can make DNA hard to re-dissolve. Especially for longer DNA, I avoid vacuum drying and airdry only briefly before re-dissolving. (Jasu)



- 7 Add your desired quantity of water. Vortex and spin down to resuspend.

Note

Beware of using water unless you are sure of what you are getting in to. The "pH" of water can vary widely (I've seen from pH 5 to pH 8.5), and depurination of DNA at low pH or degradation of RNA at high pH are possibilities. Water also typically contains trace metals, which can accelerate these reactions. I typically recommend resuspension in TE (10 mM Tris-HCl, pH 7.5, 1 mM EDTA). This makes sure your nucleic acid is at a neutral pH and the EDTA will chelate any trace metals. Since they are in such small amounts, neither the buffer nor the EDTA will affect most downstream reactions. (Kathleen)