

Apr 18, 2016

Extracting DNA from bananas

DOI

dx.doi.org/10.17504/protocols.io.esvbee6



Melissa Wilson Sayres¹

¹protocols.io



Melissa mws Wilson Sayres

Arizona State University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.esvbee6>

Protocol Citation: Melissa Wilson Sayres 2016. Extracting DNA from bananas. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.esvbee6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: March 29, 2016

Last Modified: March 16, 2018

Protocol Integer ID: 2613

Keywords: dna from banana, extracting dna, banana, spanish version of this protocol, dna

Abstract


For more information and for the Spanish version of this protocol, please see [here](#).

The protocol is based on: <http://www.scientificamerican.com/article/find-the-dna-in-a-banana-bring-science-home/>.

Guidelines

Introduction

All living things, bananas and people included, pass on information from one generation to the next using the same basic material, DNA. Within an organism, each cell contains a complete set of DNA “blueprints”. DNA contains information for how to build each cell type and for many of the features that make that organism unique. These instructions are in segments of DNA called genes. Genes and the parts of our DNA that turn genes on and off hold information for how our body develops and functions. Variants of genes, called alleles, are responsible for differences in hair color, eye color, and earlobe shape.

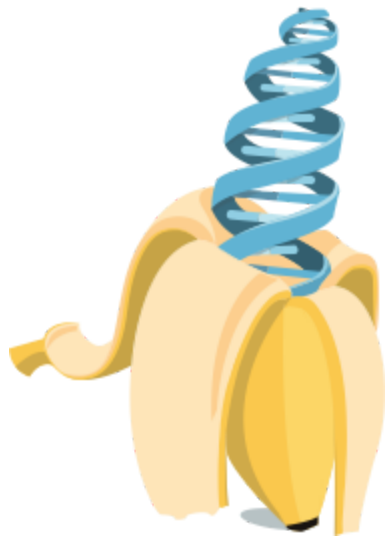



We will use common household products to break apart the cells in a banana and extract out the DNA so we can see it!

Materials

- ½ peeled ripe banana
- ½ cup hot water
- 1 tsp salt
- Resealable zip-top bag
- ½ tsp dishwashing soap
- very cold rubbing alcohol (place in freezer ahead of time)
- coffee filter
- narrow glass
- wooden stirrer

Troubleshooting



- 1 Mush the banana in the resealable bag for about a minute until it has a pudding-like consistency and all lumps are gone.
 00:01:00
- 2 Fill a cup with the hot water and salt.
- 3 Pour the saltwater mix into the bag, and close the bag. Gently mix and slosh the saltwater and mashed banana together for 30 to 45 seconds.
 00:00:30
- 4 Add dishwashing soap into the bag and gently mix the contents. You do not want the mixture to become too foamy.
- 5 Place the coffee filter in a clear glass cup, securing the top of the filter around the lip of the cup.
- 6 Pour the contents of the bag into the filter and let it sit until all of the liquid has dripped down into the cup.

Note


It is especially important to be patient here. If you try to rush this step you could break the coffee filter and mix up the banana mash with the liquid with the DNA strands in it.

- 7 Remove and discard the used coffee filter.
- 8 Tilt the glass and **slowly** add cold alcohol down the side of the cup until there is a layer that is 2.5-5cm (1-2in) thick. Keep the alcohol and liquefied banana as separate as possible.

Note

We cannot emphasize enough that this step should be done slowly. You want the cold alcohol to sit on top of the layer of liquid with the DNA in it.

- 9 Let this two-layered mixture sit for eight minutes. The longer you wait, the more defined this layer becomes. This is the DNA pieces clumping together.

 00:08:00



- 10 Stick the wooden stirrer into the cup. Spin it in place so that cloudy layer spools around it. Remove the stirrer. The substance that you see on the stirrer is DNA!