



MojoSort™ Human CD45 Nanobeads Protocol A V.1



In 1 collection

DOI

[dx.doi.org/10.17504/protocols.io.e2zb6f6](https://doi.org/10.17504/protocols.io.e2zb6f6)



Kelsey Miller¹

¹BioLegend

BioLegend

Tech. support email: tech@biolegend.com



Sam Li

BioLegend



DOI: [dx.doi.org/10.17504/protocols.io.e2zb6f6](https://doi.org/10.17504/protocols.io.e2zb6f6)

External link:

http://www.biolegend.com/media_assets/support_protocol/MojoSort_Human_CD45_Nanobeads_Protocol_12182015.pdf

Protocol Citation: Kelsey Miller . MojoSort™ Human CD45 Nanobeads Protocol A. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.e2zb6f6>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: June 03, 2016

Last Modified: August 29, 2019

Protocol Integer ID: 2873

Keywords: percentage of cd45, cd45, targeted cell, nanobead, mojosort, functional assay, cell, magnetic particle



Abstract

Note: If the percentage of CD45+ cells in your sample is less than 50%, please follow Protocol A. If it is higher than 50% then please follow protocol B.


The cells targeted by the Nanobeads are either selected or depleted by incubating your sample with the directly conjugated magnetic particles. The magnetically labeled fraction is retained by the use of a magnetic separator. After collection of the targeted cells, downstream applications include functional assays, gene expression, phenotypic characterization, etc.

Guidelines

Protocol: This procedure is optimized for the isolation of 10^7 to 2×10^8 cells per tube. If working with fewer than 10^7 cells, keep volumes as indicated for 10^7 cells. For best results, optimize the conditions to your specific cell number and tissue. Prepare fresh MojoSort™ Buffer solution by diluting the 5X concentrate with sterile distilled water.

Reagents and instruments required:

- MojoSort™ Buffer (5X) (Cat. No. 480017)
- MojoSort™ Magnet (Cat. No. 480019) or compatible magnetic separation system
- Adjustable pipettes
- 70 µm filters (one per sample)
- 5 mL (12 × 75 mm) polystyrene tubes
- Reagents for sample preparation
- Reagents and instruments (Flow cytometer) to determine yield and purity

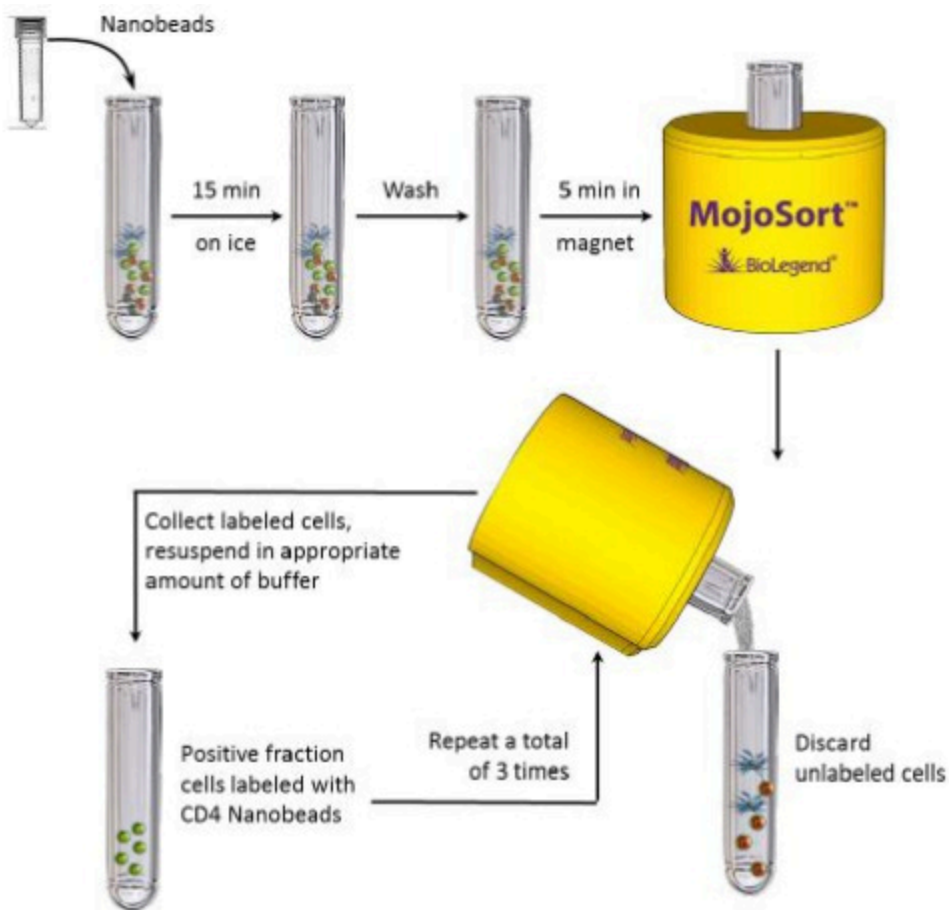





- 1 Prepare cells from your tissue of interest without lysing erythrocytes.
- 2 In the final wash of your sample preparation, resuspend the cells in MojoSort™ Buffer by adding up to 4 mL in a 5 mL (12 × 75 mm) polystyrene tube.

Note: Keep MojoSort™ Buffer on ice throughout the procedure.

- 3 Filter the cells with a 70 µm cell strainer, centrifuge at 300 x g for 5 minutes, and resuspend in an appropriate volume of MojoSort™ Buffer. Count and adjust the cell concentration to 1×10^8 cells/mL.


 00:05:00

- 4 Aliquot 100 µL of cell suspension (10^7 cells) into a new tube.
- 5 Resuspend the beads by vortexing, maximum speed, 5 touches. Add 10 µL of Nanobeads, mix well and incubate on ice for 15 minutes.
Scale up the volume accordingly if separating more cells. For example, add 100 µL for 1×10^8 cells. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.

 00:15:00

- 6 Add MojoSort™ Buffer up to 4 mL and centrifuge the cells at 300 x g for 5 minutes.
- 7 Resuspend the cells in 3 mL of MojoSort™ Buffer.
Optional: Take an aliquot before placing the tube in the magnet to monitor purity and yield

- 8 Place the tube in the magnet for 5 minutes

 00:05:00

- 9 Pour out the liquid. Resuspend the labeled cells in appropriate amount of buffer.

- 10 Repeat steps 7 – 9 on the labeled fraction 2 more times, for a total of 3 magnetic separations.

Optional: Take a small aliquot to monitor purity and yield. If desired, pool the unlabeled fractions and process simultaneously with the positive labeled cells when assessing purity and yield.