



MojoSort™ Isolation Kits Regular Protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.e2vbge6>



Kelsey Miller¹

¹BioLegend

BioLegend

Tech. support email: tech@biolegend.com



Sam Li

BioLegend

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account



DOI: <https://dx.doi.org/10.17504/protocols.io.e2vbge6>

External link:

http://www.biolegend.com/media_assets/support_protocol/MojoSort_Isolation_Kits_Regular_Protocol_V02_10192015.pdf

Protocol Citation: Kelsey Miller . MojoSort™ Isolation Kits Regular Protocol. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.e2vbge6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Created: June 03, 2016

Last Modified: August 29, 2019

Protocol Integer ID: 2869

Keywords: sample with the biotin antibody cocktail, sample with the biotin, biotin antibody cocktail, biotin, incubation with magnetic streptavidin nanobead, untouched cell, target cell, antibody cocktail, magnetic streptavidin nanobead, mojosort, functional assay, streptavidin nanobead, antibody, cell, cells of interest, isolation, magnetic separator, nanobead

Abstract

Target cells are depleted by incubating your sample with the biotin antibody cocktail followed by incubation with magnetic Streptavidin Nanobeads (Cat. No. 480015/480016). The magnetically labeled fraction is retained by the use of a magnetic separator. The untouched cells are collected. These are your cells of interest; do not discard the liquid. Some of the downstream applications include functional assays, gene expression, phenotypic characterization, etc.

Guidelines


Reagents and instruments required:

- MojoSort™ Buffer (5X) (Cat. No. 480017)
- MojoSort™ Magnet (Cat. No. 480019) or compatible magnetic separation system
- Adjustable pipettes
- 70 µm filters (one per sample)
- 5 mL (12 × 75 mm) polystyrene tubes
- Reagents for sample preparation
- Reagents and instruments (Flow cytometer) to determine yield and purity

Protocol:

This procedure is optimized for the isolation of 107 to 2×10^8 cells per tube. If working with fewer than 107 cells, keep volumes as indicated for 107 cells. For best results, optimize the conditions to your specific cell number and tissue. Prepare fresh MojoSort™ Buffer solution by diluting the 5X concentrate with sterile distilled water.

Application notes: To use this product in magnetic separation columns, a titration of the cocktail/beads should be performed. Optimal concentration for magnetic separation columns is lot-specific.

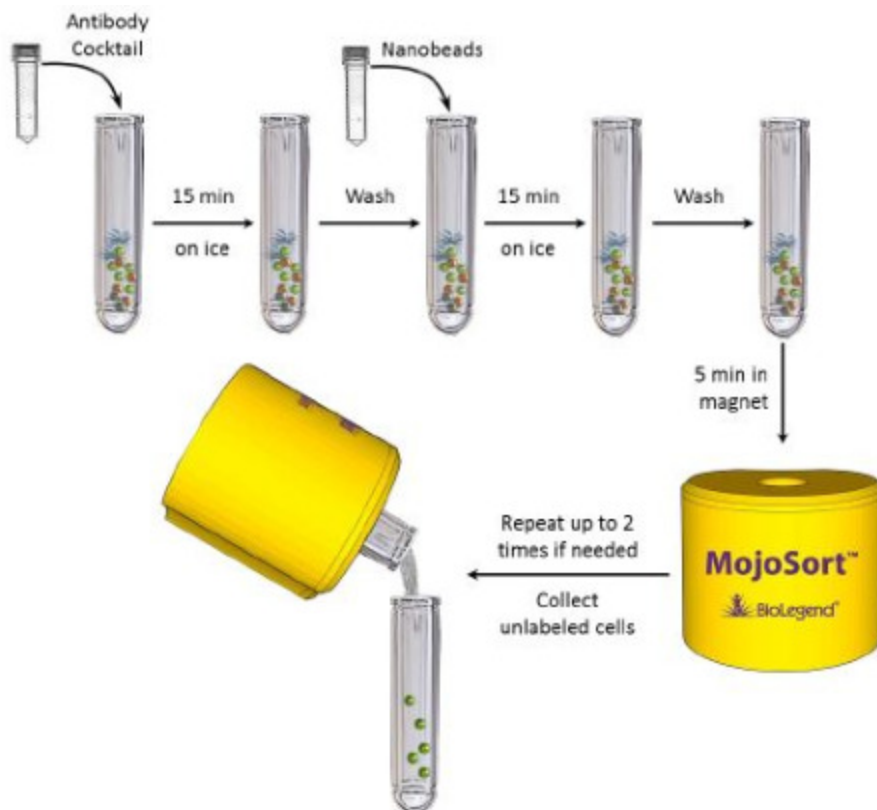



- 1 Prepare cells from your tissue of interest without lysing erythrocytes.
- 2 In the final wash of your sample preparation, resuspend the cells in MojoSort™ Buffer by adding up to 4 mL in a 5 mL (12 × 75 mm) polystyrene tube.

Note: Keep MojoSort™ Buffer on ice throughout the procedure.

- 3 Filter the cells with a 70 µm cell strainer, centrifuge at 300 x g for 5 minutes, and resuspend in an appropriate volume of MojoSort™ Buffer. Count and adjust the cell concentration to 1×10^8 cells/mL.

 00:05:00

- 4 Aliquot 100 µL of cell suspension (10^7 cells) into a new tube. Add 10 µL of the Biotin-Antibody Cocktail, mix well and incubate on ice for 15 minutes. Scale up the volume accordingly if separating more cells. For example, add 100 µL for 1×10^8 cells. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.

Optional: Keep unused cells, or take an aliquot before adding the cocktail to monitor purity and yield.

 00:15:00

- 5 Wash the cells by adding MojoSort™ Buffer up to 4 mL; centrifuge the cells at 300 x g for 5 minutes.

 00:05:00

- 6 Discard supernatant and resuspend in 100 µL of MojoSort™ Buffer.

- 7 Resuspend the beads by vortexing, maximum speed, 5 touches. Add 10 µL of Streptavidin Nanobeads. Mix well and incubate on ice for 15 minutes. Scale up the volume accordingly if separating more cells. For example, add 100 µL for 1×10^8 cells. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.

 00:15:00

- 8 Wash the cells by adding 3 mL of MojoSort™ Buffer; centrifuge at 300 x g for 5 minutes, discard supernatant.

Optional: Take an aliquot before placing the tube in the magnet to monitor purity and yield.

 00:05:00

- 9 Resuspend the cells in 3 mL of MojoSort™ Buffer.
Note: If you observe aggregates, filter the suspension. To maximize yield, you can



disrupt the aggregates by pipetting the solution up and down.

- 10 Place the tube in the magnet for 5 minutes.

 00:05:00

- 11 Pour out and collect the liquid. These are your cells of interest; **DO NOT DISCARD.**
- 12 If needed, add 3 mL of MojoSort™ Buffer and repeat steps 10 and 11 with the magnetically labeled fraction up to two times, and then pool the unlabeled fractions.

Note: Repeating the magnetic separation increases the yield, without a strong impact on the purity. The yield will typically increase about 8 – 10% with a second separation, and about 2 – 5% with a third separation. The purity may decrease 1 – 2% with each separation. Optional: Take a small aliquot before placing the tube in the magnet to monitor purity and yield.