

Aug 16, 2016

Version 2

## Mucin biotinylation - NHS-biotin reagent V.2

DOI

[dx.doi.org/10.17504/protocols.io.ex4bfqw](https://dx.doi.org/10.17504/protocols.io.ex4bfqw)

Victoria Linderberg

Thomas Crouzier Lab



Victoria Linderberg

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ex4bfqw>

**Protocol Citation:** Victoria Linderberg 2016. Mucin biotinylation - NHS-biotin reagent. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.ex4bfqw>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**Created:** May 08, 2016

**Last Modified:** March 10, 2018

**Protocol Integer ID:** 2780

**Keywords:** biotin to bovine serum mucin, mucin biotinylation, biotin reagent this protocol, biotin reagent, biotin to amine, bovine serum mucin, biotin, mucin protein, mucin, nhs molecule, protein, nh

## Abstract

This protocol describes how to attach biotin to Bovine Serum Mucin (BSM). The procedure uses a botin-NHS molecule to attach biotin to amines in the mucin proteins.

## Attachments



MAN0011206\_EZ\_NHS\_

Bi...

86KB



J. Biol. Chem.-1998-...

263KB

## Troubleshooting

## Safety warnings

- ⚠ NHS-Biotin reagents are moisture sensitive, so if stored cold, they need to be fully equilibrated to room temperature before opening the vial! Also, the biotin reagent should be prepared right before usage, since it quickly hydrolyses and becomes inactive, i.e. do not prepare stock solutions of the biotin reagent and store them!

- 1 Calculate how much NHS-Biotin that is needed for biotinylation of the desired amount of mucin (assumed to have  $M_w$  500 000 Da) and the molar ratio of biotin:mucin should be 1:2000.
- 2 Prepare Phosphate Buffered Saline (PBS) with pH 7.4.
- 3 Dissolve mucin in PBS to a concentration of 2mg/ml.
- 4 Take out the NHS-Biotin reagent and equilibrate it in room temperature before opening the vial.
- 5 Dissolve NHS-Biotin in dimethylsulfoxide (DMSO) to a concentration of 10 mM.
- 6 Add the pre-calculated volume of biotin reagent to the protein solution and incubate at 4 °C for 4 hours.
- 7 To remove unbound biotin, exchange the buffer for water by overnight dialysis.
- 8 Transfer the solution to another tube, freeze and freeze-dry.