

Feb 11, 2016

PBS / DAPI 1:2000 Dilution Mixture

DOI

<https://dx.doi.org/10.17504/protocols.io.eifbcbn>

Alan Cone¹

¹Wright State University

Ju Lab

Yeast Protocols, Tools, a...



Alan Cone

Wright State University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eifbcbn>

Protocol Citation: Alan Cone 2016. PBS / DAPI 1:2000 Dilution Mixture. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.eifbcbn>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: February 11, 2016

Last Modified: March 28, 2018

Protocol Integer ID: 2343

Keywords: dilution mixture, pb, dapi

Materials

STEP MATERIALS

 1X PBS (Phosphate-buffered saline)

 DAPI (2.5mg/mL)

 1X PBS (Phosphate-buffered saline)

 DAPI (2.5mg/mL)

Protocol materials

 DAPI (2.5mg/mL)

 1X PBS (Phosphate-buffered saline)



1 Add 1 mL 1 x PBS to a 1.5 mL microcentrifuge tube.

 1 μ L

 1X PBS (Phosphate-buffered saline)

2 Add 0.5 μ L of a 2.5 mg/mL (or a 1:2000 dilution) of DAPI to the 1 x PBS.

 DAPI (2.5mg/mL)