

Feb 04, 2016

Submixture for Step1 of Labyrinthula growth medium

DOI

<https://dx.doi.org/10.17504/protocols.io.ehmbb46>

Janina Brakel

Protist Research to Opti...

Collier Lab



Janina Brakel

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ehmbb46>

Protocol Citation: Janina Brakel 2016. Submixture for Step1 of Labyrinthula growth medium. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ehmbb46>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: February 03, 2016

Last Modified: November 12, 2017

Protocol Integer ID: 2317

Keywords: labyrinthula growth medium, submixture for step1, submixture



- 1 Add 0.6 g glucose
- 2 Add 0.06 g yeast extract
- 3 Add 0.06 g of peptone
- 4 Add 7.2 g of agar
- 5 Add 600 ml distilled water.