

May 10, 2024

AAV Craniotomy

 [Science Advances](#)

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.5jyl829q7l2w/v1>

Santiago Unda¹, Roberta Marongiu¹, Mihaela Stavarache¹, Michael G. Kaplitt¹

¹Weill Cornell Medical College



Eileen Ruth Torres

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5jyl829q7l2w/v1>

External link: <https://doi.org/10.1126/sciadv.adn6016>

Protocol Citation: Santiago Unda, Roberta Marongiu, Mihaela Stavarache, Michael G. Kaplitt 2024. AAV Craniotomy. protocols.io <https://dx.doi.org/10.17504/protocols.io.5jyl829q7l2w/v1>

**Manuscript citation:**

Kokotos AC, Antoniazzi AM, Unda SR, Ko MS, Park D, Eliezer D, Kaplitt MG, Camilli PD, Ryan TA Phosphoglycerate kinase is a central leverage point in Parkinson's disease-driven neuronal metabolic deficits. Science Advances 10(34). doi: [10.1126/sciadv.adn6016](https://doi.org/10.1126/sciadv.adn6016)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 08, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 99452

Keywords: ASAPCRN, surgery, mouse, AAV, aav craniotomy this protocol, aav craniotomy, injection of aav, aav, mouse brain, injection, stereotactic frame, hamilton syringe, brain

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: 020608

Abstract

This protocol is for injection of AAV into a mouse brain using a stereotactic frame and Hamilton syringe.

Guidelines

Note that sterile technique must be followed during this procedure to avoid health complications.

Anesthesia

- 1 Weigh mice.
- 2 Administer intraperitoneal mixture of ketamine (ButlerAnimal Health Supply) and xylazine (Lloyd Laboratories) at concentrations of 110 and 10 mg/kg of body weight, respectively.
- 3 Confirm induction of anesthesia.

Injection

- 4 Place mouse into stereotactic frame (David Kopf Instruments) and place eye lubricant.
- 5 Remove fur from skull area using either electric shaver or depilatory cream.
- 6 Sterilize skull area and then cut open with scalpel.
- 7 Adjust brain so that it is level. Locate bregma and zero X and Y axes on stereotaxic frame.
- 8 Move to injection coordinates and drill small hole into the skull.
- 9 Zero the Z axis on the stereotaxic frame.
- 10 Prepare AAV solution by diluting AAV in PBS with 500mM Mg2/Cl to a final dilution of total of 2×10^{11} genomic particles/mL.
- 11 Using a 10 μ L stereotactic syringe with a 33 G needle attached to a micro-infusion pump (World Precision Instruments), withdraw the AAV solution.
- 12 With the loaded syringe, navigate the needle to the desired X, Y, Z coordinates.



- 13 Infuse the brain with AAV at a flow rate of 0.1 - 0.4 $\mu\text{L}/\text{min}$.
- 14 To prevent reflux, after each infusion, leave the injection needle in place for 5 min, withdraw a short distance (0.3 - 0.5 mm), and then leave in the new position for an additional 2 min before complete removal of the needle.
- 15 Close skin with non-absorbable sutures.

AAV Preparation

- 16 Inject a total of 2×10^{11} genomic particles/mL of vector is injected into the appropriate coordinates.

Post-procedure

- 17 Give meloxicam 2mg/kg subcutaneously.
- 18 Perform health checks every day for the first 72 hrs post-injection.
- 19 Remove sutures 15 days post-injection.