

Oct 18, 2019

Version 2

Ablation of β -cells in Tg(ins:NTR-mCherry) transgenic zebrafish using metronidazole V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.8fkhtkw>



Sandra Rieger¹

¹University of Miami

Diabetic Complications Consortium

Tech. support email: rmcindoe@augusta.edu



Lili Liang

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8fkhtkw>

External link: <https://www.diacomp.org/shared/document.aspx?id=221&docType=Protocol>

Protocol Citation: Sandra Rieger 2019. Ablation of β -cells in Tg(*ins:NTR-mCherry*) transgenic zebrafish using metronidazole. protocols.io <https://dx.doi.org/10.17504/protocols.io.8fkhtkw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 18, 2019

Last Modified: October 18, 2019

Protocol Integer ID: 28876

Keywords: hyperglycemia, Neuropathy, transgenic zebrafish strain, hyperglycemia in larval zebrafish, zebrafish strain, nfsb gene, larval zebrafish, bacterial enzyme nitroreductase, red fluorescent protein mcherry, cell, hyperglycemia, transgenic, addition of the prodrug metronidazole, cell ablation, cell ablation system, diabetic complication, prodrug metronidazole

Abstract

Summary:

To induce hyperglycemia in larval zebrafish, utilize a genetic β -cell ablation system involving the Tg(*ins:NTR-mCherry*) transgenic zebrafish strain (courtesy of Dr. Michael Parsons, Johns Hopkins University). This strain expresses the bacterial enzyme Nitroreductase (NTR), which is encoded by the *nfsB* gene, in β -cells of the pancreas starting around 24 hours post fertilization (hpf). Because NTR is genetically fused to the red fluorescent protein mCherry, β -cells expressing NTR can be visualized with fluorescence microscopy. Upon addition of the prodrug metronidazole (MTZ) to the embryo medium, NTR converts MTZ into a toxin, which selectively ablates NTR-expressing β -cells (Curado et al., Nature Protocols 3, 948-954 (2008)). This can be monitored by the disappearance of the fluorescence under a confocal microscope.

Diabetic Complication:



Neuropathy



Materials

MATERIALS

⊗ Sea salt **Instant Ocean (Pet store)**

⊗ 1-phenyl-2-thiourea (PTU) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P7629**

⊗ Metronidazole **Merck MilliporeSigma (Sigma-Aldrich) Catalog #M3761**

⊗ Dimethyl sulfoxide (DMSO) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #276855**

⊗ Tg (ins:NTR-mCherry)

Reagent/Material	Required concentration
Tg(<i>ins:NTR-mCherry</i>)	
Sea salt	0.3 g/ L
1-phenyl-2-thiourea (PTU)	0.003 %
Metronidazole	10 mM
Dimethyl sulfoxide (DMSO)	0.5 %

Reagent Preparation:

Reagent 1: Embryo medium

Preparation: Add 0.03 % Instant Ocean salt (Pet store) into double distilled water.

Reagent 2: Embryo medium with 1-phenyl-2-thiourea (PTU) to prevent pigment formation.

Preparation: To embryo medium in a small beaker add a final of 0.003% PTU and dissolve overnight using a magnetic stir bar. Keep PTU solution in the dark.

Reagent 3: Metronidazole for β -cell ablation.

Preparation: First make a 2x MTZ stock (20 mM) by adding 0.17 g MTZ to a 50 ml Falcon tube. Next, add 49.5 ml of embryo medium containing 0.003 % PTU and 0.5 ml DMSO (makes 1 %). Dissolve MTZ by vortexing for ~ 5 minutes or until dissolved. Dilute the 2x MTZ stock 1:1 with embryo medium to make a 10 mM MTZ/ 0.5 % DMSO working solution. Store 50 % of 2x MTZ stock in the fridge overnight. Use the other half immediately. Discard the 2x MTZ stock latest after 2 days due to degradation of the MTZ.

Reagent 4: DMSO control solution.

Preparation: Make a 0.5 % DMSO solution by adding 0.25 ml DMSO to 50 ml of embryo medium.

Note:

Sigma-Aldrich, RRID:SCR_008988

Protocol:

- 1 Collect fertilized Tg (*ins:NTR-mCherry*) transgenic zebrafish embryos and incubate at 28°C for 24 hours. Remove dead embryos and replace old embryo medium with fresh embryo medium containing 0.003 % PTU to prevent pigment formation. Remove chorions from embryos according to Protocol 1 and add the 1x MTZ working solution to the embryos. Incubate in the dark (due to light-dependent degradation of MTZ) at 28°C for 24 hours. Assess ablation efficiency with confocal imaging. At 48 hpf, exchange media with fresh MTZ solution and incubate for another 24 hours. Repeat confocal analysis.

Potential Pitfalls:

- 2
 1. Ablation efficiency is low: MTZ might be old or only incompletely dissolved in the embryo medium. MTZ might alternatively be degraded if embryos were incubated in light, as MTZ is light sensitive.
 2. MTZ does not dissolve: DMSO and volume of initial solution is both crucial for dissolving the MTZ. High concentrations will not go into solution. Vortex for extended periods if it does not dissolve within 5 minutes.
 3. MTZ treatment causes larvae to appear unhealthy or die: MTZ may be old or degraded due to excessive light exposure. We found that this affects the health of larval zebrafish. DMSO concentrations, if too high (above 1 - 2 %) may also cause larvae to appear unhealthy over time (bend axis). The same is true for larvae that are incubated for more than 2 days in MTZ. If ablation is not sufficient, try to incubate for 2 days and then let larvae recover for 1 or 2 days until a second MTZ treatment.