



Feb 16, 2026

Acute Midbrain Slice Preparation

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g776dkgwz/v1>

Nicola Berretta¹, Ezia Guatteo^{1,2}, Nicola Biagio Mercuri^{3,4}

¹Experimental Neurology Laboratory, IRCCS Santa Lucia Foundation, Rome, Italy;

²Department of Medical, Human Movement and Well-Being Sciences, University of Naples 'Parthenope', Naples, Italy;

³Fondazione Santa Lucia IRCSS; ⁴Department of Systems Medicine, University of Rome Tor Vergata, Rome, Italy



Cristian González-Cabrera

Leibniz Institute für Neurologie

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g776dkgwz/v1>

Protocol Citation: Nicola Berretta, Ezia Guatteo, Nicola Biagio Mercuri 2026. Acute Midbrain Slice Preparation . **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.q26g776dkgwz/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 16, 2026

Last Modified: February 16, 2026

Protocol Integer ID: 243379

Keywords: acute midbrain slice preparation, μm horizontal midbrain slice, using chilled nmdg, chilled nmdg, protective cutting solution, graded sodium reintroduction, based protective cutting solution, sodium reintroduction, preparation, electrophysiological recording, storage in hepe

Abstract

This protocol describes preparation of 250 μm horizontal midbrain slices containing the VTA using chilled NMDG-based protective cutting solution, followed by graded sodium reintroduction and storage in HEPES-based aCSF prior to electrophysiological recording.

Materials

****NMDG-Based Cutting Solution (mM)****

92 NMDG, 2.5 KCl, 1.25 NaH_2PO_4 , 30 NaHCO_3 , 20 HEPES, 25 glucose, 2 thiourea, 5 Na-ascorbate, 3 Na-pyruvate, 0.5 CaCl_2 , 10 MgSO_4 . Equilibrate with 95% O_2 - 5% CO_2 , pH 7.3. Keep chilled.

****HEPES-Based Storage Solution (mM)****

92 NaCl, 2.5 KCl, 1.25 NaH_2PO_4 , 30 NaHCO_3 , 20 HEPES, 25 glucose, 2 thiourea, 5 Na-ascorbate, 3 Na-pyruvate, 2 CaCl_2 , 2 MgSO_4 . Equilibrate with 95% O_2 - 5% CO_2 , pH 7.3. Store at room temperature.

Step-by-Step Procedure

- 1 Anesthetize the mouse with halothane and decapitate rapidly.
- 2 Remove the brain and immediately place it into chilled carbogenated NMDG solution.
- 3 Mount the brain for horizontal slicing.
- 4 Cut 250 μ m horizontal slices containing the midbrain using a vibratome.
- 5 Transfer slices to NMDG solution at 33.0 C and incubate for 15 min.
- 6 Gradually reintroduce sodium by adding increasing volumes of 2 M NaCl solution every 5 min for 40 min total.
- 7 Transfer slices to HEPES-based aCSF at room temperature.
- 8 Allow at least 1 h recovery before recording.