

Oct 05, 2025

Add Arabinose

DOI

<https://dx.doi.org/10.17504/protocols.io.dm6gpm558gzp/v1>

igemnthuth¹

¹iGEM_NTHU_Taiwan

iGEM NTHU



igemnthuth

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.dm6gpm558gzp/v1>

Protocol Citation: igemnthuth 2025. Add Arabinose. protocols.io
<https://dx.doi.org/10.17504/protocols.io.dm6gpm558gzp/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: October 03, 2025

Last Modified: October 05, 2025

Protocol Integer ID: 228985

Keywords: synthetic replication operon, expression of the synthetic replication operon, sgrna repression system, orthogonal replication system, orthogonal replication, improving orthogonal replicon transformation efficiency, orthogonal replicon transformation efficiency, downstream operon component expression, arabinose, induced dcas9

Abstract

To control expression of the synthetic replication operon in the orthogonal replication system, this step uses an **arabinose-induced dCas9–sgRNA repression system** to reduce downstream operon component expression, thereby **improving orthogonal replicon transformation efficiency**.

Materials

- LB broth
- 10% glycerol solution (pre-chilled)
- Arabinose
- ddH₂O
- Pre-chilled electroporation cuvettes (2 mm gap)
- Pre-chilled centrifuge tubes (50 mL × 2; 1.5 mL × 1)
- Shaking incubator
- Ice-water bath and ice bucket
- Spectrophotometer (for OD600 measurement)
- Refrigerated centrifuge
- Electroporator



Pre-culture

- 1 Pick colonies that have successfully taken up the following plasmids:
 - **EcOrep operon plasmid**
 - **dCas9 + sgRNA expression plasmid** (e.g., pRT17/19/21)
 - **Mutant DNAP plasmid**
- 2 Inoculate into **15 mL LB broth** supplemented with the **corresponding antibiotics for all three plasmids**, and incubate **overnight at 37 °C, 220 rpm** (~12 h).

Main Culture and Arabinose Induction

- 3 The next day, transfer **5 mL** of the overnight culture into **200 mL** fresh LB broth (**1:40 dilution**) in a **2 L shake flask**; incubate at **37 °C, 220 rpm**.
- 4 When **OD600 ≈ 0.2**, add **arabinose** to a **final concentration of 10 mM** (e.g., add **2 mL of 1 M stock** to **200 mL** culture) and continue incubation.
- 5 Grow to **OD600 ≈ 0.4**, then immediately proceed to the following steps for electroporation pre-treatment.

Preparation of Competent Cells

- 6 Place the culture on ice to cool for **10 minutes** (gentle swirling optional).
- 7 Centrifuge at **4 °C, 4500 × g for 3 minutes**; carefully discard the supernatant.
- 8 Wash the cell pellet with **50 mL pre-chilled 10% glycerol**; repeat the wash **twice** (total **2 washes**).
- 9 Finally, **resuspend** the pellet in **1 mL pre-chilled 10% glycerol** to obtain the competent cells.
- 10 **Aliquot** into several **100 µL** tubes (for subsequent **O-replicon electroporation**). This completes the **arabinose treatment** stage.