



Oct 11, 2022

ADN PURIFICATION FROM AGAROSE GELS WITH GENELUTE GEL EXTRACTION KIT

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygx9p5dg8j/v1>

Ana Belem García González¹, Georgina Elisa Diego Macías¹, Irán Alessandra Chaparro Rodríguez¹,
Jair Alexis Gardea Sáenz¹

¹Tecnologico de Monterrey Campus Chihuahua



Georgina Elisa Diego Macías

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx9p5dg8j/v1>

Protocol Citation: Ana Belem García González, Georgina Elisa Diego Macías, Irán Alessandra Chaparro Rodríguez, Jair Alexis Gardea Sáenz 2022. ADN PURIFICATION FROM AGAROSE GELS WITH GENELUTE GEL EXTRACTION KIT . **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kxygx9p5dg8j/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 11, 2022

Last Modified: October 11, 2022

Protocol Integer ID: 71132

Keywords: dna purification from agarose gel, adn purification from agarose, adn purification, dna purification, genelute gel extraction kit, protocol for dna purification, agarose gel, purification, agarose, dna, extraction

Abstract

Protocol for DNA purification from DNA purification from agarose gels with GeneElute Gel Extraction Kit

Materials

GenElute Gel Extraction Kit (NA1111, Sigma-Aldrich):

- Column Preparation Solution
- Gel Solubilization Solution
- Wash Solution Concentrate G
- Elution Solution
- GenElute Binding Column G
- Collection Tubes, 2 mL

Isopropanol

Ethanol

Sodium Acetate 3M, pH 5.2

Thermoblock at 50-60°C.

Electrophoresis equipment for agarose gel

Tubes of 1.5 mL



- 1 Previous electrophoresis : Realization of an electrophoresis in a agarose gel of the product of ADN, charging the maximum amount of DNA per well (if necessary, use more than one well)
- 2 Extraction of the band of interest: Extract the band of interest of the gel, removing the excess of agarose
- 3 Weighting the band: Weight the piece of gel removed in a tarated tube.
- 4 Gel Solubilization: Add 3 volumes of Gel Solubilization Solution to the piece of gel, meaning, for each  100 mg of agarose gel, add  300 μL of Gel Solubilization Solution. Incubate at 50-60°C for  00:10:00 minutes or until the piece of gel has completely dissolved. Give Vortex every 2 or 3 minutes in this incubation time. 10m
- 5 Preparation of the column (This can be done alongside step 4). Add GenElute Binding Column G in a recolection tube of 2 mL. Add  500 μL of the Column Preparation Solution to each Binding Column. Centrifugate for  00:01:00 minute at 12,000 to 16,000 x g. Discard liquid 1m
- 6 Check the color of the sample: Once the gel had dissolved completely, make sure that the color is yellow before continuing. If the color is red, add  10 μL of the Sodium Acetate 3M, pH 5.2 buffer and mix. Add until the color is yellow.
- 7 Add isopropanol: Add a volume of the gel of isopropanol at 100% and mix until homogenized.
- 8 Bind ADN: Charge the sample of step 7 to the binding column prepared in step 5. Is normal a change of color from yellow to red once the sample is charged. If the volume of the sample is bigger than  700 μL , charge the sample in proportions of  700 μL . Centrifugate for one minute at 12,000-16,000 x g every time after the sample is charged. Discard the liquid obtained from the centrifugation.
- 9 Wash the column: Add  700 μL of Wash Solution to the binding column. Centrifugate for one minute at 12,000-16,000 x g. Remove the binding column of the tube and discard liquid. Colocate again the binding column in the recolection tube and centrifuge one more time for one minute to eliminate the excess of ethanol



10

Elute the DNA: Transfer the binding column to a clean, esteril tube. Add  50 μ L of the Elution Solution to the center of the membrane and incubate for  00:01:00 minute. Centrifuge for  00:01:00 minute at 12,000-16,000 x g. In case of purifying plasmid DNA or linear fragmented of DNA with a weight exceeding 3Kb, preheat the Elute Solution at 65 °C

2m