

Jun 19, 2018

Version 2

Adult Mouse Spleen Dissociation (On ice) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.q5cdy2w>

Andrew Potter¹

¹CCHMC

Human Cell Atlas Metho...



Andrew Potter

CCHMC

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q5cdy2w>

Protocol Citation: Andrew Potter 2018. Adult Mouse Spleen Dissociation (On ice). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.q5cdy2w>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: June 19, 2018

Last Modified: June 19, 2018

Protocol Integer ID: 13188

Keywords: spleen, mouse spleen into single cell, adult mouse spleen dissociation, mouse spleen, single cell, cell, mg tissue, cell size, tissue, ice, variety of cell size, dissociation

Abstract

Protocol used to dissociate adult (8-10 wk) mouse spleen into single cells. Attained >95% viability, a variety of cell sizes, and ~10 million cells from 12 mg tissue.

Guidelines

Collagenase Enzyme Mix (two tubes, 1 mL each)

7.5 mg/mL Collagenase A (Sigma, 10103578001)

7.5 mg/mL Collagenase Type 4 (Worthington, CLS-4)

100 µg/mL soybean trypsin inhibitor (Sigma, 10109886001)

125 U DNase (Applichem, A3778)

5 mM CaCl₂

740 µL DPBS (no Ca, Mg)

+12 mg chopped spleen / tube

Materials

MATERIALS

 Red Blood Cell Lysis Buffer Hybri-Max Merck MilliporeSigma (Sigma-Aldrich) Catalog #R7757

Before start

-Set centrifuges to 4° C.

-Make two tubes of 1 mL enzyme mix.

-Make ~25 mL of DPBS/0.04% BSA



- 1 Isolate whole spleen and immerse in ice-cold PBS.
- 2 Using sterile forceps, transfer spleen to petri dish on ice. Chop whole spleen coarsely for 45 sec using razor blade on petri dish, on ice until a fine paste.
 00:00:45 mince on ice
- 3 Weigh out 12 mg of minced spleen on petri dish. Using razor blade, transfer to 1.5 mL tube containing 1 mL of enzyme mix on ice.
- 4 Incubate tube on ice for 10 minutes. Triturate 10X every 2 mins and shake every min.
 00:10:00 Incubate on ice
 00:02:00 triturate 10x every 2 min
 00:01:00 Shake every min
- 5 After 10 mins of digestion, let tissue chunks settle for 1 min on ice & remove 80% of supernatant with released cells & filter using 70 μ M filter on 50 mL conical, on ice. Rinse filter with 5 mL ice-cold PBS/0.04% BSA. Leave filter and 50 mL conical on ice, it will be used for the steps as well.
 5 μ L ice-cold PBS/BSA 0.04%
- 6 Add additional 1 mL enzyme mix to tissue chunks.
- 7 Continue to triturate 10x every 2 minutes and shake every minute while incubating on ice, for 10 additional minutes (21 min total time).
 00:10:00 Incubate on ice
 00:02:00 triturate 10x every 2 min
 00:01:00 shake every min
- 8 After 21 min total time, triturate 10x and add entire volume of cell digestion to 70 μ M filter on 50 mL conical, using the same filter and tube as in the previous step. Rinse filter w/5 mL ice-cold PBS/0.04% BSA.
 5 μ L ice-cold PBS/BSA 0.04%
- 9 Transfer flow-through to 15 mL conical. Spin 650 g for five minutes at 4 °C. After spin, remove supernatant, leaving 100 μ L volume in 15 mL conical tube.
 4 °C



00:05:00 spin 650 g for 5 min

- 10 Perform RBC lysis: add 1 mL RBC lysis buffer to the 100 μ L of cells and triturate 10X. Let sit 3 minutes on ice. After incubating 3 min in RBC lysis buffer, add 10 mL ice-cold PBS/BSA 0.04% in the 15 mL conical to dilute the RBC lysis buffer. Pipet mix.

00:03:00 incubate on ice

10 μ L ice-cold PBS/BSA 0.04%

1 μ L RBC lysis buffer

- 11 Spin 650 g for 5 mins at 4 °C. Remove supernatant and re-suspend in 1 mL ice-cold PBS/BSA 0.04%.

4 °C

00:05:00 spin at 650 g for 5 min

- 12 Examine cells using hemocytometer with trypan blue. Adjust concentration to 1000 cells / μ L for 10x chromium or 100 cells / μ L for DropSeq.