



Jul 18, 2024

Agrobacterium-Mediated Transient Expression in *Nicotiana benthamiana* Leaves



DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzj7prlx1/v1>

Maryam Rahmati Ishka¹

¹Boyce Thompson Institute (BTI)

Julkowska Lab



Maryam Rahmati Ishka

Boyce Thompson Institute (BTI)

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzj7prlx1/v1>

External link: <http://n/a>

Protocol Citation: Maryam Rahmati Ishka 2024. Agrobacterium-Mediated Transient Expression in *Nicotiana benthamiana* Leaves. protocols.io <https://dx.doi.org/10.17504/protocols.io.rm7vzj7prlx1/v1>

Manuscript citation:

n/a

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 17, 2024

Last Modified: July 18, 2024

Protocol Integer ID: 103602

Keywords: Agrobacterium infiltration, Nicotiana benthamiana, plant construct delivery, transient protein expression, mediated transient expression in nicotiana benthamiana leaf, agrobacterium infiltration of tobacco leaf, nicotiana benthamiana leaf, tobacco leaf, agrobacterium, agrobacterium infiltration, mediated transient expression, labeled protein, protein

Disclaimer

n/a

Abstract

This procedure outlines the Agrobacterium infiltration of tobacco leaves for the transient expression of fluorescently labeled proteins.

Image Attribution

n/a

Guidelines

n/a

Materials

Infiltration medium: 10 mM MgCl₂, 10 mM MES (pH 5.6). Autoclave for 20 min. Add acetosyringone immediately before using it  MgCl₂ Applied Biosystems (ThermoFisher Scientific)

 20X MES Buffer Thermo Fisher Scientific Catalog #NP0002

 Acetosyringone Merck MilliporeSigma (Sigma-Aldrich) Catalog #2478-38-8

 1g Rifampicin USP Grade G-Biosciences Catalog #RC-191 to a final concentration of 200 µM.

Safety warnings

 n/a

Ethics statement

n/a

Before start

n/a

Transform constructs into *Agrobacterium tumefaciens* strains that are compatible with vectors. Strain GV3101 is typically used (GV2260 can also be used). Make glycerol stock of.

- 1 Refresh bacteria containing the construct of interest from glycerol stock on agar YEP (or any other bacterial media such as LB) medium with an appropriate antibiotic. Incubate at 28°C for two days. If *A. tumefaciens* GV3101 is used add **rifampicin**.
- 2 Inoculate **5 mL** of liquid YEP medium with appropriate **antibiotic** in a test tube. Grow at 28°C for **24 hours** with shaking at 250–300 rpm. Be sure to grow bacterial strain with **p19** as a silencing inhibitor as well!
- 3 Measure the OD₆₀₀ using 100 µL of bacterial culture.
- 4 Calculate the culture volume needed for an OD₆₀₀ of 0.2 – 0.5 per construct for a 4 mL total volume of Infiltration medium. 4 mL are used for test spot infiltration. The specific OD₆₀₀ to use depends on how well expressed your construct is in *N. benthamiana*, the general range is 0.1 – 0.5. Use an OD₆₀₀ of 0.3 for P19.
(I use 0.3 for all constructs and it works well).
- 5 Pipette the appropriate volume of each strain together in 2 mL tubes.
- 6 Centrifuge combined bacteria at 8000×g for 2 min at room temperature. Carefully remove all supernatant.
- 7 Add 200 mM **acetosyringone** to the **Infiltration medium** at a 1:1000 ratio.
Note: acetosyringone is light-sensitive, keep out of light until ready for use.
- 8 Add 2 mL of infiltration medium with **acetosyringone** to each pellet, and resuspend. Leave in the dark at room temperature for 1–2 hrs (or even longer but not overnight).
- 9 Infiltrate 4-week-old *N. benthamiana* (pre-flowering) leaf using 1 mL syringe.
- 10 Grow plants for 2–3 days after infiltration (up to 5 days, depending on expression level of constructs).



- 11 Punch **1 cm** leaf discs from leaves and observe under the microscope to visualize the fluorescence.

Protocol references

n/a