



Jun 23, 2022

Amicon NMWCO Filter Concentration and Buffer Exchange

DOI

dx.doi.org/10.17504/protocols.io.8epv59594g1b/v1

Lauren Adams¹

¹Northwestern University

Kelleher Research Group

Tech. support email: kelleher-ofc@northwestern.edu



Kelleher KRG Research Group

Northwestern University, National Resource for Translational...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8epv59594g1b/v1>

Protocol Citation: Lauren Adams 2022. Amicon NMWCO Filter Concentration and Buffer Exchange. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.8epv59594g1b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: June 23, 2022

Last Modified: June 23, 2022

Protocol Integer ID: 65205

Keywords: buffer exchange this amicon nmwco filter protocol, amicon nmwco filter concentration, amicon nmwco filter protocol, buffer exchange, assists with buffer exchange, concentration dilute elution fraction, concentration, filter, containing protein, protocol, protein

Abstract

This Amicon NMWCO filter protocol helps concentration dilute elution fractions containing protein as well as assists with buffer exchange if necessary.

Materials

Amicon NMWCO Filter (MWCO dependent on the protein target)

Centrifuge

0.5% TFA or 0.2% FA in mass spec grade H₂O

LoBind microcentrifuge tubes

Troubleshooting



- 1 Wash the filter three times with Optima grade water and centrifuging at 14,000 x g. Equilibrate the column by washing once with the same buffer as the elution buffer containing the target protein for ~10-30 minutes or until the volume has decreased.
- 2 Load the elution fraction containing the target protein into the filter and centrifuge at 14,000 x g for ~10-30 minutes or until the volume has decreased. Repeat this until the total elution volume has decreased $\leq 25 \mu\text{L}$ for LC-MS or $\leq 80 \mu\text{L}$ for I2MS. If the buffer is already acidic (TFA or FA), no additional buffer exchange is required. However, if the elution buffer is not acidic (ex: 100 mM Tris Base), buffer exchange should be performed.
- 3 Gradually incorporate larger volumes of 0.2-0.5% TFA or FA to the sample volume after sequential centrifuge cycles until the majority of the final buffer is TFA or FA. Remove sample from filter by centrifuging upside down at 14,000 x g into a clean LoBind tube. Bring the final volume up to 25 μL for LC-MS or 80 μL for I2MS with 0.2-0.5% TFA or FA before analyzing by mass spectrometry.