



Jun 24, 2023

Analysis of ER structures in Cultured Induced Neuron axons

 [Nature Cell Biology](#)

 In 3 collections

DOI

dx.doi.org/10.17504/protocols.io.yxmvm37j9l3p/v1

Melissa Hoyer¹, Harper JW¹

¹Harvard Medical School

Harper

ASAP



Melissa Hoyer

Harvard Medical School

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yxmvm37j9l3p/v1>

External link: <https://doi.org/10.1038/s41556-024-01356-4>

Protocol Citation: Melissa Hoyer, Harper JW 2023. Analysis of ER structures in Cultured Induced Neuron axons. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.yxmvm37j9l3p/v1>

**Manuscript citation:**

Melissa J. Hoyer, Cristina Capitanio, Ian R. Smith, Julia C. Paoli, Anna Bieber, Yizhi Jiang, Joao A. Paulo, Miguel A. Gonzalez-Lozano, Wolfgang Baumeister, Florian Wilfling, Brenda A. Schulman, J. Wade Harper (2024) Combinatorial selective ER-phagy remodels the ER during neurogenesis. *Nature Cell Biology* doi: [10.1038/s41556-024-01356-4](https://doi.org/10.1038/s41556-024-01356-4)

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 23, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 83960

Keywords: ASAPCRN, ER-phagy, iNeurons, axonal er accumulation within synaptic er bouton, er structures in axon, axonal er accumulation, synaptic er bouton, endoplasmic reticulum, er remodeling by autophagy, er to degradative autophagy machinery, extensive er remodeling during differentiation, deficient neuron, er structure, deficient neurons in vivodisplay, neuron, embedded receptor, phagy receptor fam134b, cultured induced neuron, degradative autophagy machinery, axon, er proteome, underlying er remodeling, autophagy, analysis of er structure, vast proteomic landscape, tractable induced neuron, phagy receptor, monitoring extensive er remodeling, receptor, selective er, linking er, proteome, protein, er remodeling

Funders Acknowledgements:

ASAP

Grant ID: ASAP-000282

Abstract

The endoplasmic reticulum (ER) has a vast proteomic landscape to preform many diverse functions including protein and lipid synthesis, calcium ion flux, and inter-organelle communication. The ER proteome is remodeled in part through membrane-embedded receptors linking ER to degradative autophagy machinery (selective ER-phagy)^{1,2}. A refined tubular ER network^{3,4} is formed in neurons within highly polarized dendrites and axons^{5,6}. Autophagy-deficient neurons *in vivodisplay* axonal ER accumulation within synaptic ER boutons,⁷ and the ER-phagy receptor FAM134B has been genetically linked with human sensory and autonomic neuropathy^{8,9}. However, mechanisms and receptor selectivity underlying ER remodeling by autophagy in neurons is limited. Here, we combine a genetically tractable induced neuron (iNeuron) system for monitoring extensive ER remodeling during differentiation. With this system, we imaged fixed iNeuron cultures, imaged these via confocal fluorescence microscopy, and quantified ER structures in axons.

Materials

| A | B | C |
|---|--|--|
| REAGENT or RESOURCE | SOURCE | IDENTIFIER |
| Cell line | | |
| NGN2 inducible embryonic stem cells | See protocol dx.doi.org/10.17504/protocols.io.br9em93e | CVCL_9773 (modified from this source line) |
| Chemicals | | |
| Dulbecco's MEM (DMEM), F12 | | |
| Phosphate Buffered Saline 1X | Corning | 21-031-CV |
| 16% Paraformaldehyde, Electron-Microscopy Grade | Electron Microscopy Science | 15710 |
| Triton-X | Sigma | T8787 |
| DAPI | Thermo Fisher Scientific | D1306 |
| E8 components | See protocol dx.doi.org/10.17504/protocols.io.bsacnaaw | |
| ND1 and ND2 components | See protocol dx.doi.org/10.17504/protocols.io.br9em93e | |
| Software | | |
| Nikon Imaging Software Elements | 5.21.3 (Build 1489) | https://cellprofiler.org/ SCR_014329 |
| Cell Profiler | CellProfiler v4.0.6 | https://cellprofiler.org/ SCR_007358 |
| Fiji | ImageJ V.2.0.0 | https://imagej.net/software/fiji/ SCR_002285 |
| Hardware | | |
| Thermo Neon™ Transfection System | Thermo Fisher Scientific | MPK5000 |



| | A | B | C |
|--|--|---|---|
| | Yokogawa CSU-X1 spinning disk confocal on a Nikon Ti-E inverted microscope | | |

Troubleshooting

Genetically modify Ngn2-inducible embryonic stem (ES) cell H9 line using Cas9

- 1 1. Genetic editing of Ngn2-inducible ES cells is done using the following protocol
"Electroporation of Cas9 protein into human pluripotent stem cells"
([dx.doi.org/10.17504/protocols.io.br87m9zn](https://doi.org/10.17504/protocols.io.br87m9zn))

Differentiation of Stable Cell ES H9 line to induced neurons (iN)

- 2 1. Differentiation to induced neurons (iN) is done by following the protocol "Neural differentiation of AAVS1-TREG3-NGN2 pluripotent stem cells"
([dx.doi.org/10.17504/protocols.io.br9em93e](https://doi.org/10.17504/protocols.io.br9em93e))

Analysis of endoplasmic reticulum (ER) in iN via confocal microscopy

- 3 Fixation and immunofluorescence is next done to label the ER, axons and dendrites of the iNs
- 3.1 1)To help get spread out iNeurons for imaging, when cells get 90 percent confluent at any point in the day 5-7 range, cells are plated into onto 35 mm-glass bottom dishes (No. 1.5, 14 mm glass diameter, MatTek) at a low confluency (approximately 1×10^5 cells per 2mL dish).
- 3.2 2)At day 20, iN are fixed. phosphate buffered saline (1xPBS) with 8% paraformaldehyde is prewarmed to 37C and then slowly added to an equal volume of media on the cells for a resulting 4% paraformaldehyde solution. iN are incubated with this solution at 37C for 15min.
- 3.3 3)Fixed iN are washed one time with room temperature 1xPBS
- 3.4 4)Fixed iN are solubilized with 0.1% triton-X in PBS for 5 min at RT.
- 3.5 5)Fixed iN are then blocked for 30 min at RT with sterile filtered blocking buffer (1% bovine serum albumin, 0.1% triton-X in PBS)
- 3.6 6)Primary antibodies for Endoplasmic reticulum (ER), axons and dendrites are added to blocking buffer at 1:300 and then spun down for 1min at 10,000 x g. The blocking buffer is completely removed and then the antibody in blocking buffer is applied to the cells (100mL applied to the center of the well where the glass coverslip is attached) for 1h at RT.

- 3.7 7) Cells are washed 4 times with PBS (5min for each wash).
- 3.8 8) Fluorescently conjugated secondary antibodies are added to blocking buffer at 1:300 and then spun down for 1min at 10,000 x g. The blocking buffer is completely removed and then the antibody in blocking buffer is applied to the cells (100mL applied to the center of the well where the glass coverslip is attached) for 1h at RT.
- 3.9 9) Cells are washed 4 times with PBS (5min for each wash) and left in 1XPBS. Dapi at 1:10,000 is added in the second 1XPBS wash.
- 4 iN are imaged via confocal fluorescence microscopy and ER, axons and dendrites of the iNs are measured.
- 4.1 1) Cells are imaged at RT using a Yokogawa CSU-X1 spinning disk confocal on a Nikon Ti-E inverted microscope at the Nikon Imaging Center in Harvard Medical School. Nikon Perfect Focus System was used to maintain cell focus over time. The microscope is equipped with a Nikon Plan Apo 40x/1.30 N.A or 100x/1.40 N.A objective lens. 488nm (100mW) and 561nm (100mW) laser lines are controlled by AOTF. All images are collected with a Hamamatsu ORCA-ER cooled CCD camera (6.45 μm^2 photodiode) with Nikon Imaging Elements.
- 4.2 2) Z series are displayed as maximum z-projections and saved using Fiji software.
- 4.3 3) The ER positive structures are detected using Cell Profiler with the same pipeline applied for each condition (see attached CellProfiler pipeline for specific thresholding details). Cell nuclei are detected using a "identify primary objects" module with 50-to-125-pixel units min and max and a global minimum cross-entropy thresholding method. Both axon and dendrite area are identified using a "identify primary objects" module with 10-to-40-pixel units min and max using default thresholding method. Each ER structure is identified using a "identify primary objects" module with 5-to-50-pixel units min and max and a global robust background thresholding method. The number of ER structures masked specifically within the axonal regions is related to the number of nuclei in each field of view to yield the number of ER structures in axons per nuclei readout. Object size and shape is measured, and these values are also exported.
- 4.4 4) Each channel z series are brightness and contrast adjusted equally and then converted to RGB for publication using FIJI software.

Protocol references



WTvATG12vpentaregionsinaxonswi...