

Dec 29, 2023

## Antibody Purification and Labeling

DOI

[dx.doi.org/10.17504/protocols.io.kxygx34rdg8j/v1](https://dx.doi.org/10.17504/protocols.io.kxygx34rdg8j/v1)

Maya Brewer<sup>1</sup>, Yuantee Zhu<sup>1</sup>, Jeff Spraggins<sup>2</sup>, Mark De Caestecker<sup>3</sup>

<sup>1</sup>Vanderbilt University Medical Center; <sup>2</sup>Vanderbilt University;

<sup>3</sup>Division of Nephrology, Vanderbilt University Medical Center

VU Biomolecular Multim...

Human BioMolecular Atl...



Maya Brewer

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx34rdg8j/v1>

**Protocol Citation:** Maya Brewer, Yuantee Zhu, Jeff Spraggins, Mark De Caestecker 2023. Antibody Purification and Labeling. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kxygx34rdg8j/v1>

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** December 26, 2023

**Last Modified:** December 29, 2023

**Protocol Integer ID:** 92732

**Keywords:** HuBMAP, BIOMIC, Antibody Purification, Antibody Labeling, Immunofluorescence, Antibody Validation, antibody purification, antibody labeling, process for antibody purification, validating antibody labeling, subsequent labeling for direct immunofluorescence, antibody, direct immunofluorescence, purification, immunofluorescence, labeling this protocol, labeling, subsequent labeling, protocol

## Abstract

This protocol describes the process for antibody purification, subsequent labeling for direct immunofluorescence, and validating antibody labeling.

## Materials

### Reagents:

Protein G High Performance SpinTrap Protein G High Performance SpinTrap (Millipore Sigma, GE28-9031-34)

AB Buffer Kit (Millipore Sigma, GE28-9030-59)

MilliporeSigma Amicon Pro Purification System, MWCO 100KDa (Fisher Scientific, ACS510012)

100 ug of Antibody

10X Phosphate Buffered Saline (Thermo Fisher Scientific, #28374)

Double Distilled Water to dilute PBS to 1X

DyLight Microscale Antibody Labeling Kit

- DyLight 488 Microscale Antibody Labeling Kit (ThermoFisher, 53025)
- DyLight 594 Microscale Antibody Labeling Kit (ThermoFisher, 53045)
- DyLight 650 Microscale Antibody Labeling Kit (ThermoFisher, 84536)
- DyLight 749/755 Microscale Antibody Labeling Kit (ThermoFisher, 84539)
- DyLight 800 Microscale Antibody Labeling Kit (ThermoFisher, 53063)

### Equipment:

- Centrifuge (swinging-bucket and fixed-angle rotors)
- Thermo Scientific Nanodrop 1000 Spectrophotometer (OR Nanodrop that reads the absorbance at 280 nm to calculate the concentration of purified proteins; contaminants/ buffers that absorb around 280 nm will affect protein concentration)
- Nanodrop must also have capability to read absorbance at 493 nm, 593 nm, 652 nm, 754 nm, and 770 nm to quantify moles of dye: moles of protein ratio.

## Troubleshooting

## Antibody Purification




- 1 This is to remove any BSA, azide, glycine that is often added by the manufacturer for stabilization. If the antibody has nothing added, skip this section.
- 2 Prepare buffers from AB buffer kit:
  - 2.1 10X binding buffer must be diluted to 1X (9 parts ddH<sub>2</sub>O and 1 part 10X binding buffer).
  - 2.2 10X elution buffer must be diluted to 1X (9 parts ddH<sub>2</sub>O and 1 part 10X elution buffer).
  - 2.3 Neutralizing buffer does not have to be diluted in any way. Prepare two collection tubes containing 30 µl of neutralizing buffer for the eluted fractions.
- 3 Resuspend medium of spintrap column
- 4 Remove the bottom cap from the column and save for later use.
- 5 Place the column in a 2 mL collection tube and centrifuge at 100 x g for 30 seconds to remove storage solution.
- 6 Add 600 µl of binding buffer to the tube
- 7 Centrifuge at 100 x g for 30 seconds
- 8 Add no more than 600 µL of the antibody to the tube
  - 8.1 Several sample applications can be made subsequently as long as the capacity of the column is not exceeded.



- 9 Secure the top and bottom cap on top the tube and incubate for 4 minutes while gently mixing
- 10 Remove the bottom cap and centrifuge the column at 100 x g for 30 seconds
- 11 Add 600 ul of binding buffer to the column to wash
- 12 centrifuge at 100 x g for 30 seconds
- 13 Add 600 ul of the binding buffer to the column for a second wash
- 14 Centrifuge at 100 x g for 30 seconds
- 15 Add 400 ul of elution buffer and mix by inverting tube several times
- 16 Place the column in one of the 2 ml collection tubes containing 30 ul of neutralizing buffer
- 17 Centrifuge at 100 x g for 30 seconds
- 18 Add 400 ul of elution buffer and mix by inverting tube several times
- 19 Place the column in the other 2 ml collection tube containing 30 ul of neutralizing buffer
- 20 Centrifuge at 100 x g for 30 seconds

## Buffer Exchange

- 21 Assemble Amicon Pro Affinity Concentrator by carefully attaching the Amicon Ultra device to the exchange device

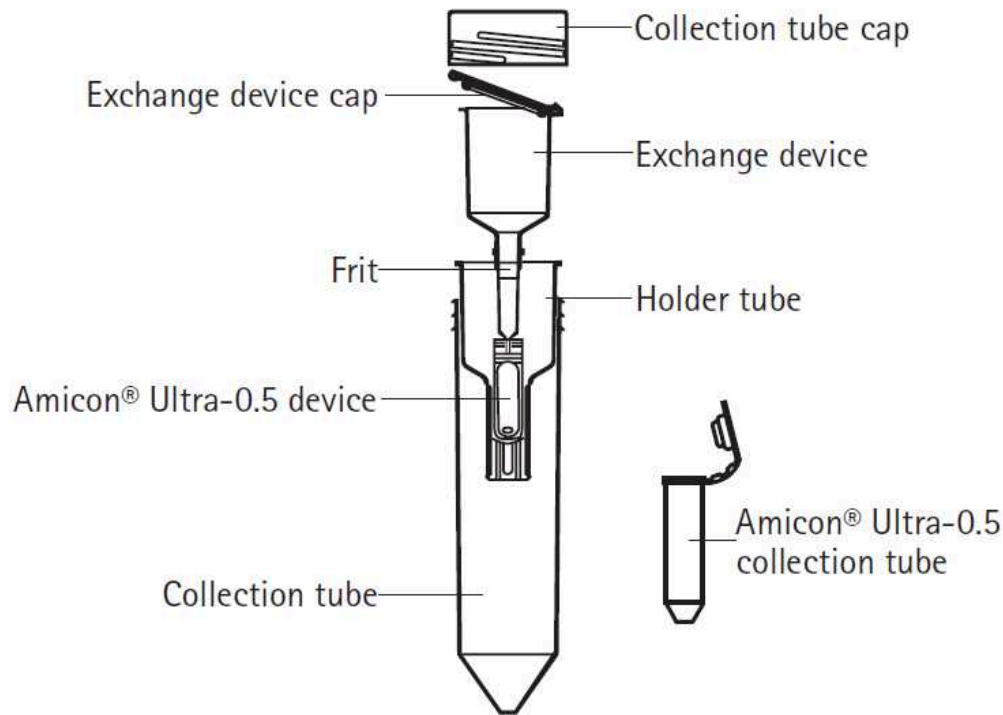

- 22 Once the concentrator is assembled, remove the collection tube cap, lift the exchange device cap, and add 1 ml of binding buffer or PBS to moisten the cellulose membrane of the Amicon Ultra device. This wash ensures that the antibody does not stick to the membrane upon its addition.
- 23 Centrifuge at 4000 x g for 3 minutes
- 24 Add 1 ml binding buffer or PBS and antibody collected in neutralizing buffer to the exchange device
- 25 Centrifuge for 15 minutes at 4000 x g in a swinging-bucket rotor
- 26 Add 2 mL of binding buffer or PBS to the device and centrifuge again at 4000 x g for 10 minutes twice
- 27 Collect purified antibody from the device by reverse spin



- 27.1 Place a collection tube on top of the Amicon Ultra-0.5 device
- 27.2 Invert the assembly and centrifuge in a fixed-angle rotor at 1000 x g for 2 minutes
- 28 Use a nanodrop to measure the purified antibody at an absorbance of 280 nm
- 28.1 Make sure that the sampling arm on the nanodrop is up
- 28.2 Using a P10 pipette, add 2  $\mu$ L of the sample onto the lower measurement pedestal
- 28.3 Lower the sampling arm and measure the absorbance of the antibody at A280. The concentration should be given in mg/mL

## Labeling Protocol

- 29 **Antibodies should be in a buffer free of any amine-containing compounds (glycine, Tris, or ammonium ions) and stabilizing proteins (bovine serum albumin). These compounds will interfere with the labeling reaction.** Labeling of antibody occurs by the DyLight dye linking to the primary amine (R-NH<sub>2</sub>) in peptides and proteins.
- 30 Follow labeling kit protocol:  
[https://www.thermofisher.com/document-connect/document-connect.html?url=https://assets.thermofisher.com/TFS-Assets%2FMSG%2Fmanuals%2FMAN0011587\\_DyLight\\_Microscale\\_Antibody\\_Label\\_UG.pdf](https://www.thermofisher.com/document-connect/document-connect.html?url=https://assets.thermofisher.com/TFS-Assets%2FMSG%2Fmanuals%2FMAN0011587_DyLight_Microscale_Antibody_Label_UG.pdf)
- 31 Collect labeled antibody from the collection tube. Aliquot then store at 4°C short term, and -20°C long term.
- 31.1 \*The conjugates can survive freeze thawing but you will need to evaluate each new antibody you use to make sure. We always evaluate one freeze/thaw cycle by repeating staining

## Validating Antibody Labeling

- 32 After labeling antibodies, we initially validate antibody labeling by staining human kidney sections with the fluorophore conjugated antibody at different dilutions, as described in the antibody validation protocol ([dx.doi.org/10.17504/protocols.io.5jyl8pwwdg2w/v1](https://doi.org/10.17504/protocols.io.5jyl8pwwdg2w/v1)). We also compare staining by indirect immunofluorescence on the same section by indirect immunofluorescence using both the conjugated and original unconjugated antibody.
- 33 If we show that the conjugated and unconjugated antibody is staining the tissue using indirect immunofluorescence but the directly conjugated fluorophore is undetectable, this suggests that the signal from the fluorophore is too low (i.e.: there may be too little dye conjugation). Conversely, if antibody staining is detectable with unconjugated antibodies, but the conjugated antibody is undetectable using both methods, it is possible too much dye is conjugated to the antibody and is inhibiting the epitope binding sites. To address this, we evaluate the dye: antibody ratio and repeat the conjugation reaction with different ratios of dye to antibody in the conjugation reaction depending on the results, as described below
- 33.1 The dye to antibody ratio measures the average number of dye molecules conjugated to each antibody molecule using serial dilutions of the conjugated antibody in PBS, and measuring the absorbance of the conjugate at A280 and the maximum excitation wavelength for the dye.
- 33.2 Ideally, the optimal dye to antibody ratio is around 2, meaning there are about 2 molecules of dye per molecule of antibody. Some dyes such as the Alexa Fluors allow higher ratios upwards to 5-8 without overlabeling the antibody. On this basis, if the dye to antibody ratio is  $>2$  and staining does not work, we repeat antibody conjugation reaction with reduced dye: antibody ratios. If the ratio of dye to antibody is  $> 2$ , then labeling is repeated after addressing the successful purification of the antibody, followed by using different ratio of dye to antibody, or decreasing the incubation time of the labeling reaction
- 33.3 Section D of this document outlines instructions for calculating the dye to antibody ratio [https://www.thermofisher.com/document-connect/document-connect.html?url=https://assets.thermofisher.com/TFS-Assets%2FMSG%2Fmanuals%2FMAN0011587\\_DyLight\\_Microscale\\_Antibody\\_Label\\_UG.pdf](https://www.thermofisher.com/document-connect/document-connect.html?url=https://assets.thermofisher.com/TFS-Assets%2FMSG%2Fmanuals%2FMAN0011587_DyLight_Microscale_Antibody_Label_UG.pdf)
- 34 If the antibody was conjugated in house, and we show that the conjugated and unconjugated antibody is staining the tissue using indirect immunofluorescence but the directly conjugated fluorophore is undetectable, this suggests that the signal from the fluorophore is too low (i.e.: there may be too little dye conjugation). Conversely, if antibody staining is detectable with unconjugated antibodies, but the conjugated antibody is undetectable using both methods, it is possible too much dye is conjugated to the antibody and is inhibiting the epitope binding sites.



- 34.1 If the antibody was conjugated commercially, and we show that the unconjugated stains the tissue while the conjugated antibody does not it is possible there is either too little or too much dye conjugated to the antibody. If this occurs consider ordering a different antibody, a different lot of the same fluorophore conjugated antibody, or a different fluorophore conjugate, and retest