

May 10, 2024

Apomorphine-induced rotations

 [Science Advances](#)

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v92yozl3e/v1>

Santiago Unda¹, Mihaela Stavarache¹, Michael G. Kaplitt¹, Roberta Marongiu¹

¹Weill Cornell Medical College



Eileen Ruth Torres

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v92yozl3e/v1>

External link: <https://doi.org/10.1126/sciadv.adn6016>

Protocol Citation: Santiago Unda, Mihaela Stavarache, Michael G. Kaplitt, Roberta Marongiu 2024. Apomorphine-induced rotations. protocols.io <https://dx.doi.org/10.17504/protocols.io.x54v92yozl3e/v1>

**Manuscript citation:**

Kokotos AC, Antoniazzi AM, Unda SR, Ko MS, Park D, Eliezer D, Kaplitt MG, Camilli PD, Ryan TA Phosphoglycerate kinase is a central leverage point in Parkinson's disease-driven neuronal metabolic deficits. Science Advances 10(34). doi: [10.1126/sciadv.adn6016](https://doi.org/10.1126/sciadv.adn6016)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 08, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 99446

Keywords: ASAPCRN, behavior, motor impairment, 6-OHDA, effects of apomorphine, apomorphine, rotational behavior, ohda lesion efficiency, induced rotation, light cycle, dark phase of light cycle, experiment, effect

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: 020608

Abstract

To test for 6-OHDA lesion efficiency, rotational behaviors are assessed while mice are under the effects of apomorphine. This test is run during the dark phase of light cycle. Experiments are performed by an examiner blinded to treatment groups.

Materials

- Apomorphine (Sigma Aldrich) dissolved in 0.9% saline/0.1% ascorbate solution
- 40cm x 40cm open field arenas
- Video camera and Ethovision or similar tracking software

Preparation

- 1 Weigh mice and habituate to the testing room for at least 1 hour prior to the start of testing.
- 2 Prepare apomorphine in 0.9% saline/0.1% ascorbate solution to dose each mouse with 0.25mg/kg via subcutaneous injection.
*Note that this solution is light-sensitive and should be shielded from light.
- 3 Clean open field arenas.
- 4 Set up tracking software on computer to record trials of  00:45:00 starting once an animal is detected within the arena. Using Ethovision, the automatic animal detection with deep learning settings is used to track center to-nose and center to-tail points in the defined arenas.

45m

Testing

- 5 Test mice in groups based on the number of open field arenas available. Before testing each group, start the tracking software. The trial time should not start until the animals are placed inside the arena.
- 6 For each group, inject each mouse with 0.25 mg/kg apomorphine via subcutaneous injection, and immediately place in the center of the corresponding open field arena. Repeat for the other mice in the same group.
- 7 Run the trial for the  00:45:00 period.
- 8 Remove mice and place back into cages.
- 9 Clean open field arenas, reset the tracking software for the next trial, and repeat steps for the rest of the groups.

45m

Analysis



- 10 Use the video recordings to analyze and count the numbers of rotations >50 degrees clockwise as well as and counterclockwise.