



Sep 22, 2024

Version 2

Arabidopsis-Microbe interaction seed treatment V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg394z1g25/v2>



Tao-Ho Chang¹

¹Plant Health Care Program, Academy of Circular Economy, National Chung Hsing University, Taichung

PSR Lab Protocols



Tao-Ho Chang

Program in Plant Health Care, Academy of Circular Economy, N...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg394z1g25/v2>

Protocol Citation: Tao-Ho Chang 2024. Arabidopsis-Microbe interaction seed treatment. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kqdg394z1g25/v2> Version created by **Tao-Ho Chang**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 05, 2023

Last Modified: September 22, 2024

Protocol Integer ID: 91876

Keywords: Arabidopsis thaliana, Seed priming, Beneficial microbe, Plant stress, Root structure architecture, microbe interaction seed treatment, treated seed, arabidopsi, bulk soil, plant, soil

Funders Acknowledgements:

National Sciece and Technology Council, Taiwan

Grant ID: 112-2313-B-005 -008 -

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

The treated seeds can grow in medium or bulk soil to determine the impact of novel materials on plants.

Guidelines

Seed treatment of arabidopsis seeds is a suitable, simple method for many plant-microbe interactions.

Materials

Arabidopsis thaliana seeds (Col-0)

50% Bleach

Distilled water

1/2 MS medium (pH = 5.7)



Safety warnings

 The only harmful chemical is 50% bleach (2% sodium hypochlorite). Please make sure wear gloves when in the seed sterile steps.

Before start

Seed priming is an important method that increases the health of the plant.



Seeds sterilisation

15m

1 The arabidopsis seeds are immersed with 50% bleach (2% Sodium hypochlorite) in 1.5 μ L tubes.

2

🌀 100 rpm, 28°C, 00:10:00

10m

3

🌀 100 rpm, 28°C, 00:00:10

10s

4 Gently remove the supernatant and leave the seeds in the tube.

5 The sterile seeds are immersed in distilled water.

6



Repeat the process of step #3 to #5 for 10 times

Seeds treatment

1d

7 The seeds were immersed in the specific concentration of treatment
1. Bacteria treatment: OD value to 0.1 and diluted 10 times.
2. Fungus treatment: The final concentration of spore suspension is 100 spores mL^{-1} .

8

🌀 100 rpm, 28°C, 01:00:00

1h

9

🌀 150 rpm, 4°C, 23:00:00

23h

Wash seeds

10s

10 The treated seeds are moved in a filter mesh (Favorgen) with a collection tube.



- 11  100 rpm, 28°C, 00:00:10 10s
- 12 Remove the filter through in collection tube
- 13 Add 500 µL of distilled water to the filter column and resuspend the seeds.
- 14  Repeat the process of step #11 to #13 for 10 times