

Mar 26, 2020 Version 1

Artificial Cerebrospinal Fluid V (ACSF.V) V.1

DOI

dx.doi.org/10.17504/protocols.io.bd8xi9xn

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Dillan Brown

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bd8xi9xn

Protocol Citation: Allen Institute for Brain Science 2020. Artificial Cerebrospinal Fluid V (ACSF.V). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bd8xi9xn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 25, 2020

Last Modified: March 26, 2020

Protocol Integer ID: 34807

Keywords: RP0205, Reagent, reagent preparation, Artificial Cerebrospinal Fluid, ACSF, surgery,

Abstract

This protocol describes how to prepare Artificial Cerebrospinal Fluid (V). Artificial Cerebrospinal Fluid (V) is used during surgical procedures to maintain tissue hydration and minimize tissue damage, adhesion formation and infection.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



RP0205_Artificial_Ce...

49KB

