



Jul 09, 2024

ATPase assay/ADP-glo kit

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl4q1x8vo5/v1>

Rania Abou El Asrar¹, Peter Vangheluwe¹

¹KU Leuven, Aligning Science Across Parkinson's (ASAP)



Rania Abou El Asrar

KU Leuven

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl4q1x8vo5/v1>

Protocol Citation: Rania Abou El Asrar, Peter Vangheluwe 2024. ATPase assay/ADP-glo kit. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.3byl4q1x8vo5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 06, 2023



Last Modified: July 09, 2024

Protocol Integer ID: 91889

Keywords: atpase assay, atpases with polyamine, atpase, activity of p5b, p5b, polyamine, assay, glo kit this protocol

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000458

Abstract

This protocol is successfully used to assess activity of P5B-ATPases with polyamines as substrate.

Materials

ADP-glo max assay (Promega, V7002)

DDM (Inalco, 1758-1350)

- 1 Prepare master mix on ice: 5 µg microsomes, 2.5 µg DDM (Inalco, 1758-1350), reaction buffer (50 mM MOPS-KOH, pH 7, 100 mM KCl, 11 mM MgCl₂ supplemented with 1 mM DTT and protease inhibitor (Sigmafast, Sigma, S8830-20TAB).
- 2 Incubate master mix for 30 minutes at room temperature in overhead shaker.
- 3 Pipet master mix (18.75 µl) into a white 96-well plate.
- 4 Prepare serial dilution of the substrate to be tested in 0.1 M MOPS-KOH, pH 7.
- 5 Pipet substrate dilution series (5 µl) into the 96-well plate.
- 6 Incubate for 5 minutes at room temperature then 5 minutes at 37 °C.
- 7 Add 5 mM pure ATP (1.25 µl) to initiate the reaction.
- 8 Incubate for 20 minutes at 37 °C followed by 10 minutes at room temperature.
- 9 Stop the reaction by addition of 25 µl ADP-glo reagent.
- 10 Incubate at room temperature for 40 minutes.
- 11 Add 50 µl/well of ADP-glo max reagent.
- 12 Incubate in the dark at room temperature for 60 minutes.



13 Measure luminescence (integration: 500 ms)