

Nov 19, 2020

Automated high throughput viral concentration from wastewater using the KingFisher Flex platform

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.bptemnje

Smruthi Karthikeyan¹, Greg Humphrey¹


¹University of California, San Diego

Coronavirus Method De...



Smruthi Karthikeyan

University of California, San Diego

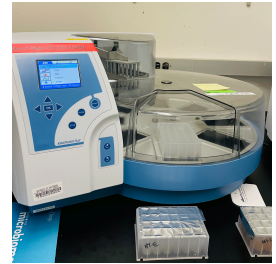

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bptemnje

External link: <https://www.medrxiv.org/content/10.1101/2020.11.16.20232900v1>

Protocol Citation: Smruthi Karthikeyan, Greg Humphrey 2020. Automated high throughput viral concentration from wastewater using the KingFisher Flex platform. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bptemnje>

Manuscript citation:

<https://www.medrxiv.org/content/10.1101/2020.11.16.20232900v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 17, 2020

Last Modified: November 19, 2020

Protocol Integer ID: 44614

Keywords: WBE, high-throughput, viral concentration, wastewater


Abstract





Large-scale wastewater surveillance has the ability to greatly augment the tracking of infection dynamics especially in communities where the prevalence rates far exceed the testing capacity. However, current methods for viral detection in wastewater are severely lacking in terms of scaling up for high throughput. In the present study, we employed an automated magnetic-bead based concentration approach for viral detection in sewage that can effectively be scaled up for processing 24 samples in a single 40-minute run. The method compared favorably to conventionally used methods for viral wastewater concentrations with a limit of detection of 8.809 viral gene copies/ml from input sample volumes as low as 10ml and can enable the processing of over 100 wastewater samples in a day.

Protocol materials

 Nanotrap Magnetic Virus Particles (10) **Ceres Nano Catalog #44202**

 MagMAX[®] Microbiome Lysis Solution **Thermo Fisher Catalog #A42361**


 MagMAX[®] Microbiome Ultra Nucleic Acid Isolation Kit, with bead plate **Thermo Fisher Catalog #A42357**

- 1 Aliquot  10 mL of raw sewage (no pre-filtering step required) into two wells of the KingFisher 24 deep-well plate (Cat# 95040470, Thermo Fisher). Each well should have  5 mL of the raw sewage sample. For instance,  5 mL of sample PL09 should be filled in A1 of plate 1 and  5 mL in A1 of plate 2.

20m



Raw, unfiltered sewage

- 2 Add  75 μ L of  Nanotrap Magnetic Virus Particles (10) Ceres Nano Catalog #44202 to each of the wells containing the raw sewage sample.

5m





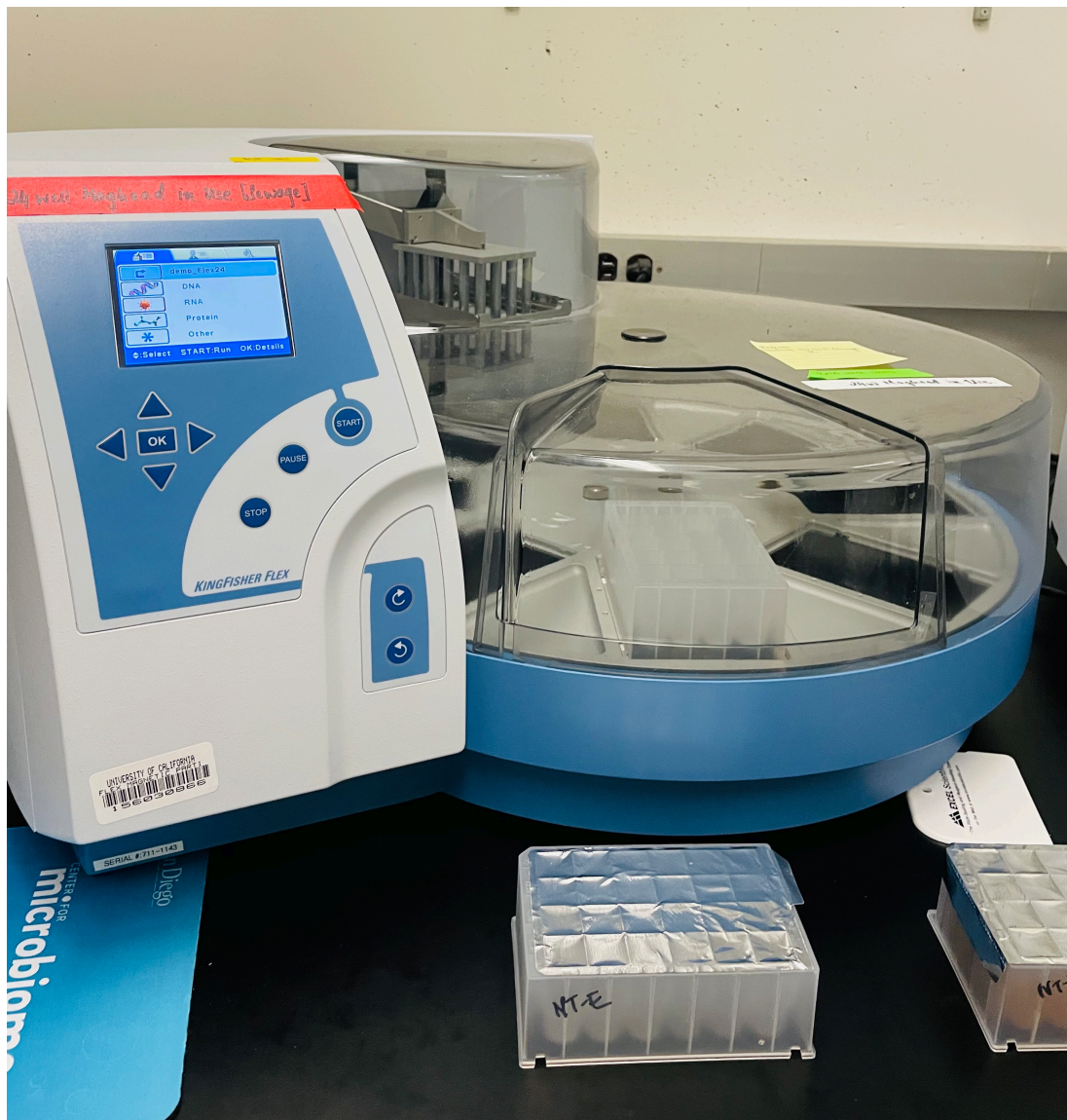
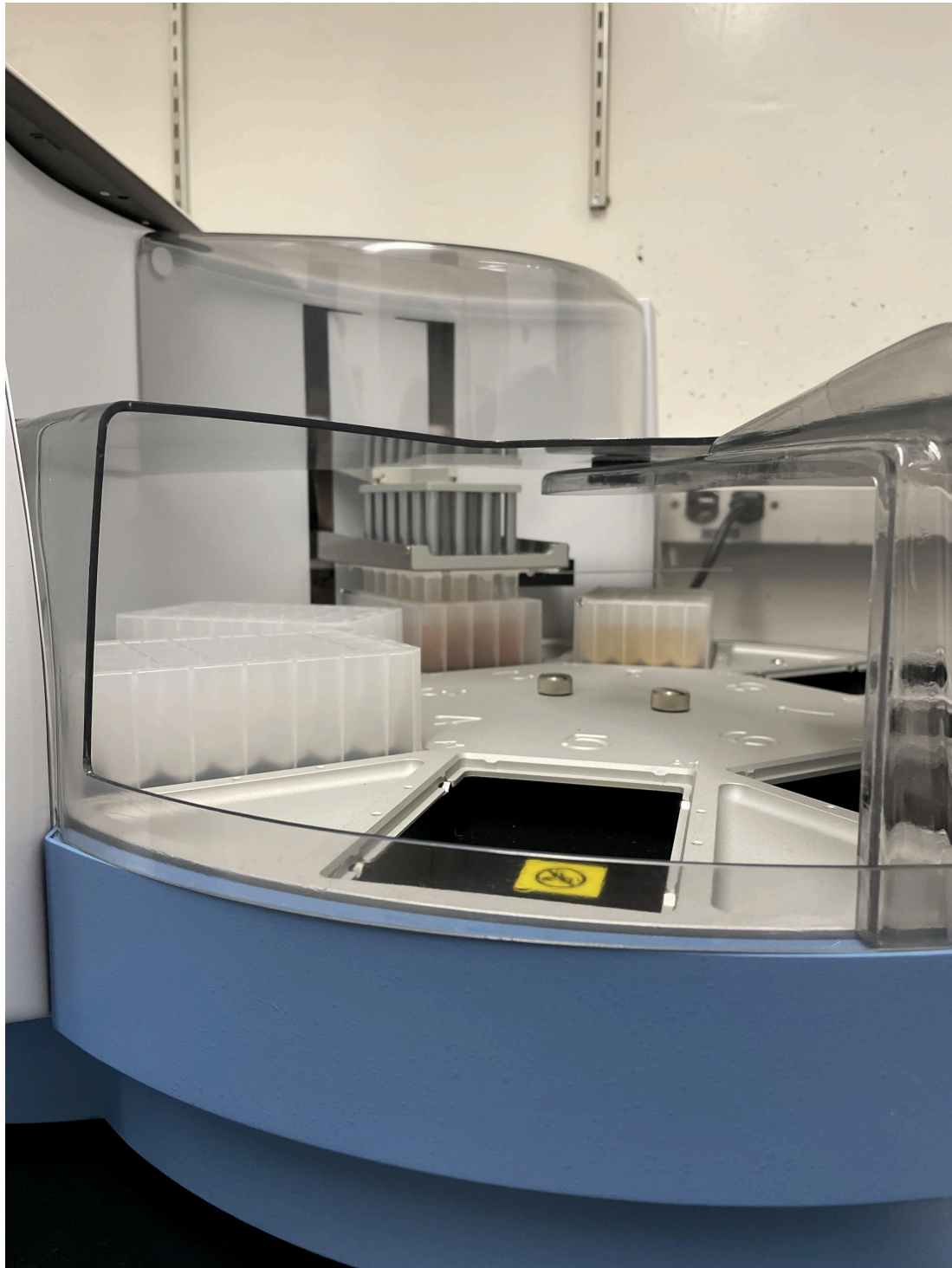


Plate 1: 5ml sample+ 75uL magnetic beads

- 3 Prepare the elution plate: Add  150 μL of 1X PBS and  800 μL of the  MagMAX[®]; Microbiome Lysis Solution **Thermo Fisher Catalog #A42361** to each of the wells of the 24 deep-well plate. 5m
- 4 Load the 3 plates into a KingFisher™ Flex **Automated Nucleic Acid Purification** system fitted with a 24 Deep-Well Head for KingFisher™ Flex Magnetic Particle Purification System (Cat. # 5400630 and 5400640).




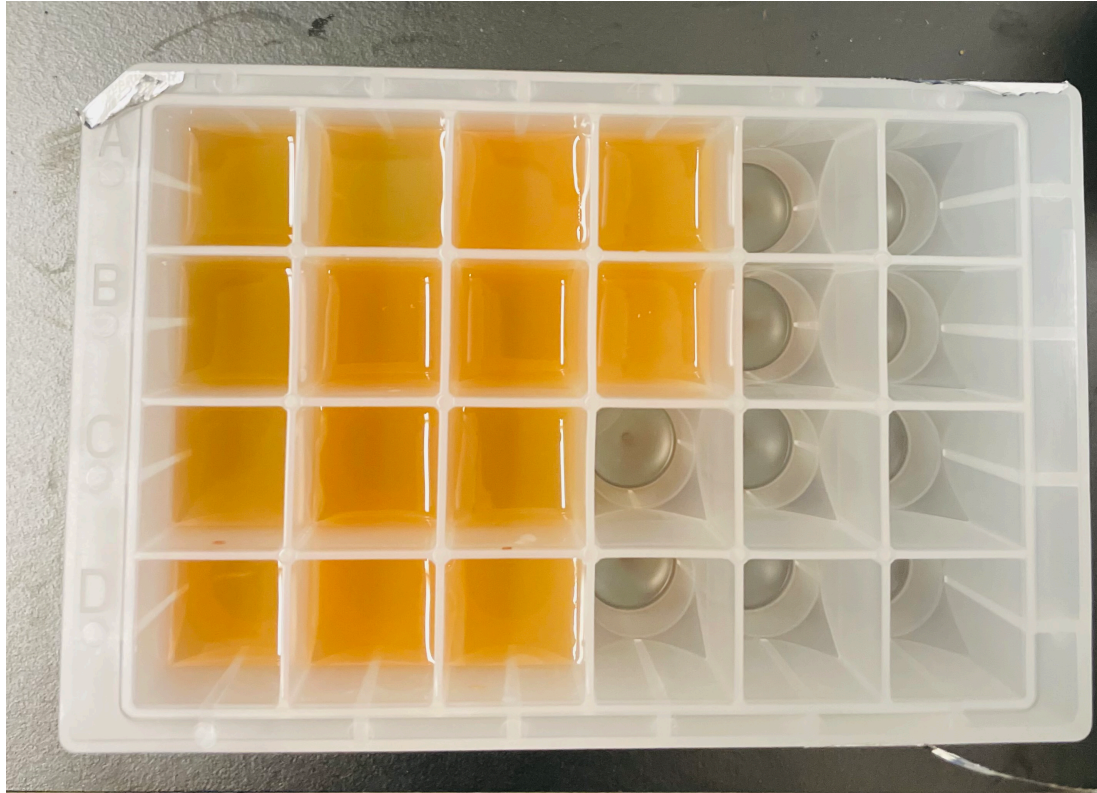






- 5 Download the program for the KingFisher Platform for the automated processing of 24 samples:  NanoTrap_Flex24_V1ck.bdz and load the program on the the equipment using the BindIt Software (Cat. # **5189009**).

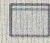

- 6 Once loaded, follow the steps on the the protocol to load the 3 plates into the system for extraction.

1h

Protocol summary			
NanoTrap_Flex24_48_Samples			
Layout	Tips	24 DW plate	
	Sample 1	24 DW plate	
	Sample 2	24 DW plate	
	Elution	24 DW plate	
	Sample 3	24 DW plate	
	Sample 4	24 DW plate	
	Elution 2	24 DW plate	
	Tips2	24 DW plate	
Steps	Bind sample 1, Drop Beads into Sample 2, Collect Remainder of Beads Sample 1, Bind Sample 2, Drop Beads into Elution, Collect Remainder of Beads Sample 2, Elution, Rinse beads into Sample 2, Collect leftover Beads from Elution, Mix1, Mix2, CollectBeads1, Mix3, Mix4, CollectBeads2, Mix5, Mix6, CollectBeads3		

Note

The protocol provided here includes an extended incubation step which can be reduced by half depending on the sample volume. The above protocol can also be increased for processing 48 samples.

- 7 The program processes the samples from the 2 sample plates and elutes the concentrated viral RNA into a single elution plate containing the lysis buffer.

Note

The Lysis buffer can vary depending on the Nucleic extraction kit used.

- 8 Nucleic acid extraction can then be performed using the



MagMAX[®] Microbiome Ultra Nucleic Acid Isolation Kit, with bead plate **Thermo Fisher Catalog #A42357**

36m



Note

The downstream extraction can be performed with any kit of choice. Here the MagMAX kit was used due to ease of integration into the KingFisher Flex system.