

Nov 08, 2022

B. burgdorferi enrichment from feeding ticks

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgqjrbovk5/v1>

Anne Sapiro¹, Jenny Zhang¹, Beth Hayes¹, Seemay Chou¹

¹University of California, San Francisco



Anne Sapiro

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgqjrbovk5/v1>

Protocol Citation: Anne Sapiro , Jenny Zhang, Beth Hayes, Seemay Chou 2022. B. burgdorferi enrichment from feeding ticks . protocols.io <https://dx.doi.org/10.17504/protocols.io.36wgqjrbovk5/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 31, 2022

Last Modified: November 08, 2022

Protocol Integer ID: 69418

Keywords: *Borrelia burgdorferi*, ticks, immunomagnetic separation, bb cells from ixodes tick, course of the tick bloodmeal, feeding tick, causative agent of lyme disease, tick bloodmeal, lyme disease, borrelia burgdorferi, ex vivo bb transcriptome, ixodes tick, global transcriptomic landscape, transmission during the tick stage, tick, encode protein, plasmids of the bb genome, full range of transcriptional change, tick stage, tick vector, transcriptional change, upregulated gene, gene expression, bb genome, majority of upregulated gene, key transcriptional regulator, dependent on key transcriptional regulator, bb for vertebrate infection, tick over the course, gene, transmitting bloodmeal

Abstract

Borrelia burgdorferi (*Bb*), the causative agent of Lyme disease, must adapt to vastly different environments as the bacterium cycles between the tick vector and a vertebrate host. During a bloodmeal, *Bb* migrates from the tick midgut to the salivary glands and changes its gene expression, priming *Bb* for vertebrate infection. These tick-specific events are dependent on key transcriptional regulators; however, the full range of transcriptional changes that occur inside of the tick over the course of several days during transmission are technically difficult to capture. We developed an experimental approach to enrich *Bb* cells from *Ixodes* ticks during a transmitting bloodmeal to longitudinally define their global transcriptomic landscape. We identified 192 genes that change expression at least twofold over the course of the tick bloodmeal, which were predominantly located on plasmids of the *Bb* genome. The majority of upregulated genes encode proteins found at the cell envelope or proteins of unknown function, including 45 upregulated outer surface lipoproteins embedded in the unusual protein-rich coat of *Bb*. The *ex vivo* *Bb* transcriptomes provide an important roadmap for investigating key determinants of *Bb* priming and transmission during the tick stage of its unique transmission cycle.

Materials

Dounce grinder (Kimble: 885303-0002) with 2 pestles
 Phosphate Buffered Saline (PBS) (3 mL/sample)
 anti-Bb antibody (Invitrogen: PA1-73004; RRID: AB_1016668) (2 µL/sample)
 Dynabeads Protein G (Invitrogen 10003D) (50 µL/sample)
 Magnet for 1.5 mL tubes (such as MagJET Separation Rack, 12 × 1.5 mL tube, Thermo Scientific MR02)
 Trizol (Invitrogen 15596018) (1.5 ml/sample)
 RNase-free non-stick 1.5 ml tubes (such as Invitrogen AM12450)

Safety warnings

 Use proper safety precautions for BSL2 agents and ticks.



Homogenize infected ticks

10m

- 1 Collect ticks for homogenization. We recommend using between 6-14 nymphal ticks depending on day of feeding. Ticks can be washed with water or surface sterilized with 1% bleach before beginning.
- 2 Add ticks and 500 μ L of PBS to the dounce grinder and homogenize using pestle A (large clearance pestle) 10 times up and down.
- 3 Remove pestle A and homogenize with pestle B (small clearance pestle) 5 times up and down.
- 4 Move homogenate to a 1.5 mL tube and increase the volume to 1 mL with PBS.
- 5 For an input sample: take out 50 μ L of the sample and add 500 μ L of Trizol. Vortex, incubate for 5 minutes at room temperature and then store at -20°C until RNA extraction.

Immunoprecipitation

1h 17m

- 6 Add 2 μ L of anti-Borrelia antibody to 1 ml (or 950 μ L) of homogenate and rotate at 4°C for 00:30:00 . 30m
- 7 [During 30 minute incubation in Step 6] Prepare Dynabeads Protein G by resuspending beads, then add 50 μ L per sample to a 1.5 mL tube. Place the tube on a magnet and remove the supernatant. Remove the tube from the magnet and resuspend the beads in 200 μ L of PBS.
- 8 After 30 minute incubation in Step 6, place the Dynabeads with PBS on the magnet and remove the supernatant. Add the homogenate + antibody mixture to the beads, resuspend, and rotate at 4°C for 00:30:00 . 30m
- 9 Place the tube on the magnet and remove the supernatant (the depleted fraction) into a new tube. 1m
- 10 After removing the depleted fraction, wash the beads 2 times with 1 mL of PBS, resuspending the beads each time. 2m



- 11 Place the tube on the magnet, remove the second PBS wash and resuspend the beads in 500 μ L Trizol. 1m
- 12 For a depleted sample, centrifuge the depleted fraction at  8000 x g, 00:07:00 to pellet the *Borrelia* and tick cells, remove ~900 μ L of supernatant and add 500 μ L of Trizol to the pellet. 7m
- 13 Incubate samples in Trizol for 5 minutes at room temperature before storing at  -20 $^{\circ}$ C until RNA extraction. 5m