

Sep 17, 2020

B-PER Lysis--CHEM 584

DOI

dx.doi.org/10.17504/protocols.io.bmg6k3ze

Ken Christensen¹

¹Brigham Young University



Ken Christensen

Brigham Young University

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bmg6k3ze

Protocol Citation: Ken Christensen 2020. B-PER Lysis--CHEM 584. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bmg6k3ze>

Manuscript citation:

<https://www.thermofisher.com/order/catalog/product/78243#/78243>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: September 17, 2020

Last Modified: September 17, 2020

Protocol Integer ID: 42238



Abstract

The Thermo Scientific B-PER Bacterial Protein Extraction Reagent enables mild extraction of proteins from bacteria (*E. coli*) without the need for mechanical disruption. The reagent may be used for soluble protein extraction and inclusion body purification from bacterial cell lysates. The B-PER Reagent with Enzymes is supplied with lysozyme and DNase I to improve the extraction efficiency of large (> 70 kDa) molecular weight proteins and proteins expressed in inclusion bodies.

Guidelines

Important Product Information

The B-PER Reagent extracts proteins from recently prepared cells and frozen cells. Extraction is typically more effective with frozen cells. The amount of reagent required depends on the amount of cell paste.

Important Product Information • The B-PER Reagent extracts proteins from recently prepared cells and frozen cells. Extraction is typically more effective with frozen cells. The amount of reagent required depends on the amount of cell paste.

B-PER Reagent effectively extracts soluble proteins from several common bacterial host strains and is especially suitable for the protease-defective expression host BL21 strains. If lysis is inefficient for a particular bacterial strain, freeze cells before extraction.



B-PER Lysis to extract soluble protein

- 1 Pellet bacterial cells by centrifugation at $5000 \times g$ in a tared tube for 00:10:00 .
- 2 Decant your spent LB and determine the mass of your cell pellet.
- 3 Freeze your pellet in the -20°C freezer overnight or quickly freeze your pellet in liquid nitrogen. When you are ready to proceed, thaw your cells at room temperature.
- 4 Add 4 mL of B-PER Reagent per gram of wet cell pellet. Pipette the suspension up and down until it is homogeneous.
- 5 Optional: Add 2 μL of lysozyme and 2 μL of DNase I per 1 mL of B-PER Reagent added.
- 6 Incubate 00:15:00 at room temperature.
- 7 Centrifuge lysate at high speed ($\sim 15,000 \times g$) for 00:05:00 to separate soluble proteins from the insoluble proteins.

Note

Note: If a large percentage of over-expressed protein remains in the pellet, the protein of interest might be expressed in inclusion bodies.