

Aug 28, 2019

Bacterial transformation

DOI

<https://dx.doi.org/10.17504/protocols.io.6taheie>



Laura Sánchez¹

¹AEGIS - Madrid iGEM Team 2019

AEGIS - Madrid iGEM 2019



Laura Sánchez

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6taheie>

External link: <https://www.addgene.org/protocols/bacterial-transformation/>

Protocol Citation: Laura Sánchez 2019. Bacterial transformation. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.6taheie>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 26, 2019

Last Modified: August 28, 2019

Protocol Integer ID: 27202

Keywords: transform, coli, heterologous DNA, plasmid, bacterial transformation transformation of competent escherichia coli culture, bacterial transformation transformation, competent escherichia coli culture, heterologous dna, external dna, introduction of external dna, cloning, heterologous protein expression, dna, piece of dna

Abstract

Transformation of competent *Escherichia coli* cultures with heterologous DNA.

This protocol allows the introduction of external DNA into competent *E. coli* cells. Typically, the piece of DNA consists of a plasmid, either for its cloning or for heterologous protein expression.

Guidelines

Work at 4°C unless otherwise indicated.

Sterile conditions are needed. Work always next to the flame.

Materials

MATERIALS

- ☒ 37°C Incubator
- ☒ Shaker incubator
- ☒ LB agar plates with the proper antibiotic(s) (eg. Kanamycin)
- ☒ Positive transformation control DNA
- ☒ Competent Cells
- ☒ SOC Media
- ☒ Ice & ice bucket
- ☒ 42°C water bath
- ☒ Microcentrifuge tubes **Eppendorf**



Safety warnings

⚠ Be careful about using disposable plastic gloves next to the flame. Burns are really painful.

Before start

Clean all the working surface with ethanol.

Make sure all the samples are placed on ice before start working.

The plates can be prepared and stored in advance.



Transformation

- 1 Take competent cells out of -80°C and thaw on ice (approximately 20-30 mins). 3m
- 2 Chill approximately 5 ng (2 µl) of DNA in a 1.5 ml microcentrifuge tube. 1m
** Please note: if draw from MoClo protocol samples, add the whole 20 µl of DNA*
- 3 Add 50 µl of competent cells to the DNA. 30s
Mix gently by pipetting up and down or flicking the tube 4–5 times to mix the cells and DNA.
Do not vortex.
- 4 Place the mixture on ice for 30 minutes. 3m
Do not mix.
- 5 Heat shock at 42°C for 45 seconds. 4s
Do not mix.
- 6 Add 950 µl of room temperature LB or SOC media to the tube. 30s
- 7 Place tube at 37°C for 60 minutes. Shake vigorously (250 rpm) or rotate. 6m
** Please note: For the duration and temperature of the heat shock step as well as for the media to be used during the recovery period, please follow the recommendations provided by the competent cells' manufacturer.*

Selection of transformants

- 8 Prepare selection plates:
 - LB + ampicillin 100 µg/mL
 - LB + kanamycin 100 µg/mL
 - LB + spectinomycin 50 µg/mL
- 9 Warm selection plates to 37°C. 1m
- 10 Spread 50–100 µl of the cells and ligation mixture onto the plates, and incubate O/N at 37°C. 1h



** Please note: Remember to add always a negative control plate (with medium without cells) to check contamination, and a positive control plate (same cells transformed with pUC) to check the efficiency of the process.*

** Tip: to ensure that all the positive results are plated, it is advisable to pellet the rest of the sample and resuspend it in a very low volume (50-100 µl), and then plate it.*

11 Check the samples for contamination. If clean, continue the protocol.

12 Choose three positive colonies of each plate.

Pick one with the inoculating loop and draw a strain in a new plate.

Do the same with the other two, using the same plate but leaving a separation between each strain.

Incubate O/N at 37°C.

** This will allow you to maintain and store in a separated plate the single clones chosen from each initial plate, and work with them.*

13 Prepare a liquid inoculum in a 15 mL Falcon tube of each of the three strains: pick a colony with the inoculating loop, resuspend it in 5 mL LB + antibiotic, and incubate 37°C O/N.

** Tip: if the chosen colonies are big enough, you may be able to do both steps 12 and 13 together. Selection of recombinants*

Selection of recombinants

14 Take a 1.5 mL sample of the Falcon tubes into a new eppendorf, and perform a miniprep.

15 Check the plasmids following the Plasmid Verification Protocol, and/or sequence them.

1h

1h

4m