

Feb 04, 2020 Version 1

BD FACS Aria II Shutdown V.1

DOI

dx.doi.org/10.17504/protocols.io.bameic3e

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Dillan Brown

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bameic3e

Protocol Citation: Allen Institute for Brain Science 2020. BD FACS Aria II Shutdown. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bameic3e>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 18, 2019

Last Modified: February 04, 2020

Protocol Integer ID: 31110

Keywords: FACS, Aria, Single-cell Sorting

Abstract

This protocol provides step by step instructions for shutting down the BD FACS Aria II and cleaning it of potentially hazardous materials.

Attachments



EQ0072_FACS_Aria_II...

40KB

