

Mar 18, 2020

Version 2

BD FACS Aria II Shutdown V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.bdvpi65n>

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bdvpi65n>

Protocol Citation: Allen Institute for Brain Science 2020. BD FACS Aria II Shutdown. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bdvpi65n>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 18, 2020

Last Modified: March 16, 2021

Protocol Integer ID: 34447

Keywords: FACS, Aria, Single-cell Sorting, bd facs aria ii shutdown, bd facs aria ii shutdown this protocol, bd facs aria ii, hazardous material, mental health, cleaning, national institute of mental health, fac

Abstract

This protocol provides step by step instructions for shutting down the BD FACS Aria II and cleaning it of potentially hazardous materials.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



EQ0072_FACS_Aria_II_...

40KB

