



Aug 27, 2024

Bioluminescence Imaging

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov191r2lr2/v1>

Santiago Unda^{1,2}, Michael G. Kaplitt^{1,2}

¹Department of Neurological Surgery, Weill Cornell Medical College, New York, NY 10065;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815, USA



Jacquelyn Haytayan

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov191r2lr2/v1>

Protocol Citation: Santiago Unda, Michael G. Kaplitt 2024. Bioluminescence Imaging. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ewov191r2lr2/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: August 27, 2024

Last Modified: September 23, 2024

Protocol Integer ID: 106521

Keywords: ASAPCRN, imaging, bioluminescence, bioluminescence stepwise imaging protocol in mice, bioluminescence imaging, bioluminescence imaging this protocol, ivis lumina series iii imaging system, bioluminescence, stepwise imaging protocol, imaging, luciferase, vagal afferent, expression in the vagal afferent, aav

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: 020608

Abstract

This protocol describes a bioluminescence stepwise imaging protocol in mice with AAV-luciferase expression in the vagal afferents using the Ivis Lumina series III imaging system.

Materials

- D-luciferin potassium salt (Xenogen)
- Isoflurane, USP (Covetrus)



Preparation and In vivo imaging

- 1 Induce anesthesia with isoflurane 2-3%.
- 2 Shave the neck, thoracic and abdominal area.
- 3 Prepare D-luciferin potassium salt (Xenogen) stock prepared in phosphate buffered saline (PBS) at 15mg/mL.
- 4 Inject D-luciferin solution (150mg/kg) intraperitoneal.
- 5 Image mice in supine position at 5-, 10- and 15-minutes post-luciferin injection.

Software protocol and analysis

- 6 Living imaging software (Ivis imaging) was configured to luminescent imaging mode with an exposure time of  00:00:00 and 5-minute delay.
- 7 Post-acquisition, the total flux of photons/radiance per second (s) in each pixel in the ROI area, in a square centimeter (cm²) of the tissue. This data is then multiplied by one steradian (sr), which result in average radiance displayed as (photons/s/cm²/sr).