

Aug 09, 2018

Version 1

Bodo saltans culture protocol V.1

DOI

dx.doi.org/10.17504/protocols.io.sh2eb8e

Fatma Gomaa¹, ZhuHong Li², Roberto Docampo², Peter Girguis¹, Virginia Edgcomb³

¹Harvard University; ²University of Georgia; ³Woods Hole Oceanographic Institution



Fatma Gomaa

Harvard University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.sh2eb8e>

Protocol Citation: Fatma Gomaa, ZhuHong Li, Roberto Docampo, Peter Girguis, Virginia Edgcomb 2018. Bodo saltans culture protocol. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.sh2eb8e>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 09, 2018



Last Modified: August 09, 2018

Protocol Integer ID: 14618

Keywords: agar medium for klebsiella, culture protocol medium recipe, paramecium medium solution, inoculation with bodo saltan, yeast extract, cerophyl powder, rye grass cerophyll, saltans food, ml of fresh medium, cerophyll to distilled water, cerophyll, bodo saltan, agar medium, ml from bodo culture, pneumoniae, fresh medium, preparation, saltan, bacterium, cold medium, atcc medium, liquid nitrogen, temperature mix, fresh bodo

Abstract

Medium recipe:

ATCC medium: 802 Sonneborn's Paramecium medium Solution 1 Rye grass Cerophyll:

Cerophyll*.....2.5 g

Distilled water.....1.0 L

Add cerophyll to distilled water and boil for 5 minutes.

Add 100 ml distilled water to compensate for evaporation.

Filter through Whatman #1 filter paper and add 0.5 g Na₂HPO₄.

Autoclave for 15 minutes at 121C.

1. *saltans* food (*K. pneumoniae*, or *E. coli*). I used only the *K. pneumoniae* so far

Agar Medium for *Klebsiella pneumoniae* ATCC-BAA-1705:

Agar.....20.0 g

Yeast extract.....4.0 g

Glucose.....0.16 g

Distilled water.....800.0 ml

Dispense in 5 ml amounts. Autoclave for 25 minutes at 121C. Slant. Bacterium, grown on solution 2, is added to solution 1 (Just add very little, few colonies) and incubated at 30C for 24 hours prior to inoculation with Bodo *saltans*.

- Cerophyll powder that works best for the *saltans* is the powder from Pines.

Culture Maintenance:

1. Prepare the bacterized *Bodo saltans* medium as described above.
2. Inoculate a T25 tissue culture flask (50ml) containing 20 to 25 ml of fresh medium with 1 to 2 ml from Bodo culture that is at or near peak density
3. Incubate horizontally at 18 to 22°C (room temperature can work fine) with cap screwed on not very tightly. 4. Subculture every 7 to 10 days.
4. I usually subculture 3 to 4 flasks every week

Cryopreservation:

Harvest and Preservation:

1. Harvest cells from a culture that is at peak density by centrifugation at 800 x g for 5 min. 2. Adjust the concentration of cells to 2×10^6 to 10^7 /mL in fresh medium (Important step, even for transfection).
2. Prepare a 20% (v/v) solution of sterile DMSO in fresh Bodo medium.
3. Add 2.0 mL of DMSO to an ice cold tube
4. Place the tube on ice and allow the DMSO to solidify (~5 min) and then add 8.0 mL of ice cold medium.
5. Invert several times to dissolve the DMSO.
6. Allow to warm to room temperature

7. Mix the cell preparation and the DMSO in equal portions. Thus, the final concentration will be 10^6 to 10^7 and 10% (v/v) DMSO. The time from the mixing of the cell preparation and DMSO stock solution before the freezing process is begun should be no less than 15 min and no longer than 30 min.
1. Dispense in 0.5 mL aliquots into 1.0 mL to 2.0 mL sterile plastic screwcapped cryules (special plastic vials for cryopreservation).
2. Place the vials in a controlled rate freezing unit. From room temperature cool at $1^{\circ}\text{C}/\text{min}$ to -40°C . If the freezing unit can compensate for the heat of fusion, maintain rate at $1^{\circ}\text{C}/\text{min}$ through the heat of fusion. At -40°C plunge into liquid nitrogen. Alternatively, place the vials in Nalgene 1°C freezing apparatus. Place the apparatus at -80°C for 1.5 to 2 hours and then plunge ampules into liquid nitrogen. (The cooling rate in this apparatus is approximately $1^{\circ}\text{C}/\text{min}$.)
3. To establish a culture from the frozen state place an ampule in a water bath set at $+35^{\circ}\text{C}$. Immerse the ampule to a level just above the surface of the frozen material. Do not agitate the ampule.
4. Immediately after thawing, do not leave in the water bath, aseptically remove the contents of the ampule and inoculate a T25 tissue culture flask containing 10 mL of Bodo medium bacterized with *Klebsiella pneumoniae* subsp. *pneumoniae* (ATCC® 700831).
5. Incubate horizontally with the cap screwed on tightly at 22°C

Attachments



Bodo saltans culture...

48KB

Troubleshooting

