


Nov 09, 2018

Version 1

Bounce PCR V.1

DOI

dx.doi.org/10.17504/protocols.io.vfne3me

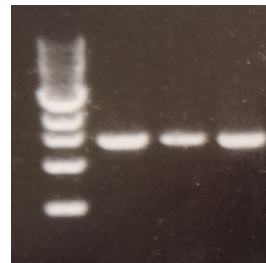

Sam Mugford¹, Saskia Hogenhout¹

¹John Innes Centre



Sam Mugford

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.vfne3me>

Protocol Citation: Sam Mugford, Saskia Hogenhout 2018. Bounce PCR. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.vfne3me>

Manuscript citation:

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 08, 2018

Last Modified: November 09, 2018

Protocol Integer ID: 17614

Keywords: PCR, Cloning, Synthetic Biology, optimisation free pcr for synthetic biology cloning application, free pcr method suitable for the rapid cloning, bounce pcr bounce pcr, bounce pcr, version of touchdown pcr, touchdown pcr, free pcr method, gene amplification, rapid cloning, molecular cloning, synthetic biology cloning application, cloning of multiple dna component, spurious priming during gene amplification, pcr, expensive alternative to traditional cloning, gene synthesis, advantage of primer, free pcr, multiple dna component, traditional cloning, biomolecular technique, synthetic biology project, cloning, most dna fragments without time, nucleic acids research, most dna fragment, increased primer, gateway recombination site, primer, target amplicon, amplification efficiency in later cycle, journal of biomolecular technique, gibson cloning, dna, early amplification cycle, amplification efficiency, vector overlaps for gibson cloning

Abstract

Bounce PCR: optimisation free PCR for synthetic biology cloning applications.

Synthetic biology projects require the cloning of multiple DNA components, and increasingly this is done through in-vitro gene synthesis. This remains an expensive alternative to traditional cloning by PCR. However, PCR often requires time-consuming optimisation. Bounce PCR is a largely optimisation-free PCR method suitable for the rapid cloning of multiple targets. Bounce PCR is a modified version of Touchdown PCR that takes advantage of primer-extension sequences -commonly used in molecular cloning - to successfully amplify most DNA fragments without time consuming optimisation. Touchdown (and related methods) employ a sequential lowering of the annealing temperature over successive cycles to minimise non-specific primer-binding during the early amplification cycles, but maximise amplification efficiency in later cycles when the target amplicon is more abundant (Don et al. 1991; Rowther et al. 2012). However, the lower annealing temperature in later cycles may still allow for non-specific amplification. Bounce PCR takes advantage of the increased primer melting temperature caused by the addition to the template of primer-extension sequences (for example: restriction enzyme recognition sites, Gateway recombination sites, or vector overlaps for Gibson cloning) to maintain the specificity in later cycles- without sacrificing efficiency- by increasing the annealing temperature again after an initial round of touchdown.

Don, R. H., Cox, P. T., Wainwright, B. J., Baker, K., & Mattick, J. S. (1991). 'Touchdown' PCR to circumvent spurious priming during gene amplification. *Nucleic acids research*, 19(14), 4008.

Rowther, F. B., Kardooni, H., & Warr, T. (2012). TOUCH-UP gradient amplification method. *Journal of biomolecular techniques : JBT*, 23(1), 1-3.



Guidelines

This protocol describes only the thermal cycling parameters of the PCR, and should work with your preferred DNA polymerase and protocol for setting up the reaction mixture. The advantage of Bounce PCR relies on the presence of primer-extension sequences, if your primers are an exact match to your template, without 5' extension sequences, then it is unlikely to offer an advantage over standard Touchdown PCR.

Materials

PCR mixture in suitable reaction tube.

Thermal cycler.

Troubleshooting



- 1 Initial denaturation: 94 degree C for 2 minutes
- 2 Denaturation 94 degree C for 30 seconds
- 3 Annealing: 60 degree C for 40 seconds (decreasing by 1 degree C per cycle for the first 15 cycles ("Touchdown"), then increasing by 0.6 degree C per cycle for the next 25 cycles ("Bounce"))
- 4 Extension: 72 degree C for 1 min per kb of target length
- 5 Repeat steps 2-4 for 40 cycles, varying the temperature of step 3 as described
- 6 Final extension step 72 degree C for 6 min