

Oct 14, 2021

Version 2

BPHL SARS-CoV-2 Tiled Amplicon Illumina Sequencing V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.by3tpynn>

Jason Blanton¹

¹Florida Department of Health, Bureau of Public Health Laboratories



Jason Blanton

Florida Department of Health, Bureau of Public Health Labora...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.by3tpynn>

Protocol Citation: Jason Blanton 2021. BPHL SARS-CoV-2 Tiled Amplicon Illumina Sequencing . protocols.io

<https://dx.doi.org/10.17504/protocols.io.by3tpynn> Version created by [Jason Blanton](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 14, 2021

Last Modified: October 14, 2021

Protocol Integer ID: 54099

Keywords: SARS-CoV-2, next generation sequencing, tiled amplicon, amplicon generation, tiled amplicon approach, sequencing workflow, amplicon approach, nextseq, bphl sar, using artic v3 primer, iseq, gohl protocol3, amplicon illumina, artic v3 primer, library pooling, miseq

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

The use of reagents, equipment, or services in this protocol is not an endorsement either stated or implied by the Florida Department of Health.

Abstract

This protocol details the Florida Department of Health's Bureau of Public Health Laboratories' (BPHL) wet lab portion of our SARS-CoV-2 next generation sequencing workflow. The method is a tiled amplicon approach using ARTIC V3 primers. The amplicon generation was adapted from the Matteson protocol¹. The library preparation is Illumina NexteraXT. Library pooling and normalization were adapted from the Gohl protocol³.

This protocol is for loading a MiSeq, but we have had equal success running on iSeqs and NextSeqs as well. Up to 96 libraries can be run on a MiSeq and up to 384 on a NextSeq.

Protocol materials

 SuperScript™ IV VILO™ Master Mix Thermo Fisher Scientific Catalog #11756500

RNA Extraction

- 1 Extract RNA from positive COVID-19 clinical specimens with the KingFisher Flex instrument using the Applied Biosystems™MagMAX™ Viral/Pathogen II (MVP II) Nucleic Acid Isolation Kit and its associated protocol.

cDNA Generation

- 2 cDNA from RNA (from any extraction method) is produced using  SuperScript™ IV VILO™ Master Mix **Thermo Fisher Scientific Catalog #11756500** with the following ratios, per sample:

 4 µL SuperScript IV VILO Master Mix

 6 µL Nuclease-free water

 10 µL Viral RNA

- 2.1 Tightly seal reaction wells, mix reaction components with plate mixer and spin down

- 2.2 Run the following thermal cycler protocol:

 25 °C 10 minutes

 50 °C 10 minutes

 85 °C 5 minutes

 4 °C ∞

- 2.3 Store cDNA at -20°C

ARTIC Amplicon Generation

- 3 ARTIC amplicons are produced by preparing two PCR reactions per sample (primer pool 1 is one reaction, primer pool 2 is the other). Per sample, combine below reagents in the listed ratios:



🧪 12.5 μ L Q5 HI-FIDEL 2X MASTER MIX 500 rxn

🧪 1 μ L IDT ARTIC V3 Primer Pool (20uM)

🧪 9 μ L nuclease-free water

🧪 2.5 μ L cDNA

3.1 Tightly seal reaction wells, mix reaction components with plate mixer and spin down

3.2 Run the following thermal cycler protocol:

🔥 98 °C 30 seconds

🔥 95 °C 15 seconds

🔥 64 °C 5 minutes

total of 35 cycles of steps 2 and 3

🔥 4 °C ∞

3.3 Combine PCR-amplified DNA from primer pool 1 and 2 together and dilute to 0.2-0.6 ng/ μ L (for Illumina)

3.4 Proceed to Illumina library prep method of choice (NexteraXT, Flex, Illumina compatible)

Library Quantification & Pooling

4 Quantify the DNA concentration of each clean library using the Qubit High Sensitivity dsDNA kit

4.1 Pool equal concentrations (e.g., 1-10 ng) of each library. Total volume does not matter

4.2 Concentrate using 0.7x AMPureXP beads (ex. for 240 μ L add 168 μ L of beads)

4.3 Allow binding at room temperature for at least 5 minutes before clearing with magnet

4.4 Wash beads 2x with 80% EtOH while still on magnet

- 4.5 Remove all EtOH and allow to pellet to dry for 5 minutes
- 4.6 Remove tube from magnet and add 20 μ L Tris-HCl pH 8.0 to pellet. Slowly pipette mix
- 4.7 Incubate at room temperature for at least 5 minutes before clearing with magnet
- 4.8 Check DNA fragment distributions of the pooled sample. Peak fragment size from 400 bp tiled amplicons with proper ligated adaptors should be $\sim$ 500nt
- 4.9 Quantify the DNA concentration of the pooled library using the Qubit High Sensitivity DNA kit

Note: At least 0.76 ng/ μ L is required to achieve 2 nM for library pooling. Libraries will need to be concentrated or re-amplified if less than this amount.
- 4.10 Convert DNA libraries from ng/ μ L to moles:
concentration [nM] = concentration [ng/ μ L] / ((ave.library size x 660)/1,000,000)

Example: if average size of library is 580 bp and concentration is 2.5 ng/ μ L:
 $(580 \times 660) / 1,000,000 = 0.382$
 $2.5 / 0.382 = 6.5$ nM
- 4.11 Dilute the pooled library to 2 nM in 10 mM Tris pH 8.0

Final Dilution and Loading

- 5 Add 10 μ L of a 2 nM library to 10 μ L of freshly prepared 0.2 N NaOH
- 5.1 Mix by flicking, spin down, incubate at room temperature for 5 minutes
- 5.2 Add 980 μ L prechilled HT1 to the tube containing denatured library. The result is 1 ml of a 20 pM denatured library
- 5.3 Add 240 μ L of the 20 pM denatured library, 348 μ L prechilled HT1, and 12 μ L of 20 pM denatured PhiX. Invert to mix, spin down



- 5.4 The result is 600 μ L of an 8 pM denatured library with 5% PhiX ready to be loaded on a MiSeq v3 cartridge

References

- 6
1. Nate Matteson, Nathan Grubaugh, Karthik Gangavarapu, Josh Quick, Nick Loman, Kristian Andersen 2020. PrimalSeq: Generation of tiled virus amplicons for MiSeq sequencing. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bez7jf9n>
 2. Grubaugh, ND. et al. An amplicon-based sequencing framework for accurately measuring intrahost virus diversity using PrimalSeq and iVar. *Genome Biology* 20,8 (2019) <https://genomebiology.biomedcentral.com/articles/10.1186/s13059-018-1618-7>
 3. Daryl Gohl 2020. SARS-CoV-2 Tailed Amplicon Illumina Sequencing. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bipikdke>
 4. Gohl DM, Garbe J, Grady P, Daniel J, Watson RHB, Auch B, Nelson A, Yohe S, Beckman KB, A rapid, cost-effective tailed amplicon method for sequencing SARS-CoV-2. *BMC Genomics* doi: [10.1186/s12864-020-07283-6](https://doi.org/10.1186/s12864-020-07283-6)