

Dec 02, 2024

## 🌐 Bulk RNAseq extraction

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6d1okvqe/v1>

Kimberly Paquette<sup>1</sup>, Andrew Pyman<sup>2</sup>, Jose Bras<sup>3</sup>


<sup>1</sup>Center for Alzheimer's and Related Dementias, National Institute on Aging, Bethesda, Maryland, USA;

<sup>2</sup>Van Andel Institute, Grand Rapids, Michigan, USA; <sup>3</sup>Regeneron Pharmaceuticals, Tarrytown, New York, USA



Kimberly Paquette

Center for Alzheimer's and Related Dementias, National Insti...

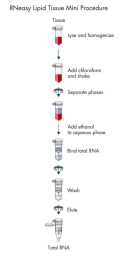

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.bp2l6d1okvqe/v1>

**Protocol Citation:** Kimberly Paquette, Andrew Pyman, Jose Bras 2024. Bulk RNAseq extraction. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.bp2l6d1okvqe/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** December 02, 2024

**Last Modified:** December 02, 2024

**Protocol Integer ID:** 113448

**Keywords:** ASAPCRN, RNA , Brain, Extraction, bulk rnaseq extraction rna extraction from frozen brain, bulk rnaseq extraction rna extraction, rneasy lipid tissue mini kit, frozen brain, using qiagen, extraction, rna extraction, using qiagen rneasy, qiagen rneasy, rna, qiagen

**Funders Acknowledgements:**

ASAP

Grant ID: ASAP-024407

## Abstract

RNA extraction from frozen brain using Qiagen's RNeasy Lipid Tissue Mini kit.

## Attachments



HB-0144-  
004\_HB\_RNY\_L...  
254KB



## Materials

|  | Reagent                   | Manufacturer             | Cat no. |
|--|---------------------------|--------------------------|---------|
|  | RNeasy Lipid Tissue Mini  | Qiagen                   | 74804   |
|  | Chloroform                | Millipore Sigma          | C2432   |
|  | Pure Ethanol              | Millipore Sigma          | E7023   |
|  | Qubit RNA Broad Range kit | Thermo Fisher Scientific | Q10210  |
|  | TissueRuptor II           | Qiagen                   | 9002755 |
|  | TissueRuptor II probes    | Qiagen                   | 990890  |
|  | RNase Zap                 | Thermo Fisher Scientific | AM9780  |
|  | RNase-Free DNase set      | Qiagen                   | 79254   |



## Cutting Brain Tissue

- 1 Manually cut 40-50mg of frozen brain tissue on dry ice. Weigh and place back in  -80 °C freezer until ready for extraction.

Note: do not let tissue thaw. Keep frozen at all times.

## RNA extraction

- 2 Using the Qiagen RNeasy Lipid Tissue Mini kit, prepare reagents and materials for extraction in a chemical hood. 

Note: use an RNase free environment and clean with RNase Zap.

- 3 Follow extraction steps outlined in Qiagen RNeasy Lipid Tissue Mini kit using the Tissue Ruptor II to homogenize frozen brain tissue.

- 3.1 Include on column DNase digestion step.

Note: keep extracted RNA on wet ice.

## Quality Control

- 4 Using the Qubit RNA Broad Range assay kit, measure the concentration of extracted RNA using  2 µL of sample.

- 5 Aliquot desired amount of sample to continue with for sequencing (  >500 ng ).

- 5.1 Optional: sequence with Illumina's TruSeq stranded riboreduced total RNA at the recommended 50M reads/sample.

## Storing RNA

- 6 Snap freeze each RNA sample with liquid nitrogen and store in  -80 °C freezer indefinitely.