

Nov 17, 2025

Calcium Imaging Protocol for hiPSC-derived Dopaminergic neurons

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ld65d7g5b/v1>

Judith Kreutzmann¹

¹Karolinska Institute

SOX6 mDA differentiation



Roberto Garcia Swinburn

Karolinska Institute Stockholm

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ld65d7g5b/v1>

Protocol Citation: Judith Kreutzmann 2025. Calcium Imaging Protocol for hiPSC-derived Dopaminergic neurons. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n92ld65d7g5b/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 09, 2025

Last Modified: November 17, 2025

Protocol Integer ID: 226859

Keywords: calcium imaging protocol for hipsc, calcium imaging protocol, imaging protocol to analyse calcium dynamic, dopaminergic neuron, derived dopaminergic neuron, analyse calcium dynamic, imaging protocol

Abstract

This is the imaging protocol to analyse calcium dynamics *in vitro* KOLF2.1J differentiated midbrain dopaminergic neurons

Protocol materials

 Fluo-4 Calcium Imaging Kit **Thermo Fisher Catalog #F10489**

 Pluronic® F-127 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P2443**



Preparation of Loading Solution

- 1 Prepare the dye-loading solution in Krebs-Ringer buffer:
 - 5 μ M Fluo-4 AM  Fluo-4 Calcium Imaging Kit **Thermo Fisher Catalog #F10489**
 - 0.625% Pluronic F-127
 -  Pluronic® F-127 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P2443**Mix thoroughly to ensure dye solubilization.



Dye Loading

30m

- 2 Add loading solution to the 35 mm culture dishes containing live cells.
- 3 Incubate at 37°C for 30 minutes **in the dark**.  37 °C

30m

Washing

- 4 After incubation, **wash cells twice** with fresh Krebs-Ringer buffer to remove excess dye.
- 5 Maintain cells in **fresh Krebs-Ringer buffer** at **37°C** during imaging.  37 °C

Calcium Imaging

- 6 Perform imaging using a Nikon CrEST X-Light V3 **spinning disk confocal microscope**:
 - Objective: 20x Air lens
 - Excitation wavelength: 480 nm
 - Sampling frequency: 1 Hz (one frame per second)Use **Nikon NIS-Elements software** to control the microscope and acquire time-lapse fluorescence data.



Data Analysis

- 7 Export image sequences from NIS-Elements.

- 8 Analyze calcium signals using the following tools:
- FIJI (ImageJ) for image preprocessing and intensity extraction.

Software

ImageJ/Fiji	NAME
Windows 7	OS
National Institutes of Health	DEVELOPER
http://wsr.imagej.net/distros/win/ij152-win-java8.zip	REPOSITORY
https://imagej.nih.gov/ij/	SOURCE LINK

- MATLAB (R2020b) for custom analysis scripts.

Software

MATLAB	NAME
Microsoft Windows 10	OS
Mathworks	DEVELOPER
Company Website	REPOSITORY
https://www.mathworks.com/products/matlab.html	SOURCE LINK

- FluoroSNNAP (GitHub: <https://github.com/tapan-patel/FluoroSNNAP>) for spike detection and activity analysis.