

Jul 09, 2025

Version 2

Calf-Intestinal Alkaline Phosphatase Treatment in situ-Killinger 2024 V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvoky7bl4o/v2>

Bryan Killinger¹, Solji Choi²

¹Rush University; ²Rush University Medical Center



Bryan Killinger

Rush University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvoky7bl4o/v2>

Protocol Citation: Bryan Killinger, Solji Choi 2025. Calf-Intestinal Alkaline Phosphatase Treatment in situ-Killinger 2024. protocols.io <https://dx.doi.org/10.17504/protocols.io.5qpvoky7bl4o/v2> Version created by **Bryan Killinger**

Manuscript citation:

Choi SG, Tittle T, Garcia-Prada D, Kordower JH, Melki R, Killinger BA. Alpha-synuclein aggregates are phosphatase resistant. bioRxiv [Preprint]. 2024 Apr 9:2023.11.20.567854. doi: 10.1101/2023.11.20.567854. PMID: 38645137; PMCID: PMC11030248.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: July 09, 2025

Last Modified: July 09, 2025

Protocol Integer ID: 222154

Keywords: sensitivity immunodetection of aggregated pser129, physiological from aggregated pser129, blocking buffer component bovine serum album, intestinal alkaline phosphatase treatment, physiological ps129 detection, intestinal alkaline phosphatase treatment in situ, phosphatase, buffer component bovine serum album, selective dephosphorylation, low level protease, aggregated pser129, success with mp biomedical, synuclein, mp biomedical, protease, sensitivity immunodetection, serum

Funders Acknowledgements:

NINDS

Grant ID: 1R01NS128467

Michael J. Fox Foundation

Grant ID: ASAP-024442

Michael J. Fox Foundation

Grant ID: MJFF-022480

Abstract

Alpha-synuclein phosphorylated at serine 129 (PSER129) occurs in two pools, non-aggregated (physiological) and aggregated (disease). This protocol allows for the selective dephosphorylation of non-aggregated PSER129 and enhances the specificity and sensitivity immunodetection of aggregated PSER129. Thus, this protocol can be used to differentiate physiological from aggregated PSER129.

For subsequent detection of non-aggregated-(physiological PS129) we have observed that the blocking buffer component bovine serum album (BSA) fraction V from several sources interferes with physiological PS129 detection. We have had success with MP Biomedicals AlbumiNZ Protease-reduced BSA, as well as omitting BSA from blocking buffers. We have not determined the source of interference, but it may include low level proteases or phosphatases.

Materials

Dilution media:

	A	B
	Nacl	150 mM
	Tris-HCl, pH 7.4	50 mM
	Triton-X100	0.5%

CIAP buffer:

	A	B
	Nacl	100 mM
	Tris- Hcl	50 mM
	Mgcl2, pH 7.9	10 mM

-  Alkaline Phosphatase Calf Intestinal (HC), 1,000u **Promega Catalog #M2825**



Day 1

1d 1h 10m

- 1 Wash free-floating tissue (3 × 10 minutes) in dilution media.



- Dilution media:

	A	B
	NaCl	150 mM
	Tris-HCl pH 7.4	50 mM
	Triton-X100	0.5%

- 1.1 Wash free-floating tissue 00:10:00 in dilution media (1/3).

10m



- 1.2 Wash free-floating tissue 00:10:00 in dilution media (2/3).

10m



- 1.3 Wash free-floating tissue 00:10:00 in dilution media (3/3).

10m



- 2 Incubate the samples with 1% Triton X-100 in DM for 00:10:00 .

10m



- 3 Wash in DM 00:10:00 .

10m



- 4 Wash the tissues in CIAP buffer (2×10 minutes).



- CIAP buffer:

	A	B
	NaCl	100 mM
	Tris- HCl	50 mM



A	B
MgCl ₂ , pH 7.9	10 mM

Autoclave and store Room temperature .

4.1 Wash the tissues in CIAP buffer 00:10:00 (1/2).

10m



4.2 Wash the tissues in CIAP buffer 00:10:00 (2/2).

10m



5 Incubate the tissues with CIAP at a dilution of 1:333 for 24:00:00 at 37 °C on a shaker.

1d



- CIAP concentration per bottle: 20 u/μl (Promega, Cat.# M2825).

- In 500 μL CIAP buffer, add 1.5 μL CIAP (30 units).

Day 2

3h 20m

6 Wash in DM (2 × 10 minutes).



6.1 Wash in DM 00:10:00 (1/2).

10m



6.2 Wash in DM 00:10:00 (2/2).

10m



7 Heat water bath to 80 °C - 85 °C 01:30:00 before the antigen retrieval step.

1h 30m

8 Place the dish containing sodium citrate buffer in the water bath and heat it for 00:10:00 .

10m



- Sodium Citrate Buffer, pH 6.0 (1L): 2.94 g Sodium citrate-Trisodium salt (Dihydrate) in 1000 mL DI water. pH 6.0. Add 0.5 mL Tween-20. Mix well.

9 Wash the tissues in sodium citrate buffer 00:10:00 .

10m



10 Incubate the tissues in the heated sodium citrate buffer for 00:30:00 .

30m



11 Cool the dish containing tissues to Room temperature (at least 00:20:00).

20m

12 Wash in DM for 10 min x 2 times.



12.1 Wash in DM for 00:10:00 (1/2).

10m



12.2 Wash in DM for 00:10:00 (2/2) .

10m



13 Tissues are now ready for downstream assays.