

Jun 20, 2018

Version 1

Carbon isotope descrimination - part I - trapline - v0.1 V.1

DOI

dx.doi.org/10.17504/protocols.io.qywdxxe

Stephen Driever¹, Steven J Burgess¹

¹University of Illinois at Urbana-Champaign



Steven J Burgess

University of Illinois at Urbana-Champaign

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.qywdxxe>

Protocol Citation: Stephen Driever, Steven J Burgess 2018. Carbon isotope descrimination - part I - trapline - v0.1. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.qywdxxe>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: June 14, 2018



Last Modified: June 20, 2018

Protocol Integer ID: 13046

Keywords: Photosynthesis, Plant physiology, carbon isotope discrimination, carbon isotope discrimination during photosynthesis, measurement of carbon isotope discrimination, photosynthesis, trapline, gas for measurement, collecting gas

Abstract

Protocol for collecting gas for measurement of carbon isotope discrimination during photosynthesis.

See the following references for more background:

H. Griffiths (1993) **Photosynthesis and Production in a Changing Environment** pp 181-1927

Farquhar et al. (1989) Annu. Rev. Plant Physiol. Plant Mol. Bioi. 1989. 40:503-37

Troubleshooting

Setting up IRGA

- 1 Set the following environmental variables in the LICOR LI-6800

CO2S	400 ppm
Txchg	25 °C
H2OR	10-14 mbar (aim for Tleaf 25 °C)
PAR	2000 (100% Red)
Flow	250 ml min ⁻¹
Fan speed	14500 rpm
deltaP	0.1 kPa

Note

Control temperature on block (Txchg), not leaf otherwise photosynthetic rate will start to oscillate due to constant adjustment of stomata.

TLeaf should ideally be within +/- 0.03oC of 25oC as measurement is temperature sensitive. 25oC is the value that C3 photosynthetic calculations are parameterized for. You can manipulate with humidity or Txchg on reference


Preparing plant for measurement

- 2
 - Clamp leaf in cuvette, under a bright ambient light.
 - Adjust H2OR so that Tleaf is around 25 °C. This will need adjusting when leaf is reaching stability in the cuvette.



Note

For tobacco a 4 week old plant prior to flowering is in the ideal state (health example below).

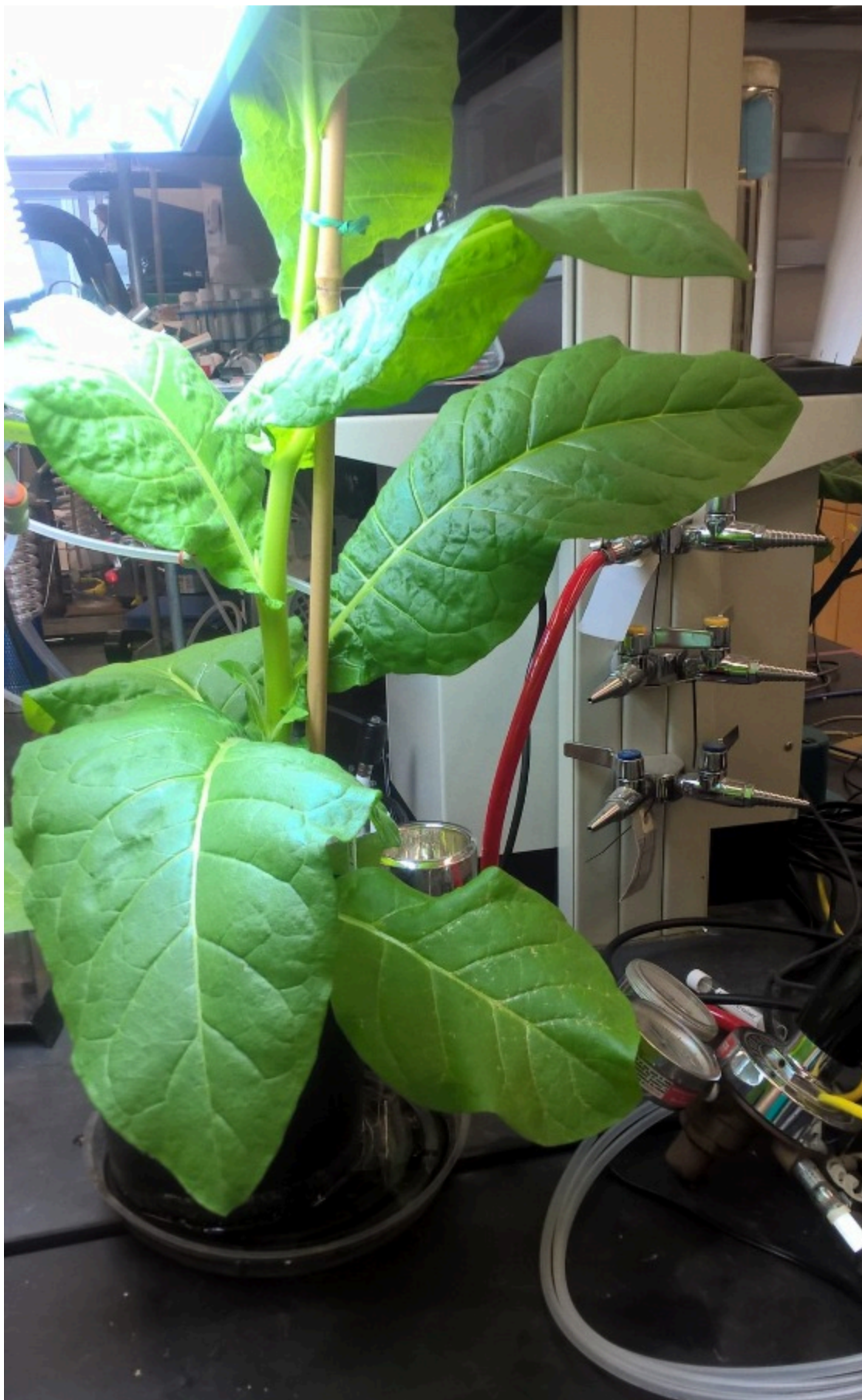



Make sure not to stress the plant by introducing tension when clamping.

Illuminate from above (at least $200\mu\text{E}$), this makes sure the surrounding leaf tissue is photosynthesising to reduce chance of diffusion of metabolites influencing measurements, this is more important for homobaric leaves such as tobacco.

It helps to let the plant acclimate in the lab for $\sim 24\text{h}$ prior to measurement.




Wait with measurements and CO_2 trapping until complete steady state has been reached.

- Assimilation rate (A) and steady state fluorescence do not change over time (10 min window)
- Stomatal conductance (g_s) does not change over time (10 min window)

This can take $\sim 2\text{h}$ depending on the plant. Check regularly (\sim every 15 mins after the first half hour)

Note that mesophyll conductance can be altered by a whole range of factors (cell size, leaf expansion, age of leaf etc etc) so care should be taken in performing comparisons and should try to keep to plants at a consistent developmental stage and growth conditions.

 01:00:00

- 3 When plant has reached a steady state switch inlet gas to 2% oxygen. Use a flow controller to moderate gas use, setting flow to ~ 6 .



Change the constant value in LICOR

Note

2% O₂ is used to suppress photorespiration simplifying downstream calculations of carbon flux. Note this will change assimilation and photosynthesis values which will take time to stabilize again, you may witness oscillations in A and F which will dampen down over time.




When there is no flow the value should be ~0.5 or less



If the cylinder is running low order a replacement, always keep a backup cylinder to avoid running out of gas



 01:00:00

Setting up preline

4 Turn on cold finger

Note

This will be used to cool down acetone to below -70°C as required for sublimation of frozen CO_2




There is only an ON/OFF setting.

- 5 Place cold finger in dewar of acetone and wait to cool down ($>-70^{\circ}\text{C}$), it should turn milky white

Note



Cover the top of the dewar with foli to prevent ice forming and falling into the acetone. Do not rest the finger on the bottom of the dewar otherwise an ice lump forms and it will stick

 01:00:00

- 6 Check temperature of acetone is below -70°C

Note



- 7 Fill two dewars 3/4 full with LN2

Note




These will be used to condense gases for sample collection. This can be done whilst waiting for the acetone to cool down



- 8 Open glass valves and wait until vacuum gauge reads around 5.5 mbar on each prep line. Adjust by opening/closing pump valves if required.

Note

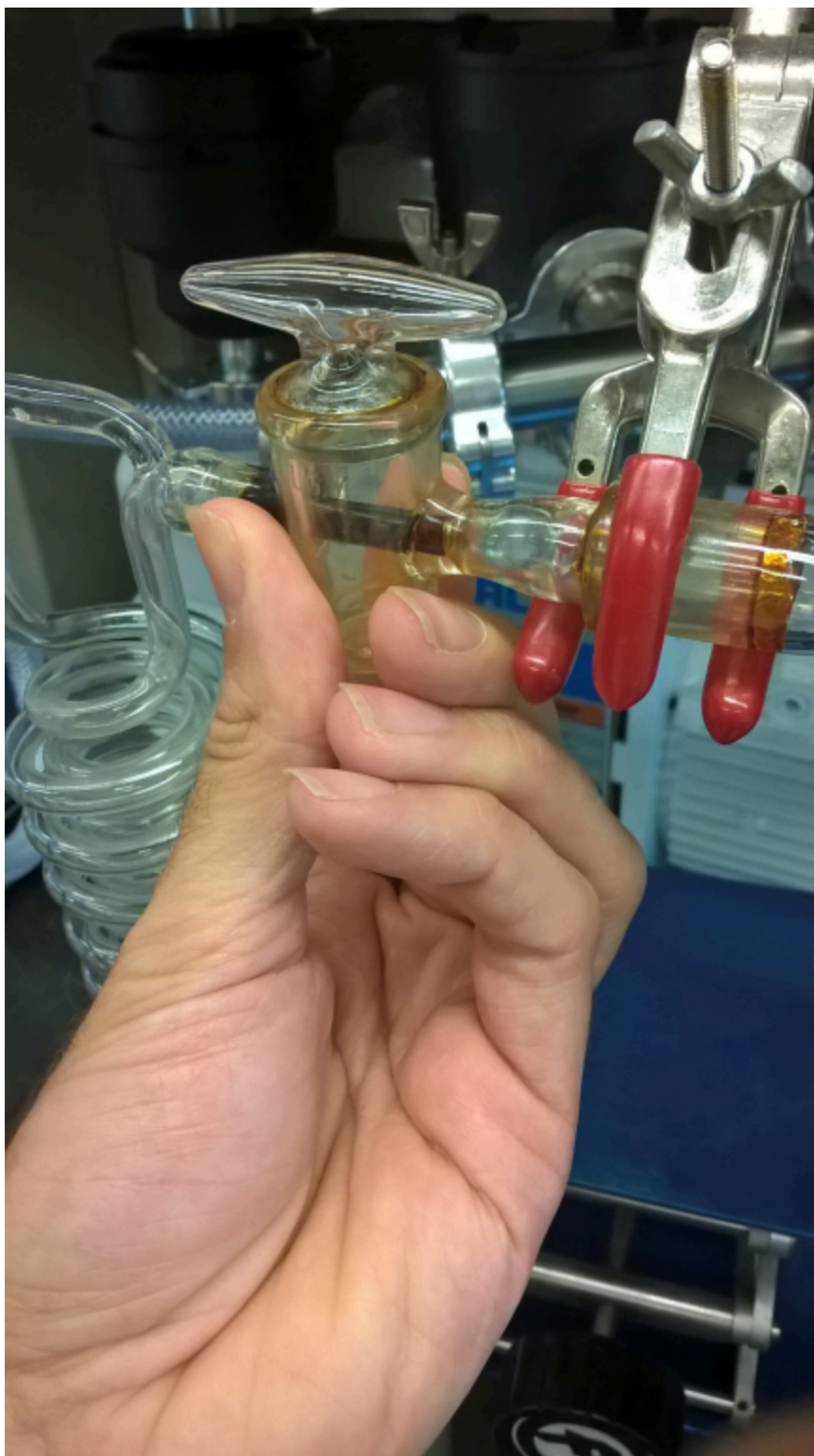

Provide support from below when opening/closing valve. Apply downward pressure while twisting to minimize the chance of introducing leaks (this is the most common source in the system).



Adjust the vacuum pressure by opening the valve to the pump (red arrow). This ensures that approximately the same amount of gas is collected for reference and sample and ensures consistency between measurements



- 9 Check that both collection tubes are open
- 10 Check if leaf is still in stable state (both A and gs)



Start Measurement


11 Match LI-6800 manually

Check settings

- automatch OFF
- Fluorescence OFF

12 Immerse both traps in LN2 (2 dewars)


Note

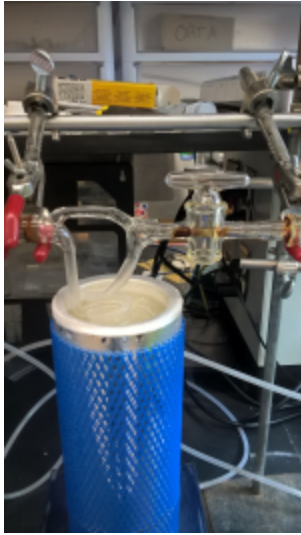

The distillation tube should ideally be covered close to the top



13 Set timer and start logging every minute min for 5 min (LI-6800, autoprogram)

 05:00:00


Note



- 14 Take fluorescence measurement (F_m' 's)
- 15 Close the glass valves, watch the vacuum drop to low 10^{-2} or high 10^{-3} E.

Note

Twist perpendicular to the flow direction otherwise there can be leaks, check the pressure gauge

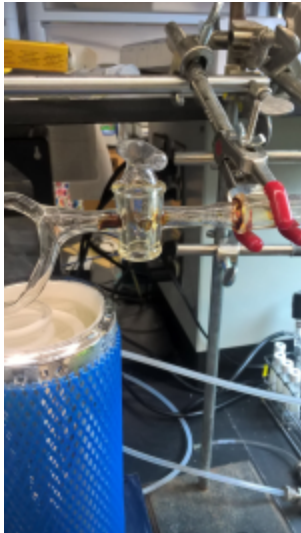


- 16 Close the pump valves.

Note




- 17 For each prep line (one after the other) make sure the collection tubes are open.

Note

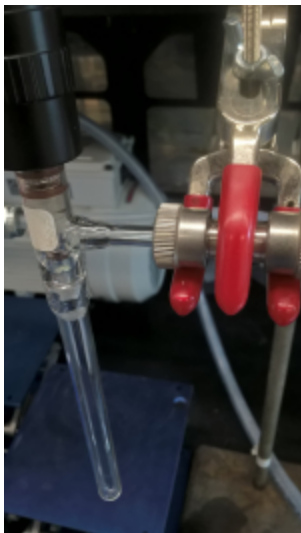

- 18 In a single motion remove distillation tube from LN2 and submerge in cold acetone to sublimate gases


Note

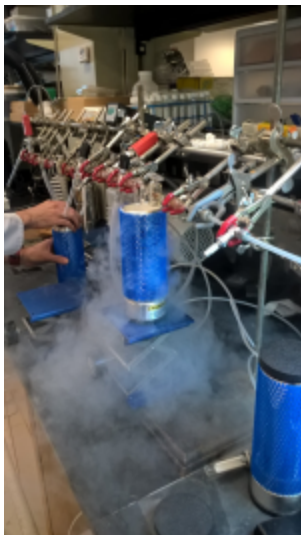

The pressure in the prep line should rise as gas is released

- 19 Quickly submerge the collection tube in LN2 to condense the released gas

Note

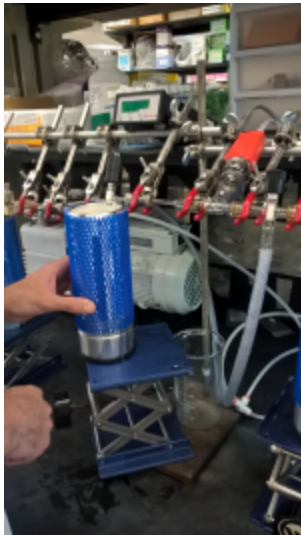

The pressure should fall again as gas is drawn down into the collection tube. By following this procedure it is possible to collect gases and remove the water vapour from samples which would interfere with measurements



Collection tube should be dipped in LN2, covering at least the bottom 2cm

20 Close collection tube

21 Remove acetone and liquid N dewars

Note

When the vacuum stabilizes all gas has been condensed again, this is the point when you can close tube and remove from LN2



Collect drips of acetone from distillation tube in beaker



- 22 Remove collection tube from prep line and leave to thaw in rack
Put new (empty and open) collection tube in place and open pump valve

Note



Opening the pump valve cleans out the prep line and residual air in the new collection tube

- 23 Close low O₂ cylinder and adjust constant value in LICOR