

May 08, 2019



Case - Plasma Insulin Measurement by ELISA

DOI

dx.doi.org/10.17504/protocols.io.yesftee



Henri Brunengraber¹

¹Case Western Reserve University

Mouse Metabolic Phenotyping Centers Tech. support email: info@mmpc.org



Lili Liang

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account





DOI: https://dx.doi.org/10.17504/protocols.io.yesftee

External link: https://mmpc.org/shared/document.aspx?id=270&docType=Protocol

Protocol Citation: Henri Brunengraber 2019. Case - Plasma Insulin Measurement by ELISA. protocols.io

https://dx.doi.org/10.17504/protocols.io.yesftee



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 21, 2019

Last Modified: May 08, 2019

Protocol Integer ID: 20658

Keywords: Mouse plasma, Insulin ELISA kit, mouse insulin elisa kit, plasma insulin measurement by elisa summary, plasma insulin measurement, radioactive quantification of insulin, insulin, mouse plasma, elisa summary

Abstract

Summary:

Mouse Insulin ELISA kit is used for the non radioactive quantification of insulin in mouse plasma.

Materials

MATERIALS

Mouse Sensitive Insulin ELISA Mercodia Catalog #10-1247-01

Mouse Ultrasensitive Insulin ELISA Mercodia Catalog #10-1249-01

Reagents and Materials:

Reagent/Material	Quantity Required	Vendor	Stock Number
Mouse Sensitive Insulin ELISA	10ul serum/plasma per well	Mercodia	nr 10-1247-01
Mouse Ultrasensitive Insulin ELISA	25ul Serum/plasma per well	Mercodia	10-1249-01

Note:

Mouse Sensitive Insulin ELISA, Cite this, (Mercodia Cat# 10-1247-01, RRID:AB_2783837)

Troubleshooting



1 **Protocol:**

- 1. Plasma or serum are isolated from mice fasted or after glucose injection. Heparin or EDTA coated tubes are acceptable for plasma fractionation.
- 2. Insulin is measured according to the manufacturer's directions.

2 **Manufacturer's Protocol below:**

MOUSE INSULIN ELISA KIT

96-Well Plate (Cat. #10-1247-01)



3 Principle of procedure (provided by manufacturer)

Mercodia Mouse Insulin ELISA is a solid phase two-site enzyme immunoassay. It is based on the direct sandwich technique in which two monoclonal antibodies are directed against separate antigenic determinants on the insulin molecule. During incubation insulin in the sample reacts with peroxidase-conjugated anti-insulin antibodies and antiinsulin antibodies bound to microplate wells. A simple washing step removes unbound enzyme labeledantibody. The bound conjugate is detected by reaction with 3,3',5,5'tetramethylbenzidine. The reaction is stopped by adding acid to give a colorimetric endpoint that is read spectrophotometrically. This assay is a Sandwich ELISA based, sequentially, on:

- 1) capture of insulin molecules from samples to the wells of a microtiter plate coated by pre-titered amount of a monoclonal mouse anti-rat insulin antibodies and the binding of biotinylated polyclonal antibodies to the captured insulin,
- 2) wash away of unbound materials from samples,
- 3) binding of horseradish peroxidase to the immobilized biotinylated antibodies,
- 4) wash away of free enzyme conjugates, and
- 5) quantification of immobilized antibody- enzyme conjugates by monitoring horseradish peroxidase activities in the presence of the substrate 3,3',5,5'-tetramethylbenzidine. The enzyme activity is measured spectrophotometrically by the increased absorbency at 450 nm, corrected from the absorbency at 590nm, after acidification of formed products. Since the increase in absorbency is directly proportional to the amount of captured insulin in the unknown sample, the latter can be derived by interpolation from a reference curve generated in the same assay with reference standards of known concentrations of rat insulin.