


Dec 07, 2023

🌐 CD8 Cell Density In Substantia Nigra And Cerebral Peduncle Image Analysis

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvo3roxv4o/v1>

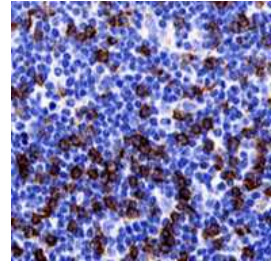

Hemanth Ramesh Nelvagal^{1,2}, Toby Curless^{1,2}, Zane Jaunmuktane^{1,3,2}

¹Department of Clinical and Movement Neurosciences, UCL Institute of Neurology, Queen Square, London WC1N 3BG, UK;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20 815, USA;

³Queen Square Brain Bank for Neurological Disorders (QSBB), 1 Wakefield Street, London WC1N 1PJ, UK



Hemanth Ramesh Nelvagal

University College London, University of London

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvo3roxv4o/v1>

Protocol Citation: Hemanth Ramesh Nelvagal, Toby Curless, Zane Jaunmuktane 2023. CD8 Cell Density In Substantia Nigra And Cerebral Peduncle Image Analysis. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5qpvo3roxv4o/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 07, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 86040

Keywords: ASAPCRN, Human Brain, Substantia Nigra, Cerebral Peduncle, CD8, Parkinson's disease, QuPath, NZConnect, Image Analysis, Pathology, cerebral peduncle image analysis qupath, cd8 cell density in substantia nigra, cd8 cell density, cd8 density in the substantia nigra, cd8 density, bioimage analysis software, stained brain section, cerebral peduncle, digital pathology, brain section, image analysis, whole slide image analysis, brain

Funders Acknowledgements:

The Michael J. Fox Foundation for Parkinson's Research (MJFF) and the Aligning Science Across Parkinson's (ASAP) Initiative
Grant ID: ASAP-000478

Abstract

QuPath is a bioimage analysis software designed for digital pathology and whole slide image analysis. This protocol describes how to measure CD8 density in the substantia nigra and cerebral peduncle using haematoxylin and DAB-stained brain sections.

Materials

- CD8 IHC-stained sections
- NZConnect (Hamamatsu)
- StarDist
- QuPath

Annotation

- 1 Manually annotate the substantia nigra and cerebral peduncle on NZConnect (Hamamatsu), a web-based whole-slide image (WSI) viewer.
- 2 Download the annotations using a Python script then import into QuPath [\[1\]](#) using a Groovy script.

QuPath Deconvolution and CD8 Density Measurement

- 3 In QuPath, set the colour deconvolution to facilitate the detection of haematoxylin and DAB staining on CD3 IHC-stained sections.
- 4 Segment all cell nuclei using StarDist [\[2\]](#) via the QuPath StarDist extension[\[3\]](#), follow with an object classifier to classify CD8-positive cells.
- 5 Calculate CD8-positive cell density was calculated by the number of CD8-positive cells divided by the area of the region of interest (CD8-positive cells per mm²).

Note

References

[1] Bankhead, P., Loughrey, M.B., Fernández, J.A. et al. QuPath: Open source software for digital pathology image analysis. *Sci Rep* **7**, 16878 (2017). <https://doi.org/10.1038/s41598-017-17204-5>

[2] Schmidt, U., Weigert, M., Broaddus, C., & Myers, G. (2018, September). Cell detection with star-convex polygons. In *International Conference on Medical Image Computing and Computer-Assisted Intervention* (pp. 265-273). Springer, Cham. <https://arxiv.org/abs/1806.03535>

[3] StarDist extension for **QuPath** <https://github.com/qupath/qupath-extension-stardist>



Protocol references

- [1] Bankhead, P., Loughrey, M.B., Fernández, J.A. *et al.* QuPath: Open source software for digital pathology image analysis. *Sci Rep* **7**, 16878 (2017). <https://doi.org/10.1038/s41598-017-17204-5>
- [2] Schmidt, U., Weigert, M., Broaddus, C., & Myers, G. (2018, September). Cell detection with star-convex polygons. In *International Conference on Medical Image Computing and Computer-Assisted Intervention* (pp. 265-273). Springer, Cham. <https://arxiv.org/abs/1806.03535>
- [3] StarDist extension for **QuPath**
<https://github.com/qupath/qupath-extension-stardist>