

Feb 13, 2026

Cell apoptosis assay

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ld1jzn15b/v1>

Guangli Suo¹

¹CAS Key Laboratory of Nano-Bio Interface, Suzhou Institute of Nano-Tech and Nano-Bionics, Chinese Academy of Sciences, Jiangsu, 215123, China.

MTS



Guangli Suo

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ld1jzn15b/v1>

Protocol Citation: Guangli Suo 2026. Cell apoptosis assay. protocols.io
<https://dx.doi.org/10.17504/protocols.io.n92ld1jzn15b/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: February 13, 2026

Last Modified: February 13, 2026

Protocol Integer ID: 243183

Keywords: cell apoptosi, quantitative assessment of apoptotic cell population, apoptosi, apoptotic cell population, flow cytometry, using fluorescein isothiocyanate, fluorescein isothiocyanate, propidium iodide, assay

Abstract

Quantitative assessment of apoptotic cell populations by flow cytometry using fluorescein isothiocyanate (FITC)-conjugated annexin V and propidium iodide (PI) dual staining.



Cell apoptosis assay

- 1 The apoptosis of cells cultured on MWC-chips was assessed using Annexin V-FITC and propidium iodide (PI) double staining method (Vazyme). For cell lines, BRC cell lines (MDA-MB-231) or normal human foreskin fibroblasts (HFF) cell lines were seeded in the MWC-chips and cultured in Sel-Medium or normal cultural condition for 3 days. Subsequently, cells were harvested, trypsinized and stained with Annexin V-FITC and PI. For patient-derived cells, clinical BRC and adjacent tissues were dissociated into single cells which were then seeded onto MWC-chips for 2 days of culture in Sel-Medium followed by 3 days of culture in SG-Medium according to the MTSs generation protocol. The resulting MTSs and RCs were collected, trypsinized and stained with Annexin V-FITC and PI. Finally, the stained cells were analyzed by flow cytometry (CytoFLEX S, Coulter Beckman). Cells negative for PI but positive for Annexin V were considered apoptotic, while cells positive for both PI and Annexin V were classified as necrotic.