

Feb 13, 2019

Cell dissociation from airway biopsies with cold-active protease for single-cell RNA-seq

DOI

dx.doi.org/10.17504/protocols.io.x3efqje

Laure-Emmanuelle Zaragosi¹, Pascal Barbry¹


¹Université Côte d'Azur, CNRS, IPMC, 06560 Valbonne, France

Human Cell Atlas Metho...



Laure-Emmanuelle Zaragosi

Institut de Pharmacologie Moléculaire et Cellulaire

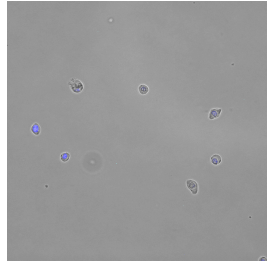

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.x3efqje

Protocol Citation: Laure-Emmanuelle Zaragosi, Pascal Barbry 2019. Cell dissociation from airway biopsies with cold-active protease for single-cell RNA-seq. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x3efqje>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 12, 2019

Last Modified: February 13, 2019

Protocol Integer ID: 20294

Keywords: bronchial biopsy, airway epithelium, single-cell, dissociation, cold-active protease



Abstract

This protocol provides details on the cell dissociation that should be performed to obtain single-cell suspensions from airway biopsies.

Biopsies may come from tracheal, bronchial or nasal epithelium.

Cell dissociation is performed at 4°C to avoid gene expression alterations and maximize viability.

The typical cell number recovery is 40 000 cells for one biopsy.

Cell suspensions are suitable for single-cell RNA-sequencing protocols.

Guidelines

Storage Conditions of Reagents

Reagent	Storage Condition
HBSS	4°C
20 mM EDTA	room temperature
BSA (Sigma, A8806)	4°C
Protease from <i>Bacillus Licheniformis</i> (Sigma, P5380)	Store 100 µL aliquots (100 mg/mL) in DPBS at -80°C
Hoechst 33342 (10 mg/mL)	4°C
NucGreen™ Dead 488 ReadYProbes™ Reagent	room temperature

Required Equipment

Equipment	Supplier	Catalog no.
Countess II	Thermo	AMQ AF10



FL auto mate d cell count er	Fishe r Scien tific	00
---	------------------------------	----

The protocol workflow is as follows:

1. Perform airway biopsies in the desired zone
2. Dissociation: mince with scalpel then triturate on ice in dissociation buffer
3. Remove red blood cells if necessary
4. Prepare cells for Chromium/DropSeq

All steps should be performed on ice or at 4°C

Use wide-bore 1 mL pipet tip for all biopsy transfers.



Materials

MATERIALS

✕ EDTA

✕ 23G Needles **Catalog #4657667**

✕ Protease from *Bacillus Licheniformis* **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P5380**

✕ Quick-Read 10 Chamber Slide **Globe Scientific Catalog #3805**

✕ Countess™ Cell Counting Chamber Slides **Catalog #C10314**

✕ DPBS no calcium, no magnesium **Invitrogen - Thermo Fisher Catalog #14190136**

✕ 21G needle **VWR International (Avantor) Catalog #BD-305165**

✕ HBSS **Gibco - Thermo Fisher Scientific Catalog #14060040**

STEP MATERIALS

✕ Quick-Read 10 Chamber Slide **Globe Scientific Catalog #3805**

✕ Ammonium Chloride Solution 100 mL **STEMCELL Technologies Inc. Catalog #7800**

✕ Flowmi cell strainer **Catalog #H13680-0040**


✕ Hoechst 33342, Trihydrochloride, Trihydrate - 10 mg/mL Solution in Water **Thermo Fisher Scientific Catalog #H3570**

✕ NucGreen™ Dead 488 ReadyProbes™ Reagent **Thermo Fisher Scientific Catalog #R37109**

Protocol materials

 23G Needles **Catalog #4657667**

 Flowmi cell strainer **Catalog #H13680-0040**

 Hoechst 33342, Trihydrochloride, Trihydrate - 10 mg/mL Solution in Water **Thermo Fisher Scientific Catalog #H3570**

 Protease from Bacillus Licheniformis **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P5380**

 DPBS no calcium, no magnesium **Invitrogen - Thermo Fisher Catalog #14190136**

 HBSS **Gibco - Thermo Fisher Scientific Catalog #14060040**

 Quick-Read 10 Chamber Slide **Globe Scientific Catalog #3805**

 EDTA

 NucGreen™ Dead 488 ReadyProbes™ Reagent **Thermo Fisher Scientific Catalog #R37109**

 Quick-Read 10 Chamber Slide **Globe Scientific Catalog #3805**


 Countess™ Cell Counting Chamber Slides **Catalog #C10314**

 21G needle **VWR International (Avantor) Catalog #BD-305165**

 Ammonium Chloride Solution 100 mL **STEMCELL Technologies Inc. Catalog #7800**


 Ammonium Chloride Solution 100 mL **STEMCELL Technologies Inc. Catalog #7800**

 NucGreen™ Dead 488 ReadyProbes™ Reagent **Thermo Fisher Scientific Catalog #R37109**

 Hoechst 33342, Trihydrochloride, Trihydrate - 10 mg/mL Solution in Water **Thermo Fisher Scientific Catalog #H3570**

 Flowmi cell strainer **Catalog #H13680-0040**

Safety warnings

-  Samples coming from patients with undetermined viral status should be process in cell culture rooms with the appropriate safety level.



Before start

Prepare *Bacillus Licheniformis* enzyme mix just prior to starting dissociation:

	Volume (μl)	Reagent	Final concentration
	850	DPBS	1X
	50	20 mM EDTA	0.5 mM
	100	Protease from <i>B. Licheniformis</i> (100 mg/mL)	10 mg/mL

Prepare Inactivation buffer:

Make stock of 10% BSA in HBSS and store at -20 °C.

To make HBSS/BSA 2% aliquot 40 mL of HBSS in 50 mL conical and pipet in 10 mL of 10% BSA stock.

Prepare Wash buffer:

To make HBSS/BSA 1% aliquot 20 mL of HBSS in 50 mL conical and pipet in 20 mL of HBSS/BSA 2%.

Prepare Resuspension buffer:

To make HBSS/BSA 0.05% aliquot 1 mL of HBSS/BSA 2% in 50 mL conical and pipet in 39 mL of HBSS.

Prepare cell staining reagent:

- HBSS: 500 μL
- Hoechst 33342 (10 mg/mL): 1 μL
- NucGreen™ Dead 488 ReadyProbes™ Reagent: 1 drop



- 1 Perform bronchial biopsy at the desired level of the airways (to be performed by a medical doctor)

Equipment

Biopsy forceps

NAME

Medi-Globe

BRAND

GBF-21-18-120

SKU

- 2 Put the biopsy in 1 mL DPBS in a well of a 6-well plate, observe aspect, and then transfer into 1 mL of ice-cold dissociation buffer in a 1.5 mL eppendorf tube. Use wide-bore 1 mL pipet tip for all biopsy transfers.

 1 mL

Note

PREPARATION OF DISSOCIATION MIX (Fresh at each experiment)

Ingredients:

- PBS
- Protease from *Bacillus Licheniformis* (100 mg/mL stock solution in PBS)
- EDTA 10 mM

For 1 mL of dissociation mix add:

- 850 microlitres of PBS
- 100 microlitres of protease (Final concentration: 10 mg/mL)
- 50 microlitres of EDTA (Final concentration: 0.5 mM)

 4 °C

Expected result




- 3 If transportation or storage is necessary: place tube on ice or in polystyrene box containing ice packs. Biopsies can be stored for 60 min in dissociation buffer.
- 4 Carefully remove biopsy from the dissociation buffer, with a wide-bore 1 mL pipet tip and place in a 100 mm petri dish, taking as little dissociation buffer as possible. Mince with a scalpel equipped of a 10 blade. Drag the biopsy out of the liquid and mince very carefully into the smallest possible pieces. With a wide-bore pipet tip, transfer back the minced biopsy with a small volume of dissociation buffer. Rinse the petri dish with dissociation buffer, at the location of biopsy mincing to recover as many cells as possible.
- 5 Incubate cells on ice for 90 to 120 min after mincing, with gentle trituration with needles 5 times every 5 min. Use needle with decreasing sizes from 21G to 23G.

 01:30:00 Incubation

 00:05:00 Trituration

 4 °C

- 6 Inactivate protease by adding 200 μ L of Inactivation buffer (HBSS containing 2% BSA)


 200 μ L Inactivation buffer

Note

Prepare Inactivation buffer:

HBSS : 40 mL

10% BSA stock: 10 mL

- 7 Spin at 400g for 5 min at 4°C
- 8 Discard supernatant leaving 10 μ L of residual liquid on the pellet.
- 9 Resuspend in 100 μ L of wash buffer (HBSS + 1% BSA)
 100 μ L wash buffer

Note**Prepare Wash buffer:**

HBSS : 20 mL

HBSS/BSA 2%: 20 mL

- 10 Observe cells under an inverted microscope to evaluate red blood cells (RBC) content. RBC content is better evaluated using an automated cell counter such as Countess, after addition of Hoechst 33342 to an aliquot of the cell suspension to discriminate nucleated cells from non-nucleated cells.

- 11 If RBC content is lower than 50%, go directly to step 18.
Perform RBC lysis: add 900 μ L of Ammonium Chloride 0.8% to 100 μ L of cell suspension (9 volumes).

Ammonium Chloride Solution 100 mL **STEMCELL Technologies Inc. Catalog #7800**900 μ L Ammonium Chloride 9 vol. for 1 cell vol.

- 12 Incubate on ice for 10 min.



00:10:00 RBC lysis



4 °C

- 13 Add 200 μ L of Inactivation buffer

200 μ L Inactivation buffer

- 14 Spin at 400g for 5 min at 4°C

- 15 Discard supernatant leaving 10 μ L of residual liquid on the pellet.

- 16 Resuspend in 100 μ L of wash buffer and monitor correct RBC lysis under microscope





100 μ L wash buffer

- 17 Add 1 mL of wash buffer



1 mL wash buffer





- 18 Spin at 400g for 5 min at 4°C
- 19 Discard supernatant leaving 10 µL of residual liquid on the pellet.
- 20 Resuspend in 500 µL of wash buffer
 500 µL wash buffer
- 21 Filter cell suspension through Flowmi cell strainer
 Flowmi cell strainer **Catalog #H13680-0040**
<https://www.youtube.com/watch?v=taS1BuTnvds>
- 22 Add 500 µL of wash buffer to filtered cells.
 500 µL wash buffer
- 23 Discard supernatant leaving 10 µL of residual liquid on the pellet.
- 24 Resuspend in 40 µL of resuspension buffer (HBSS + 0.05% BSA).
 40 µL Resuspension buffer

Note

Prepare Resuspension buffer:

HBSS : 39 mL


HBSS/BSA 2%: 1 mL

- 25 Mix 10 µL of cells with 10 µL of cell counting solution (HBSS with 20 µg/mL Hoechst 33342 and NucGreen). Incubate for 1 min at room temperature.
 Hoechst 33342, Trihydrochloride, Trihydrate - 10 mg/mL Solution in Water **Thermo Fisher Scientific Catalog #H3570**
 NucGreen™ Dead 488 ReadyProbes™ Reagent **Thermo Fisher Scientific Catalog #R37109**

Note

Preparation of cell staining reagent:

- HBSS: 500 μ L
- Hoechst 33342 (10 mg/mL): 1 μ L
- NucGreen™ Dead 488 ReadyProbes™ Reagent: 1 drop

 00:01:00

- 26 Count with Countess automated cell counter using both sides of chambers. Monitor the percentage of nucleated cells (Hoechst +) and dead cells (GFP+).

Equipment

new equipment

NAME

Thermo Fisher Scientific

BRAND


AMQAF1000

SKU


Countess™ II FL Automated Cell Counter with Dapi and GFP cubes

SPECIFICATIONS

Expected result

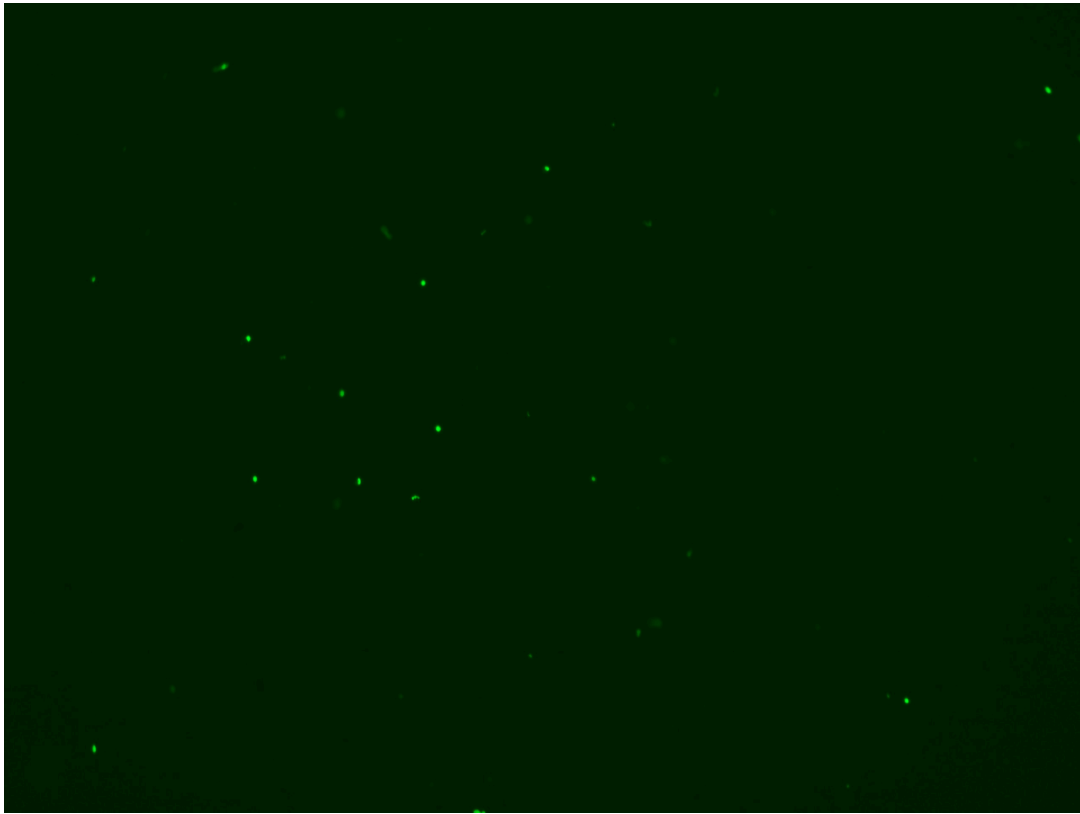


Countess GFP image after NucGreen and Hoescht33342 staining

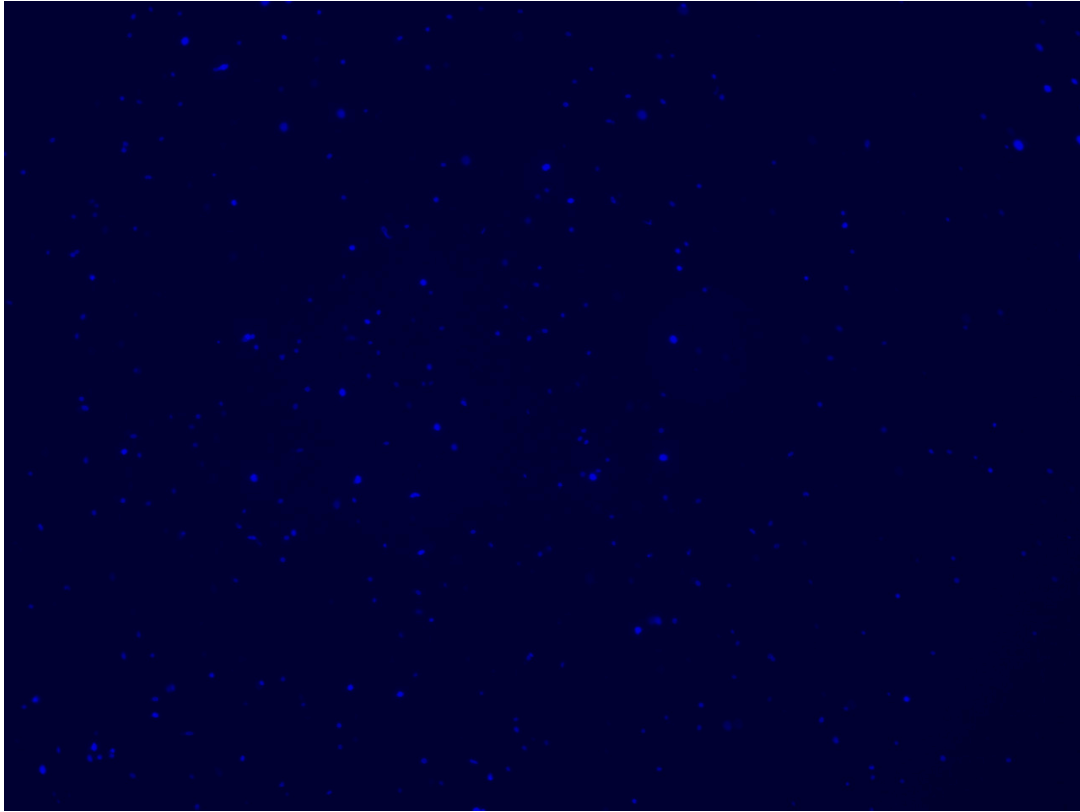

Countess Dapi image after NucGreen and Hoescht33342 staining




Countess report after NucGreen and Hoescht33342 staining

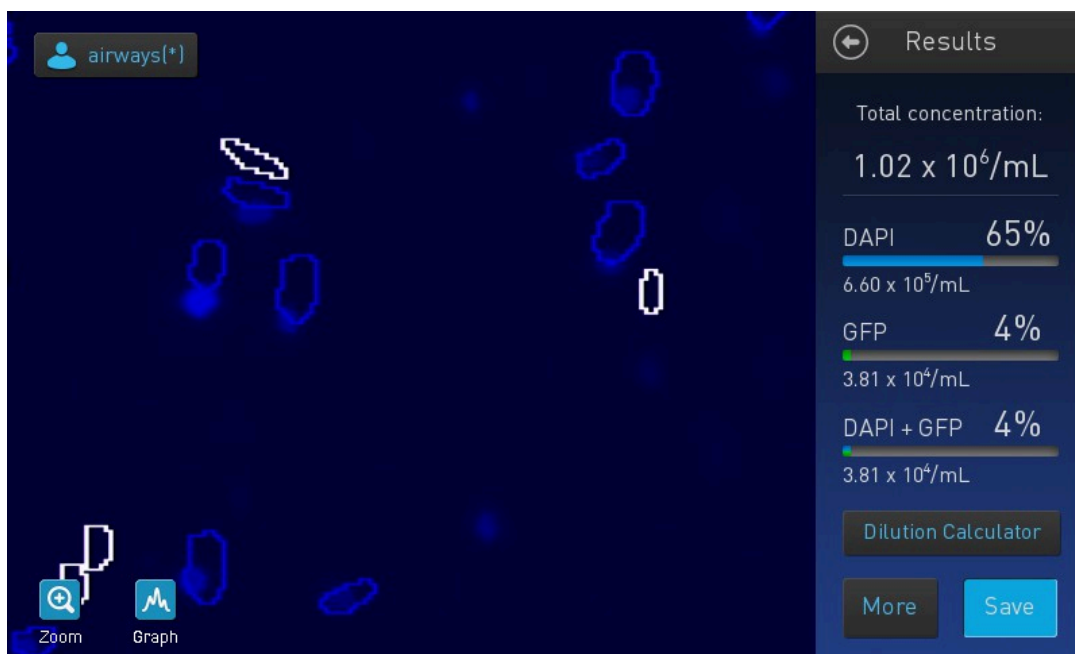

Countess report after NucGreen and Hoescht33342 staining

Expected result

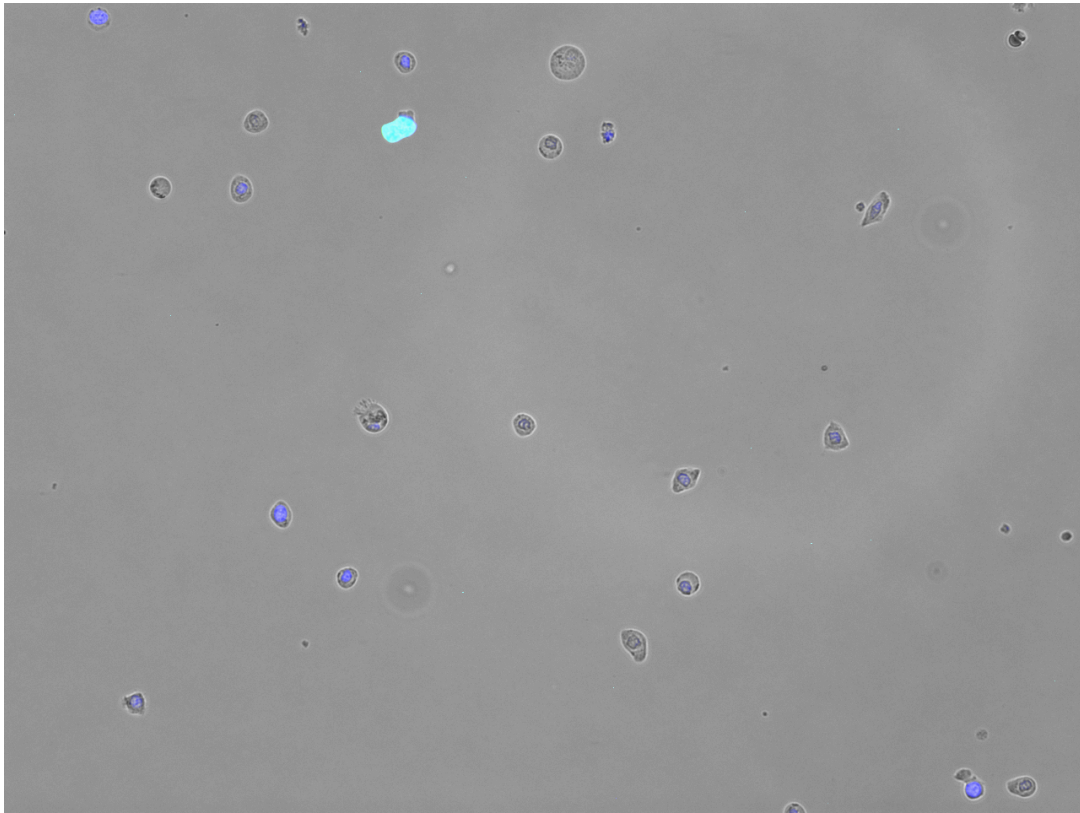

Fluorescence image after NucGreen and Hoescht33342 staining

- 27 Adjust concentration to a range of 700 to 1000 cells/ μ L (with wash buffer) for 10X Chromium. Monitor final cell concentration.