

🛡️ Cell Fixation and Permeabilization Protocol using 70% Ethanol V.3

DOI

dx.doi.org/10.17504/protocols.io.bacziax6



Sam Li¹

¹BioLegend

BioLegend

Tech. support email: tech@biolegend.com



Sam Li

BioLegend



DOI: dx.doi.org/10.17504/protocols.io.bacziax6

External link: <https://www.biolegend.com/protocols/cell-fixation-and-permeabilization-protocol-using-70-ethanol/4252/>

Protocol Citation: Sam Li . Cell Fixation and Permeabilization Protocol using 70% Ethanol. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bacziax6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: December 11, 2019


Last Modified: December 11, 2019

Protocol Integer ID: 30841

Keywords: fix perm, ethanol, flow cytometry

Materials

MATERIALS

 Cell Staining Buffer **BioLegend Catalog #420201**



- 1 Prepare 70% Ethanol and chill to -20°C. Tip: Do not freeze ethanol for long-term storage.
- 2 Prepare target cells of interest and wash 2X with PBS, centrifuging at 350xg for 5 minutes
- 3 Discard supernatant and loosen the cell pellet by vortexing.
- 4 Add 3ml cold 70% ethanol drop by drop to the cell pellet while vortexing.
- 5 Continue vortexing for 30 seconds and then incubate at -20°C for 1 hour.
- 6 Wash 2X with BioLegend Cell Staining Buffer (Cat. [420201](#)) and resuspend cells at 0.5- 1.0×10^7 cells/ml.
- 7 Use 100µl cell suspension/staining tube.
Note: Please note that certain markers or fluors may not survive ethanol fixation. Protein-based fluors, like PE and APC, tend to have more difficulty, while synthetic fluors, like Brilliant Violet™, tend to have a higher chance of surviving the process.