

Apr 15, 2025

Version 2

Chloroform-free DNA Extraction - Ammonium Acetate Precipitation Method V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9pekzg3e/v2>

NEOF - NERC Environmental Omics Facility VF¹

¹University of Sheffield

NEOF - NERC Environmental Omics Facility VF: School of Biosciences, Western Bank, University of Sheffield, Sheffield, S10 2TN, UK



NEOF - NERC Environmental Omics Facility VF

University of Sheffield

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9pekzg3e/v2>

Protocol Citation: NEOF - NERC Environmental Omics Facility VF 2025. Chloroform-free DNA Extraction - Ammonium Acetate Precipitation Method. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x54v9pekzg3e/v2> Version created by **NEOF - NERC Environmental Omics Facility VF**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: April 15, 2025

Last Modified: April 15, 2025

Protocol Integer ID: 126730

Keywords: dna extraction, cheap, cost-effective, non hazardous, ammonium acetate, blood, tissue, feather, hair, insect, quick, effective dna extraction method, chloroform, ammonium acetate precipitation method, extraction

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISKThe protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services. We accept no responsibility or liability for materials damaged or destroyed nor costs incurred through the use of this protocol

Abstract

A chloroform-free, cost-effective DNA extraction method for a variety of sample types.



Materials

Digsol: To make 500ml

- 20ml 0.5M EDTA (pH 8.0)
- 3.425g NaCl
- 25ml 1M Tris-HCl (pH 8.0)
- 430ml ddH₂O

Autoclave then add 25ml of 20% SDS.

Low TE (Tris_{10mM}, EDTA_{0.1mM}): To make 500ml

- 5ml 1M Tris-HCl (pH 8.0)
- 100µl 0.5M EDTA (pH8.0)
- 495ml ddH₂O

Autoclave.

Protocol materials

⊗ Proteinase K, 2mL **Qiagen Catalog #19131**

⊗ Ammonium Acetate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A1542-500G**

⊗ 1M DTT **Merck MilliporeSigma (Sigma-Aldrich) Catalog #43816**

Safety warnings

- ! Use GLP and wear protective equipment. Avoid contact with skin.
All chemicals can be disposed of down the sink with copious amounts of water to dilute the ethanol down to >20% of the waste.

If inhaled: If unconscious, place in recovery position and seek medical advice. Keep the respiratory tract clear. If symptoms persist, call a physician.

In case of skin contact: Wash off immediately with soap and plenty of water while removing all contaminated clothes and shoes. If symptoms persist, call a physician.

In case of eye contact: Remove contact lenses. Protect unharmed eye. Rinse thoroughly with plenty of water for at least 15 minutes and consult a physician.

If swallowed: If accidentally swallowed, obtain immediate medical attention. Rinse mouth with water. Never give anything by mouth to an unconscious person.



- 1 Add  250 μL Digsol buffer (see Materials for recipe) and  10 μL  Proteinase K, 2mL **Qiagen Catalog #19131** to a labelled 1.5ml tube. 1m
- 2 **Tissue:** Cut into small pieces ($<1\text{cm}^2$) with a sterile razor blade on a sterile glass plate before adding to the 1.5mL tube. 5m
Blood in ethanol: Centrifuge blood sample at 13,000rpm for about 1 min (to pellet sample).
Remove sample from ethanol with toothpick and blot onto tissue. When almost dry, transfer the toothpick into the 1.5mL tube and jiggle to dislodge the blood. Remove toothpick and place in disinfectant.
Swab in ethanol: Dry of ethanol and place into the 1.5mL tube. Snap off the end of the swab so the lid may close.
Feather: Cut the calamus of 1-3 feathers into small pieces with a sterile razor blade on a sterile glass plate before adding to the tube. If feathers are very small and blood spots are present on the tips, feathers can be added whole.
Hair: Place hair(s) into the 1.5mL tube preferably with the largest root and add  5 μL  1M DTT **Merck MilliporeSigma (Sigma-Aldrich) Catalog #43816**
Insect: Degut and add to 1.5mL tube. If the insect is very small, starve for 48hrs before freezing. Add  5 μL  1M DTT **Merck MilliporeSigma (Sigma-Aldrich) Catalog #43816** . Crush with a pestle or add a metal lysing bead and place on TissueLyser (Qiagen) for 2min, 30/s.
- 3 Vortex and place samples in a rotating oven at  55 $^{\circ}\text{C}$ for  03:00:00 or  Overnight for maximum digestion. 3h
- 4 Add  300 μL 4M  Ammonium Acetate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A1542-500G** . 1m
- 5 Vortex several times over a period of at least  00:15:00 at  Room temperature . 15m
- 6  13000 rpm, 00:10:00 10m



- 7 Aspirate the supernatant (clear liquid containing the DNA) into a new labelled 1.5ml tube. Discard tube containing the pelleted protein debris. 1m
- 8 Add  1 mL 100% ethanol and invert the tube gently several times to precipitate DNA. 1m
- 9  13000 rpm, 00:10:00 10m
- 10 Pour off ethanol in a smooth motion, taking care not to lose the DNA pellet. 1m
- 11 Add  500 μ L 70% ethanol and invert gently several times to clean the pellet. 1m
- 12  15000 rpm, 00:05:00 5m
- 13 Pour off ethanol in a smooth movement or using a pipette gently draw off the supernatant if fear of losing the pellet. Stand tubes upside-down on clean tissue until dry (approx. **30-60 minutes**). This can be sped up by using the heat of a lamp from above. 30m
- 14 Once fully dry add approx.  100 μ L Low TE (see Materials for recipe). Add less if a very tiny pellet or no pellet is observed. Flick sample to dislodge pellet. 1m
- 15 Place tubes in a thermomixer or shaking oven at  50 °C for  00:30:00 to dissolve the pellet. If the pellet has not completely dissolved add more Low TE. 30m
- 16 Store at  -20 °C (long term) or  4 °C (short term).



Protocol references

<https://onlinelibrary.wiley.com/doi/epdf/10.1046/j.0962-1083.2001.01355.x>

Richardson DS, Jury FL, Blaakmeer K, Komdeur J, Burke T (2001) Parentage assignment and extra-group paternity in a cooperative breeder: the Seychelles warbler (*Acrocephalus sechellensis*). *Molecular Ecology***10**, 2263-2273.

<https://onlinelibrary.wiley.com/doi/abs/10.1034/j.1600-048X.2000.310208.x>

The Evolution of Cooperative and Pair Breeding in Thornbills *Acanthiza* (Pardalotidae)

James A. Nicholls, Michael C. Double, David M. Rowell and Robert D. Magrath

Journal of Avian Biology

Vol. 31, No. 2 (Jun., 2000), pp. 165-176