



Jan 18, 2024

Version 2

Cleavage Under Targets and Release Using Nuclease (CUT&RUN) V.2

 [PLOS One](#)

✓ Peer-reviewed method

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9p8ppg3e/v2>

Geneva Miller¹, Lindsey M. Rollosso¹, Carrie Saada¹, Serenity J. Wade¹, Danae Schulz¹

¹Harvey Mudd College, Claremont, CA

PLOS ONE Lab Protocols

Spotlight series



Danae Schulz

Harvey Mudd College

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9p8ppg3e/v2>

External link: <https://doi.org/10.1371/journal.pone.0292784>

Protocol Citation: Geneva Miller, Lindsey M. Rollosso, Carrie Saada, Serenity J. Wade, Danae Schulz 2024. Cleavage Under Targets and Release Using Nuclease (CUT&RUN). **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.x54v9p8ppg3e/v2> Version created by **Danae Schulz**

Manuscript citation:

Miller G, Rollosso LM, Saada C, Wade SJ, Schulz D (2023) Adaptation of CUT&RUN for use in African trypanosomes. PLOS ONE 18(11): e0292784. <https://doi.org/10.1371/journal.pone.0292784>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 10, 2024

Last Modified: January 18, 2024

Protocol Integer ID: 93219

Keywords: Trypanosoma brucei, CUT&RUN, genomic occupancy, mapping genomic protein binding sites, interest in the protozoan parasite trypanosoma brucei, protozoan parasite trypanosoma brucei, protocol for parasite, epitope tag for the protein, parasite, cleavage under target, epitope tag, using nuclease, native protein, protein, nuclease

Abstract

This Cleavage Under Targets and Release Using Nuclease (CUT&RUN) protocol produces genomic occupancy data for a protein of interest in the protozoan parasite *Trypanosoma brucei*. The data produced is analyzed in a similar way as that produced by ChIP-seq. While we describe the protocol for parasites carrying an epitope tag for the protein of interest, antibodies against the native protein could be used for the same purpose.

The **last step** contains a supplemental video with extra context and tips, as part of the protocols.io Spotlight series, featuring conversations with protocol authors.

Attachments



[720-1571.docx](#)

27KB



Guidelines

Guidelines

Pay extra attention to time-sensitive incubation periods, particularly the DNA cleavage step. We find that working through the protocol rather speedily produces the best results, so don't dawdle.

Additional Notes

We keep our centrifuges at  10 °C to make sure the samples don't freeze if the temperature drifts slightly. All spin steps of the protocol can be performed anywhere between  4 °C and  10 °C .

While we have not rigorously tested whether increasing the exposure time to saponin influences permeabilization, we recommend that the saponin permeabilization step be completed in smaller batches of samples if large numbers of samples (>10) are being processed simultaneously.



Materials

Materials

1.  Spermidine trihydrochloride **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S2501**
2.  Saponin **Merck Millipore (EMD Millipore) Catalog #558255**
3.  Sodium Chloride **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S9888**
4.  Calcium chloride **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C4901**
5.  Tris Hydrochloride (Tris-HCl) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #RES3098T-B7**
6.  cOmplete™ EDTA-free Protease Inhibitor Cocktail **Merck MilliporeSigma (Sigma-Aldrich) Catalog #11873580001**
7.  Ultrapure 0.5M EDTA pH 8.0 **Invitrogen - Thermo Fisher Catalog #15575020**
8.  Bioworld EGTA Buffer 0.5M pH 8.0 100ml **Fisher Scientific Catalog #50-255-957**
9.  mouse anti-rabbit IgG-PE **Santa Cruz Biotechnology Catalog #sc-3753**
10.  Anti-HA antibody produced in rabbit **Merck MilliporeSigma (Sigma-Aldrich) Catalog #H6908** , primary antibody will vary depending on your system
11. Rabbit anti-H3 (gift)
12.  Sodium dodecyl sulfate solution **Merck MilliporeSigma (Sigma-Aldrich) Catalog #428018**
13.  Proteinase K **Thermo Fisher Scientific Catalog #EO0491**
14.  Monarch RNase A – 1 ml (2×0.5ml) **New England Biolabs Catalog #T3018L**
15.  Ampure XP beads **Beckman Coulter Catalog #A63881**
16.  CUT&RUN pAG-MNase and Spike-In DNA **Cell Signaling Technology Catalog #40366**

Equipment

1. Microcentrifuge (Eppendorf 5424R)
2. Swinging bucket centrifuge (Eppendorf 5920R)
3. 1.7ml Olympus microtubes (Genesee 22-281)
4. Olympus 50ml Centrifuge Tubes (Genesee 21-108)
5. Roto-Mini Plus Variable Speed Rotator with tube holders, 115V (ThermoFisher Scientific, 1159P34)
6. Novocyte Flow Cytometer 2000

Buffers

NP-S Buffer with 0.1% Saponin



	A	B
	Individual Components	
	spermidine	0.5 mM
	Saponin	0.1% (vol/vol)
	NaCl	50 mM
	Tris-Cl (pH 7.5)	10 mM
	Store at 4°C	
	Add protease inhibitors just before use.	

NP-S Buffer no detergent

	A	B
	Individual Components	
	spermidine	0.5 mM
	NaCl	50 mM
	Tris-Cl (pH 7.5)	10 mM
	Store at 4°C	
	Add protease inhibitors just before use.	

Antibody Buffer**NP-S Buffer** with 2mM EDTA**2X Stop Buffer**

	A	B
	EDTA	20mM
	EGTA	40mM
	Add 50 pg of yeast spike-in DNA to each reaction, or 50 pg/100 µl of 2X Stop Buffer.	



Before start

Before Start

We use 50-75 million bloodstream parasites per sample. Parasites are cultured in HMI9 media with incubation at  37 °C and 5% CO₂. Cultures should be prepared in advance so that sufficient numbers of parasites are available for each sample.

The protocol works best when everything is kept cold prior to cutting with the protein A-MNase fusion protein. We recommend keeping buffers chilled on ice and pre-cooling centrifuges to  4-10 °C . Protease inhibitors should be added to the NP-S buffer just before commencing the experiment. The amount of **2X Stop Buffer** required for the experiment should be calculated and yeast spike-in DNA should be added prior to starting (see 2X Stop Buffer recipe below).



Prepare cells

1 Count parasite cultures with hemocytometer or other preferred counting method.

2 Spin down cells in centrifuge at  2800 x g, 00:10:00 .

10m

Note

This spin step can be performed at  Room temperature or at  10 °C .



2.1 Remove supernatant and resuspend in small amount of remaining media (~  100 µL).

3 If needed, combine samples from multiple Eppendorf tubes so that each final tube has 75 million cells, and spin again at  2800 x g, 10°C, 00:04:00 in microcentrifuge. Remove supernatant.

4m



Permeabilize cells

4 Wash all samples with  1 mL **NP-S Buffer with 0.1% Saponin.**



5 Spin at  4600 x g, 10°C, 00:04:00 . Remove sup.

4m



Primary Antibody Binding

6 Resuspend each sample in  100 µL **NP-S Buffer with 0.1% Saponin.**

7 Add EDTA to  2 millimolar (mM) final ( 4 µL of  0.05 Mass Percent EDTA/ tube).



8 Add  5 µg α-HA antibody (or other antibody against protein of interest for experimental sample) OR add  5 µg α-H3 antibody (for control sample, might require





titration) OR add  5 μ g α -IgG antibody (for control sample).

9 Incubate for  00:45:00 on tube rotator at  Room temperature .

45m



10 Add  1 mL NP-S Buffer no detergent to each sample.



11 Spin at  4600 x g, 10°C, 00:04:00 . Remove supernatant.

4m



12 Add  1 mL NP-S Buffer no detergent to each sample.



13 Take out aliquot of 1 million cells for flow cytometry analysis (~  14 μ L if starting with 75 million cells). Set aside  On ice .

14 Spin the remainder of the sample at  4600 x g, 10°C, 00:04:00 . Remove supernatant. (total 2 washes).

4m



pAG-MNase binding

15 Resuspend each sample in  100 μ L of buffer **NP-S Buffer no detergent**.

16 Add  1.5 μ L of pAG-MNase enzyme to each sample.



17 Incubate for  00:45:00 at  Room temperature on rotator.

45m



FACS Sample preparation

18 Add  1 mL **NP-S Buffer no detergent** to 1 million cell aliquot prepared above.



**Note**

We also routinely use HMI-9 media for the flow cytometry antibody incubation and washes.

- 19 Spin at  4600 x g, 00:04:00 . 4m 
- 20 Resuspend in  100 μ L NP-S Buffer no detergent.
- 21 Stain with α -rabbit IgG PE at 1:200 for  00:15:00 at  Room temperature . 15m
- 22 Wash. 
- 22.1 Wash in  1 mL NP-S Buffer no detergent (or HMI-9) at  7000 rpm, 00:04:00 . 4m
- (1/2)
- 22.2 Wash in  1 mL NP-S Buffer no detergent (or HMI-9) at  7000 rpm, 00:04:00 . 4m
- (2/2)
- 23 Resuspend in  300 μ L NP-S Buffer no detergent (or HMI-9).
- 24 Transfer sample into flow cytometry tube.
- 25 Analyze on flow cytometer.

pAG-MNase wash

- 26 

**Note**

continued after 45m incubation above.

Add  1 mL **NP-S Buffer no detergent** to each sample.

27 Spin at  4600 x g, 10°C, 00:04:00 . Remove supernatant.

4m



28 Add  1 mL **NP-S Buffer no detergent** to each sample.



29 Spin at  4600 x g, 10°C, 00:04:00 . Remove supernatant.

4m

**Targeted Digestion Preparation**

30 Make sure to prepare enough **2X Stopbuffer** with spike in control.

Note

This is the section where timing is very important.

Targeted Digestion

31 Resuspend each sample in  100 µL buffer **NP-S Buffer no detergent**.

32 Incubate at  Room temperature for  00:05:00 .

5m



33 Add  2 µL  100 millimolar (mM) CaCl_2 to all samples (final concentration = 2mM), mix by flicking.



34 Incubate all samples at  25 °C for  00:05:00 .

5m



**Note**

We recommend using a heat block as ambient temperature in the lab can vary.

35 Add  100 μL 2X Stop buffer to each sample and mix by flicking.

**Chromatin Release**

36 Incubate  00:10:00 at  37 °C to release CUT&RUN fragments from the insoluble nuclear chromatin.

10m



37 Spin at  4600 x g, 10°C, 00:04:00 . Remove supernatant into new tube. SAVE SUPERNATANT.

4m

**DNA Extraction**

10m

38 To all samples add  2 μL of 10% SDS (final concentration = 0.1%),  3.3 μL of  10 μL proteinase K (165 $\mu\text{g}/\text{ml}$), and  1.33 μL of  1 μL RNase A (6.5 $\mu\text{g}/\text{ml}$).



39 Mix by gentle flicking and incubate for  00:10:00 at  70 °C .

10m



40 Purify using Ampure XP beads at 1.8X or phenol chloroform extraction.

Note

Following DNA purification, high-throughput sequencing libraries can be prepared using the preferred method of the research lab.

Spotlight video

41

<https://www.youtube.com/embed/8ray2eO8zJE?si=AFat8OzneyCF5ZQK>



Protocol references

Citations

Ashby E, Paddock L, Betts HL, Liao J, Miller G, Porter A, Rollosso LM, Saada C, Tang E, Wade SJ, Hardin J, Schulz D. 2022. Genomic occupancy of the bromodomain protein Bdf3 is dynamic during differentiation of African trypanosomes from bloodstream to procyclic forms. *mSphere* 7:e00023-22.

Skene PJ, Henikoff S. 2017. An efficient targeted nuclease strategy for high-resolution mapping of DNA binding sites. *eLife* 6:e21856.